

# Application News

## No. B71

### MALDI-TOF Mass Spectrometry

## Molecular Weight Measurement of Glycoproteins Using a Benchtop MALDI-TOF Mass Spectrometer

MALDI-TOF mass spectrometers are a mass spectrometer type that is used in a wide range of fields likewise LCMS, in terms of high through-put and high sensitivity. These instruments are being utilized more and more for simple molecular weight measurement and profiling of synthesized products and high-molecular compounds. This is because instruments of this type have several features: singly-charged ions are generated so molecular weights can be recognized easily, the applicable mass range is wide, and there are many solvent options because the sample is dried before measurement.

On the other hand, due to changes in social conditions in these several years, government offices and private enterprises strongly request the reduction of costs for both introduction and running of instruments used for such applications. The benchtop MALDI-TOF mass spectrometer "MALDI-8020" is a new instrument that can sufficiently meet such market needs. The noteworthy point of this instrument is that it has a shorter flight tube, which is the key feature of its small size, while retaining the performance equal to or higher than that of a conventional model.

Molecular weight measurement of proteins modulated by post-translation modification using a MALDI-TOF mass spectrometer is the most basic application in the fields of biochemistry and molecular biology. Molecular weights of glycoproteins, in particular, are observed as a molecular weight distribution of multiple components because they are heterogeneous due to the difference in their sugar chain structures.

This article introduces an example of measuring molecular weights of such glycoproteins using the benchtop MALDI-TOF mass spectrometer "MALDI-8020".

Y. Yamazaki

### ■ Benchtop MALDI-TOF MS: MALDI-8020

The MALDI-8020 (Fig. 1) is a compact-design, minimal-space linear mode-only MALDI-TOF mass spectrometer. Its performance in linear mode (positive ion) is comparable to the same mode of a conventional MALDI-TOF mass spectrometer. Equipped with a 200 Hz solid-state laser and a load-lock chamber mechanism that enables a target plate change while maintaining the degree of vacuum at the measuring position, the instrument ensures rapid measurements.

### ■ Example of Glycoprotein Molecular Weight Measurement

Fig. 2 shows the amino acid sequence of Ribonuclease B. It is known that due to the N-linked glycosylation of the high mannose type, this protein is detected with its main component being a modification approx. 1217 greater than its intrinsic mass. Bovine derived ribonuclease B (5 pmol) was mixed with ferulic acid (10 mg/mL, 50 % acetonitrile/0.1 % formic acid solution), and analyzed in linear mode using MALDI-8020. The result is shown in Fig. 3. The singly-charged protonated molecule of the detected ribonuclease B is detected at  $m/z$  14890, which is consistent with the theoretical value (14899.3). In addition, the distribution due to heterogeneous sugar chains is detected as a mass difference (theoretical value: 162) derived from hexose. As indicated above, the resolving power, mass accuracy, etc., of the linear mode of the MALDI-8020 are comparable to those of a larger conventional MALDI-TOF mass spectrometer in the same mode.



Fig. 1 Benchtop MALDI-TOF MS: MALDI-8020

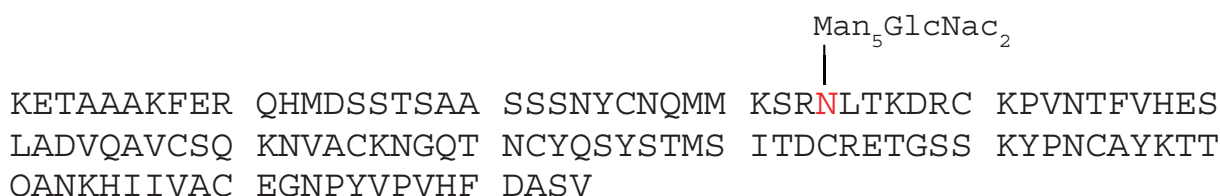


Fig. 2 Amino Acid Sequence and Glycosylation of Ribonuclease B

## Conclusion

This example demonstrate that the benchtop MALDI-TOF mass spectrometer "MALDI-8020" can be used for measuring molecular weights of glycoproteins and detecting the distribution originating from glycosylation at the same level as larger-sized MALDI-TOF mass spectrometers.

In recent years, MALDI-TOF mass spectrometers have evolved into high specification instruments and have increased in size and complexity. This results in over-specifications against the needs of measurement in linear mode, which occupies the most of MALDI applications. Size and initial/running costs also pose a barrier for introducing the instrument for applications requiring only modest performance.

The MALDI-8020, the world's smallest linear-only MALDI-TOF mass spectrometer, satisfies the needs of protein analysis in linear mode, and its future dissemination in biochemistry and molecular biology is expected.

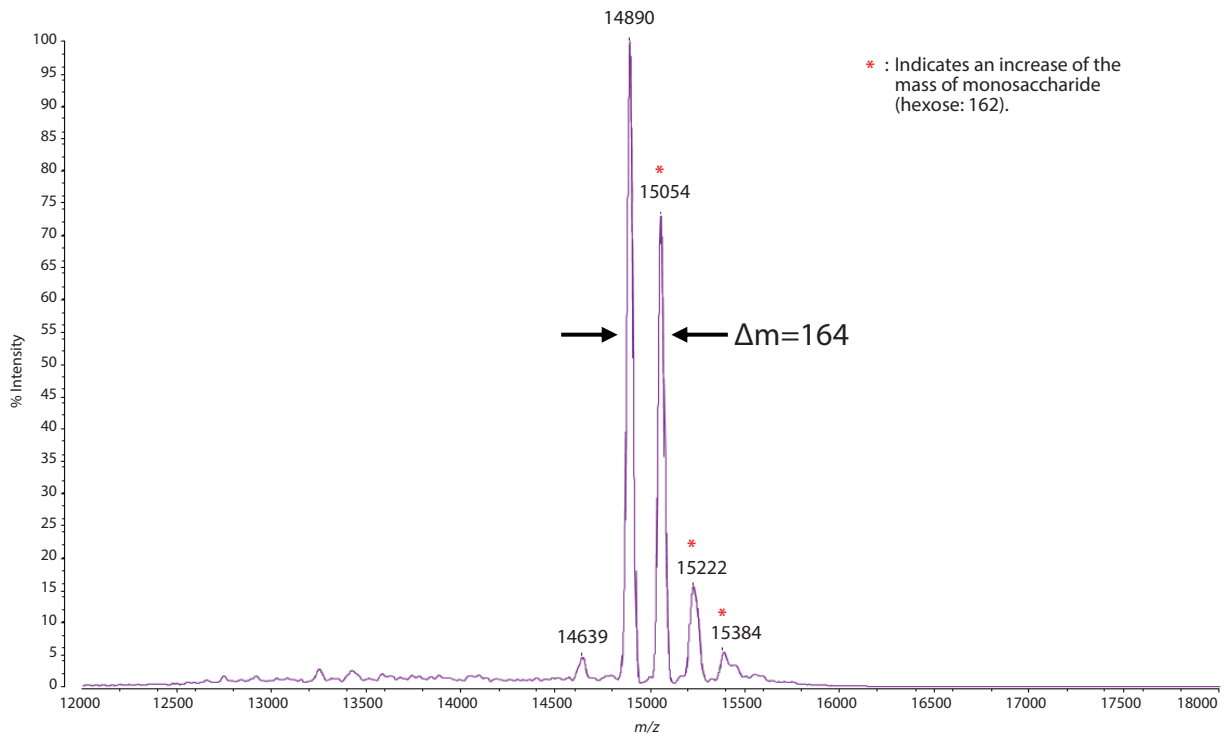


Fig. 3 Mass Spectrum of Ribonuclease B

First Edition: Nov. 2017



For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. Shimadzu disclaims any proprietary interest in trademarks and trade names used in this publication other than its own. See <http://www.shimadzu.com/about/trademarks/index.html> for details.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

Shimadzu Corporation

[www.shimadzu.com/an/](http://www.shimadzu.com/an/)