


# Settings for copying MS images

MS images can be copied from the “MS image” panel


The screenshot displays the software interface with several panels. The 'MS image' panel on the right shows a heatmap with a white square highlighting a specific region. The 'Data Matrix Table' in the center shows a table of data points. The 'MS image list' panel at the bottom right shows a list of MS images. The 'Analysis Parameters' panel at the bottom left shows various settings.

MS	MS	MS	MS	MS	MS	MS	MS	MS	MS
1	796.52130	796.52130	796.52130	796.52130	796.52130	796.52130	796.52130	796.52130	796.52130
2	796.52130	796.52130	796.52130	796.52130	796.52130	796.52130	796.52130	796.52130	796.52130
3	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
4	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
5	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
6	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
7	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
8	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
9	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
10	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
11	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
12	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
13	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
14	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
15	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
16	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
17	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
18	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
19	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445
20	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445	797.03445

The contents to be copied can be selected

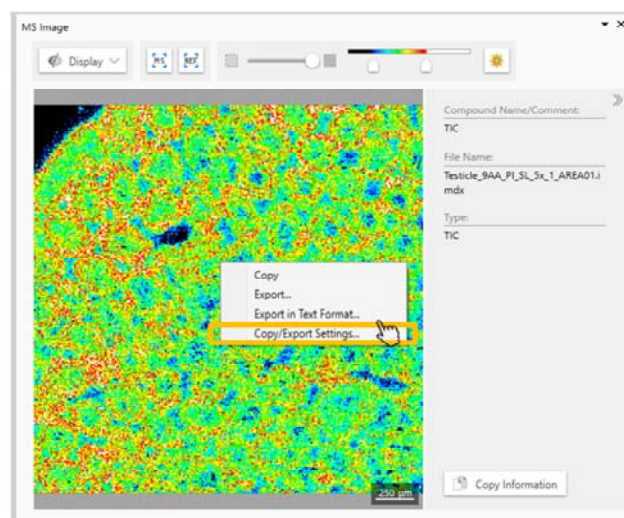

## Various settings can be applied

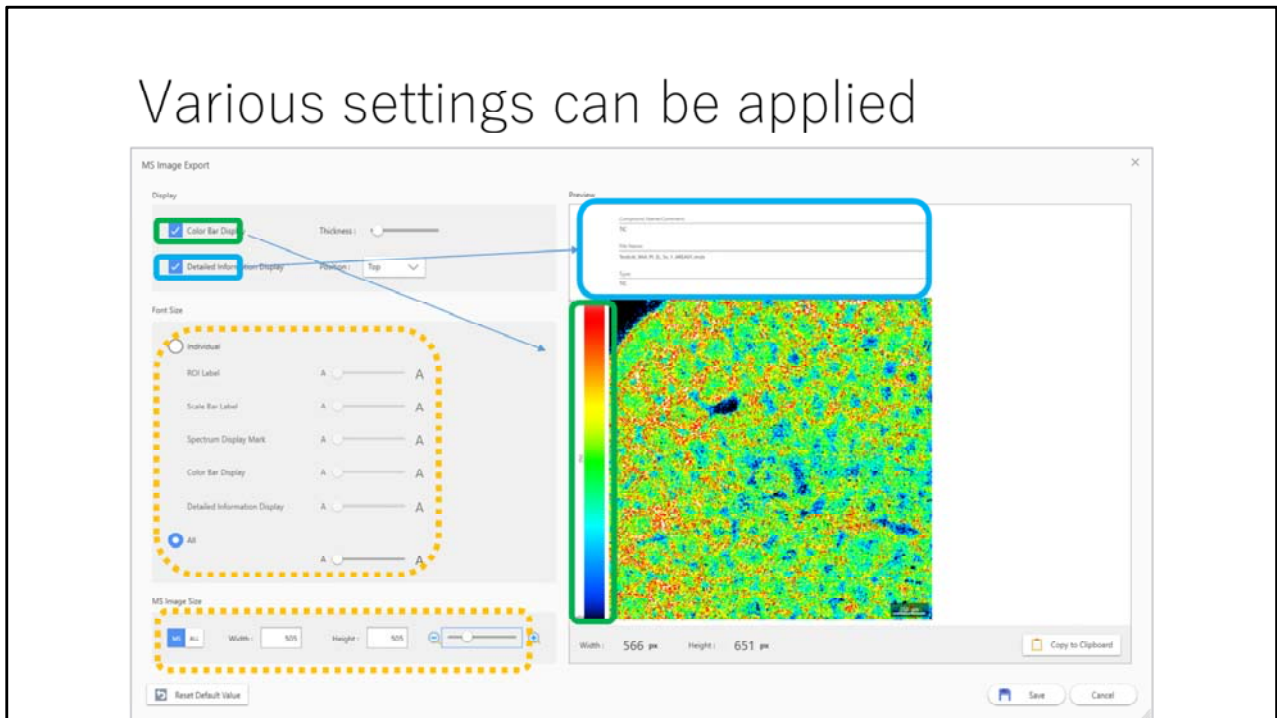

Choose whether to display the colour bar or detailed information.  
The font size of the text and the overall size of the image can also be adjusted.

For the MS image size, the pixel count can be changed and a region can be selected to swap with either the MS image or the whole reference image.  
The scale bar can be displayed on the main image.