

Application Note


No. 48

Life Science

Comprehensive Measurement of Metabolites Using GC-MS/MS and LC-MS/MS

- An Application to the Research of the Intestinal Environment -

Takanari Hattori^{*1,*2}, Akihiro Kunisawa^{*1,*2}, Shuichi Kawana^{*1}, Shin-ichi Kawano^{*1,*2}, Yoshihiro Hayakawa^{*1},
Junko Iida^{*1,*2}, Eiichiro Fukusaki^{*2,*3}, Mitsuharu Matsumoto^{*4}

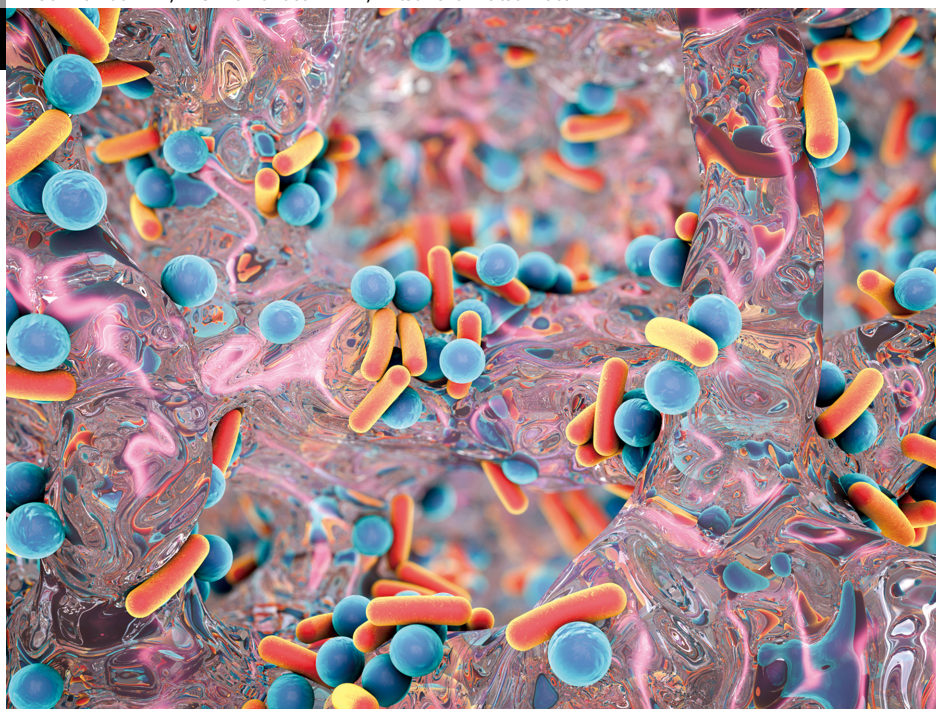

Life Science

1. Introduction

The human intestine harbors over 100 species of bacteria with a total of approximately 100 trillion bacterial cells, creating a complex intestinal microbiota. Studies in recent years have revealed that intestinal bacteria may be beneficiary to the host by contributing to health preservation and promotion. This has raised the attention on researches regarding the intestinal microbiota. In the field of medicine, researches are investigating the relationship of the intestinal microbiota with diseases such as colorectal cancer as well as the effects on drug efficacy and the immune system. In the field of foods, there are studies aiming to verify the health effects of lactic acids in functional foods. As a means to reveal the effects of the intestinal microbiota on the host, next-generation sequencing is used to analyze the species and distribution of bacteria in the microbiota. Another method gaining attention in recent years is metabolomic analysis using a mass spectrometer for a comprehensive analysis of the metabolites produced by the intestinal microbiota. This is because it is considered that some of the metabolites produced by the intestinal microbiota are absorbed by the intestinal tract, taken into

the bloodstream, and then circulated through the entire body.

Metabolomic analysis using a mass spectrometer generally employs a gas chromatograph mass spectrometer (GC-MS) or a liquid chromatography mass spectrometer (LC-MS) to comprehensively measure the metabolites (the metabolome) within a sample. As indicated in Fig. 1, it is necessary to select either a GC-MS or an LC-MS as appropriate for the target compound and the measurement purpose. A GC-MS features excellent robustness, and although derivatization is required for hydrophilic metabolites such as amino acids, organic acids, and sugars, it can comprehensively measure several hundred compounds in one measurement. An LC-MS on the other hand does not require derivatization. Capable of efficiently measuring specific metabolites (up to 100 compounds), it is optimal for the routine measurement of specific compounds.

This article introduces a method for the comprehensive measurement of the primary metabolites in mouse feces using GC-MS/MS and LC-MS/MS.

*1 Analytical & Measuring Instruments Division, Shimadzu Corporation

*2 Osaka University and Shimadzu Analytical Innovation Research Laboratory

*3 Graduate School of Engineering, Osaka University

*4 Research Laboratories, Kyodo Milk Industry Co., Ltd.

2. Pretreatment⁽¹⁾

The pretreatment procedure for extracting metabolites from mouse feces is shown in Fig. 2. Fresh feces were collected from male C57BL/6J mice reared in a normal environment. Taking 50 mg of the feces, we added 450 μ L of phosphate-buffered saline. After mixing, the supernatant was centrifuged and subjected to ultrafiltration. For the measurement of primary metabolites using GC-MS/MS, the resulting filtrate is subjected to derivatization and is then used as the measurement sample. Details of the derivatization method are described in the "Pretreatment Procedure Handbook for Metabolites Analysis" (C146-E323). For the measurement of primary metabolites using LC-MS/MS, the measurement sample was prepared by diluting the filtrate by a factor of 10.

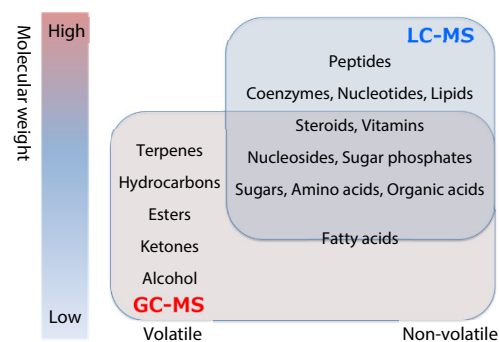


Fig. 1 Target Compounds for GC-MS and LC-MS

3. Analytical Conditions

GC-MS/MS analysis was done using the MRM method included in the Smart Metabolites Database™ together with the GCMS-TQ™ 8040. The Smart Metabolites Database contains the information of MRM transitions of 475 compounds, mainly of metabolites in biological samples, and allows the simultaneous measurement of 475 metabolites. Details are described in Application Data Sheet No. 104. For LC-MS/MS analysis, we used the ion pairing LC-MS/MS method and non ion-pairing LC-MS/MS method included in the LC/MS/MS Method Package for Primary Metabolites (ver. 2) together with the LCMS-8040 and the LCMS-8050. The ion pairing LC-MS/MS method makes possible the simultaneous measurement of 55 metabolites that are highly important for metabolomic analyses in the field of life sciences. This includes those of the glycolytic system, TCA cycle, pentose-phosphate pathway, amino acids and nucleotides. In addition, the non ion-pairing LC-MS/MS method enables the simultaneous measurement of 97 other metabolites, such as organic acids, that cannot be measured with the ion pairing LC-MS/MS method. Details are described in Application Data Sheet Nos. 42 and 49.

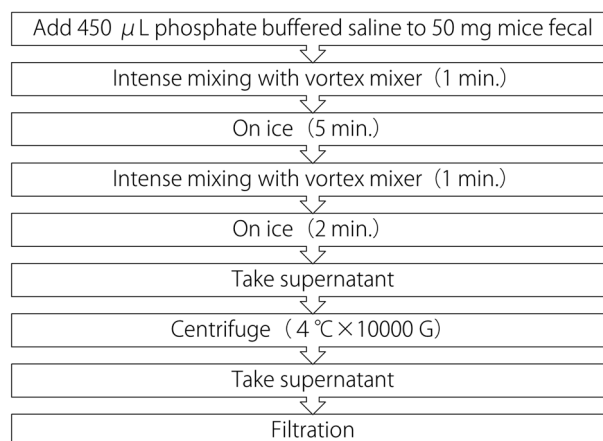


Fig. 2 Pretreatment Procedure of Mouse Feces

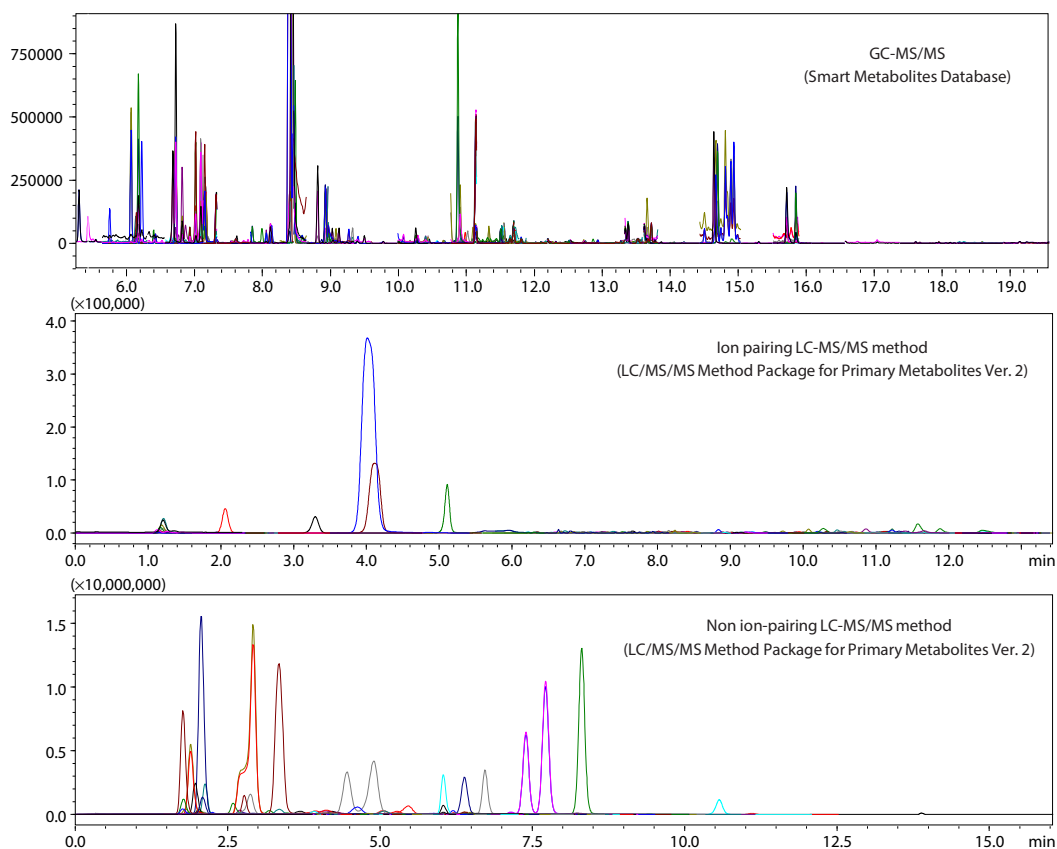


Fig. 3 MRM Chromatograms of Mouse Fecal Extracts

4. Analysis Results

The MRM chromatograms obtained from measuring mouse fecal extracts using GC-MS/MS and LC-MS/MS are shown in Fig. 3. The detected compounds and the repeatability of peak area ratios with respect to internal standard substances are listed in Table 1. GC-MS/MS analysis detected 100 compounds, mainly of short-chain fatty acids, organic acids, and sugars. Sugars are difficult to measure using LC or LC-MS/MS, but 17 compounds were detected by GC-MS/MS. The

ion pairing LC-MS/MS method detected 17 compounds, mainly which are amino acids. The non ion-pairing LC-MS/MS method detected a total of 75 components, including amino acids, nucleotides, nucleosides, and organic acids of the TCA cycle. As indicated in Fig. 4, using both GC-MS/MS and LC-MS/MS is highly effective for the comprehensive measurement of metabolites in feces.

Table 1 Metabolites Detected from Mouse Fecal Extracts and Repeatabilities

| Category | Compounds | Repeatability (%RSD, n=6) | | |
|--|--------------------------------------|---------------------------|---------------------|-------------------------|
| | | GC-MS/MS | LC-MS/MS (Ion pair) | LC-MS/MS (Non ion-pair) |
| Amino acids | Alanine | 40.4 | 3.0 | 3.0 |
| | Asparagine | - | 9.7 | 1.4 |
| | Aspartic acid | - | 1.1 | 1.7 |
| | Asymmetric dimethylarginine | - | - | 1.5 |
| | Cystine | - | - | 7.2 |
| | Glutamic acid | - | 0.5 | 1.6 |
| | Glutamine | - | 3.1 | 1.0 |
| | Glycine | 1.0 | 5.9 | 4.0 |
| | Histidine | - | 18.0 | 4.7 |
| | Isoleucine | - | - | 1.7 |
| | Leucine | 16.6 | - | 2.4 |
| | Lysine | - | 5.0 | 0.7 |
| | Methioninesulfoxide | - | - | 2.1 |
| | Phenylalanine | 10.0 | 0.6 | 1.0 |
| | Proline | - | - | 2.3 |
| | Serine | - | 9.3 | 1.9 |
| | Symmetric dimethylarginine | - | - | 4.3 |
| | Threonine | - | 3.5 | 2.4 |
| | Tryptophan | - | - | 1.2 |
| | Tyrosine | - | 2.4 | 0.9 |
| | Valine | - | - | 4.8 |
| Amino acid related metabolites | 3-Methyl-2-oxovaleric acid | 17.2 | - | - |
| | 4-Aminobenzoic acid | 8.7 | - | - |
| | 4-Aminobutyric acid | - | - | 3.9 |
| | 4-Hydroxyproline | - | - | 3.8 |
| | 5-Oxoproline | 1.4 | - | - |
| | Dimethylglycine | - | - | 5.5 |
| | N-Acetyls erine | 5.3 | - | - |
| | Pantothenic acid | 7.4 | - | 2.6 |
| Nucleic acid related metabolites | Sarcosine | 2.2 | - | - |
| | Adenine | - | - | 2.1 |
| | Adenosine | - | - | 9.6 |
| | Adenosine 3',5'-cyclic monophosphate | - | - | 14.8 |
| | Adenosine monophosphate | - | - | 3.6 |
| | Allantoin | - | - | 4.2 |
| | Cytidine | - | - | 1.2 |
| | Cytidine monophosphate | - | - | 1.3 |
| | Cytosine | - | - | 9.2 |
| | Guanine | - | - | 5.8 |
| | Guanosine | - | - | 2.1 |
| | Guanosine monophosphate | - | - | 6.4 |
| | Hypoxanthine | - | - | 1.4 |
| | Inosine | - | - | 1.5 |
| | Thymidine | - | - | 10.1 |
| | Thymidine monophosphate | - | 6.6 | 13.1 |
| | Thymine | - | - | 6.9 |
| | Uracil | - | - | 3.1 |
| | Uric acid | - | - | 7.0 |
| | Uridine | 11.9 | - | 3.7 |
| | Uridine diphosphate | - | 8.1 | - |
| Sugars | Xanthine | - | - | 7.6 |
| | Arabinose | 3.9 | - | - |
| | Fructose | 3.2 | - | - |
| | Fucose | 3.9 | - | - |
| | Galactose | 3.3 | - | - |
| | Glucose | 3.0 | - | - |
| | Lactose | 9.4 | - | - |
| | Lyxose | 3.4 | - | - |
| | Maltose | 19.2 | - | - |
| | Mannose | 5.7 | - | - |
| | Psicose | 5.0 | - | - |
| | Rhamnose | 4.2 | - | - |
| | Ribulose | 4.5 | - | - |
| | Sorbose | 6.1 | - | - |
| | Sucrose | 7.0 | - | - |
| | Tagatose | 11.1 | - | - |
| | Xylose | 6.0 | - | - |
| | Xylulose | 5.2 | - | - |
| TCA cycle | 2-Ketoglutaric acid | - | - | 2.9 |
| | Aconitic acid | - | - | 2.4 |
| | Citric acid | 4.0 | - | 6.4 |
| | Fumaric acid | 3.0 | - | - |
| | Isocitric acid | - | - | 5.0 |
| | Malic acid | - | - | 2.8 |
| | Malic acid | 3.2 | - | - |
| | Pyruvic acid | 8.8 | - | - |
| Glycolysis | Succinic acid | 1.0 | - | 2.2 |
| | Dihydroxyacetone phosphate | 7.3 | - | - |
| | Glycerol 3-phosphate | 11.4 | - | - |
| Pentose phosphate pathway | Lactic acid | 1.9 | - | 5.0 |
| | 6-Phosphogluconic acid | - | 5.6 | - |
| | Glyceric acid | 0.9 | - | - |
| Methylation and Transsulfuration cycle | Ribose | 3.5 | - | - |
| | Cystathionine | - | - | 5.1 |
| | Cysteine | - | - | 7.1 |
| Urea cycle | Methionine | - | 1.6 | 2.5 |
| | S-Adenosylhomocysteine | - | - | 8.9 |
| | Arginine | - | 5.0 | 2.6 |
| | Citrulline | - | - | 2.1 |
| | Ornithine | - | - | 2.8 |

| Category | Compounds | Repeatability (%RSD, n=6) | | |
|----------|---------------------------------|---------------------------|---------------------|-------------------------|
| | | GC-MS/MS | LC-MS/MS (Ion pair) | LC-MS/MS (Non ion-pair) |
| Others | 1,6-Anhydroglucose | 8.0 | - | - |
| | 1-Hexadecanol | 2.6 | - | - |
| | 2-Aminobutyric acid | - | - | 4.8 |
| | 2-Aminoisobutyric acid | 7.2 | - | - |
| | 2-Hydroxybutyric acid | 4.1 | - | - |
| | 2-Hydroxyglutaric acid | 1.9 | - | - |
| | 2-Hydroxyisocaproic acid | 8.1 | - | - |
| | 2-Hydroxyisovaleric acid | 7.2 | - | - |
| | 3-Hydroxybutyric acid | 2.0 | - | - |
| | 3-Hydroxyglutaric acid | 0.6 | - | - |
| | 3-Hydroxyisobutyric acid | 12.4 | - | - |
| | 3-Hydroxyisovaleric acid | 4.2 | - | - |
| | 3-Hydroxyphenylacetic acid | 4.7 | - | - |
| | 3-Hydroxypropionic acid | 1.6 | - | - |
| | 3-Methoxy-4-hydroxybenzoic acid | 6.7 | - | - |
| | 3-Phenylacetic acid | 4.4 | - | - |
| | 4-Hydroxybenzoic acid | 4.9 | - | - |
| | 4-Hydroxyphenylacetic acid | 5.8 | - | - |
| | 4-Hydroxyphenylacetic acid | 7.6 | - | - |
| | Acetoacetic acid | 4.0 | - | - |
| | Acetylcamitine | - | - | 3.0 |
| | Adipic acid | 7.2 | - | - |
| | Arabitol | 5.7 | - | - |
| | Argininosuccinic acid | - | - | 14.1 |
| | Azelic acid | 5.0 | - | - |
| | Benzoic acid | 17.4 | - | - |
| | Caproic acid | 2.8 | - | - |
| | Carnitine | - | - | 6.9 |
| | Carnosine | - | - | 2.0 |
| | Catechol | 13.1 | - | - |
| | Cholesterol | 14.7 | - | - |
| | Cholic acid | - | - | 2.6 |
| | Choline | - | - | 1.6 |
| | Creatine | - | - | 2.5 |
| | Creatinine | - | - | 3.0 |
| | Decanoic acid | 8.5 | - | - |
| | Dopamine | - | - | 6.4 |
| | Elaidic acid | 8.8 | - | - |
| | FMN | - | - | 8.0 |
| | Galacturonic acid | 5.6 | - | - |
| | Glucuronic acid | 7.1 | - | - |
| | Glutaric acid | 2.7 | - | - |
| | Glycolic acid | 2.5 | - | - |
| | Glyoxylic acid | 5.7 | - | - |
| | Histamine | - | - | 2.6 |
| | Hydroquinone | 3.7 | - | - |
| | Inositol | 1.1 | - | - |
| | Lauric acid | 16.7 | - | - |
| | Malonic acid | 3.8 | - | - |
| | Mannitol | 3.8 | - | - |
| | Margaric acid | 6.0 | - | - |
| | meso-Erythritol | 2.6 | - | - |
| | Methylsuccinic acid | 5.7 | - | - |
| | Monostearin | 38.8 | - | - |
| | Myristic acid | 3.9 | - | - |
| | N-Acetylmannosamine | 8.9 | - | - |
| | Nicotinic acid | 2.2 | - | 3.4 |
| | Nonanoic acid | 5.6 | - | - |
| | Norepinephrine | - | - | 1.7 |
| | Octadecanol | 4.1 | - | - |
| | Octanoic acid | 4.2 | - | - |
| | Oleic acid | 7.4 | - | - |
| | Palmitic acid | 1.4 | - | - |
| | Phenylacetic acid | 1.4 | - | - |
| | Phosphoric acid | 1.1 | - | - |
| | Protocatechuic acid | 9.9 | - | - |
| | Pyridoxine | 10.6 | - | - |
| | Ribitol | 11.5 | - | - |
| | Sebacic acid | 7.0 | - | - |
| | Serotonin | - | - | 2.4 |
| | Stearic acid | 1.4 | - | - |
| | Suberic acid | 4.2 | - | - |
| | Taurocholic acid | - | - | 4.1 |
| | Threitol | 3.3 | - | - |
| | Threonic acid | 8.1 | - | - |
| | Triethanolamine | 5.2 | - | - |
| | Xylitol | 7.0 | - | - |

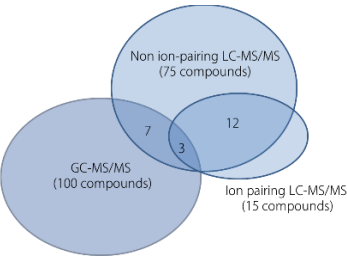


Fig. 4 The Number of Metabolites Detected from Mouse Fecal Extracts

5. A Comparison of Young Mice and Old Mice

As an example application, we analyzed and compared the feces of male C57BL/6J mice reared in a normal environment but of different age in weeks (10 weeks old and 70 weeks old). A total of 66 metabolites were detected. In order to compare the content of metabolites in both mice, we used the Shimadzu Multi-omics Data Analysis Pack to create graphs of peak area ratios of the 66 detected metabolites with respect to internal standard substances. The results as indicated in Fig. 5 suggest that there is a difference between the two groups with regard to various metabolites including amino acids such as arginine, aspartate, and lysine, nucleic acid metabolites such as adenine and adenosine, and cholic acid. In terms of cholic acid, results indicated that the content tends to be lower in the feces of old mice compared to young mice. Cholic acid results from the deconjugation of

taurocholic acid. Probable causes of the low cholic acid content in the feces of old mice may be that the secretion of bile acids or the number of bacteria which cause deconjugation decreased due to age. Regarding arginine, the content tended to be lower in the feces of young mice compared to old mice. Arginine in the intestinal lumen is absorbed and used for the synthesis of polyamines which are indispensable for cell growth. This suggests that the arginine content may be low in young mice because they are actively absorbing arginine from the intestinal lumen and synthesizing polyamines. As demonstrated so far, this method can also be considered to be effective in investigating the factors of aging using feces.



Fig. 5 A Comparison of the Individual Metabolite Content Detected in the Fecal Extracts of Young Mice and Old Mice

6. Conclusion

We introduced a comprehensive measurement of the metabolites in mouse feces using GC-MS/MS and LC-MS/MS. Since the compounds that can be measured differ between GC-MS/MS and LC-MS/MS, using both analysis methods enables a comprehensive measurement of the metabolites. In addition, by using the MRM database for GC-MS/MS and the method package for LC-MS/MS, comprehensive

measurement can be done easily even by those who are not very familiar with measurement. The immense data obtained by the measurements can be then visualized using the Shimadzu Multi-omics Data Analysis Pack, which will facilitate the examination of the results. These measurement and analysis methods are expected to be highly effective for metabolomic analyses.

References

- (1) M. Matsumoto, et al., Scientific Reports 2, 223 (2012).

Smart Metabolites Database and GCMS-TQ are trademarks of Shimadzu Corporation.



Shimadzu Corporation

www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. Shimadzu disclaims any proprietary interest in trademarks and trade names used in this publication other than its own. See <http://www.shimadzu.com/about/trademarks/index.html> for details.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.