

# Application News

## No. G290A

### Gas Chromatograph

## Analysis of Residual Solvents in drug products using Nexis GC-2030 combined with HS-20 head space sampler - USP <467> Residual Solvents Procedure A -

Residual solvents in pharmaceuticals are defined as volatile organic compounds used in or generated from the manufacture of drug substances, pharmaceutical additives, or drug products. They are strictly controlled according to risk classifications from Class 1 to Class 3, which are based on the risk to human health.

Headspace GC methods specified in the USP (U.S. Pharmacopeia), General Chapters <467> Residual Solvents, are commonly used for analysis of residual solvents. This Application News presents data obtained using the Shimadzu HS-20 Headspace Sampler and Nexis GC-2030 Gas Chromatograph, from Class 1 and Class 2 standard solutions, in accordance with Water-Soluble Articles, Procedure A, in USP <467> Residual Solvents.

E. Kobayashi, T. Murata



Fig. 1 Nexis GC-2030 + HS-20

### Class 1

Fig.2 shows the Class 1 standard solution chromatogram. Procedure A requires that the S/N ratio obtained for 1,1,1-Trichloroethane in this chromatogram be 5 or higher. As shown, the S/N ratio was 220. Even for carbon tetrachloride, which had the lowest sensitivity level, the S/N was 20.

Table 3 indicates the S/N ratio of each peaks and the repeatability of the peak area (n=6).

Table 3 S/N ratio and Repeatability of Class1

No.	Compounds	S/N ratio	%RSD (n=6)
1	1,1-Dichloroethane	320	2.8
2	1,1,1-Trichloroethane	220	2.3
3	Carbon tetrachloride	20	2.9
4	Benzene	170	2.5
5	1,2-Dichloroethane	60	3.4

### Instruments and Analytical Conditions

Table 1 GC Method for USP 467 Procedure A

Model	: Nexis GC-2030
Detector	: FID-2030
Headspace Sampler	: HS-20
Column	: SH-Rxi-624 Sil MS (0.32 mm I.D. x 30 m, d.f. = 1.8 µm)
Column Temperature	: 40 °C (20 min) - 10 °C /min - 240 °C (20 min) Total 60 min
Injection Mode	: Split 1 : 5
Carrier Gas Controller	: Constant Linear Velocity (He)
Linear Velocity	: 35 cm/sec
Detector Temperature	: 250 °C
Injection Volume	: 1 mL

Table 2 HS-20 Method for USP 467 Procedure A

Oven Temperature	: 80 °C
Sample Line Temperature	: 110 °C
Transfer Line Temperature	: 120 °C
Vial Stirring	: Off
Vial Volume	: 20 mL
Vial Heat-retention Time	: 60 min
Vial Pressurization Time	: 1 min
Vial Pressure	: 75 kPa
Loading Time	: 1 min
Needle Flush Time	: 5 min

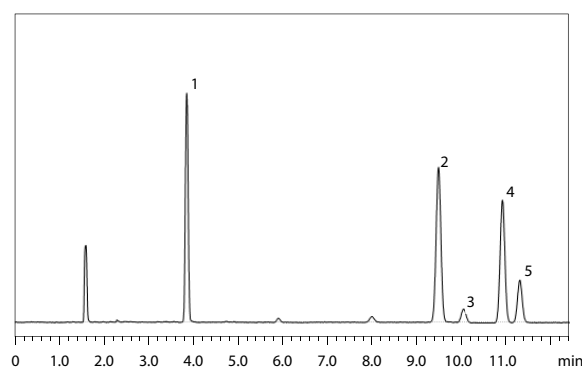


Fig. 2 Chromatogram of WATER-SOLUBLE ARTICLES Class1 Standard Solution by Procedure A

■ Class2

Due to the large number of components in the Class 2 standard solution, it was separated into two mixtures: A and B. Respective measurement results are shown in Fig.3 and Fig.4. Procedure A requires that the resolution for acetonitrile and methylene chloride in

the Class 2 standard solution Mixture A chromatogram be 1.0 or greater.

Fig.3 shows that, using the Restek SH-Rxi-624SiIMS low-bleed column, the specified peaks are completely separated, with a resolution of 2.4.

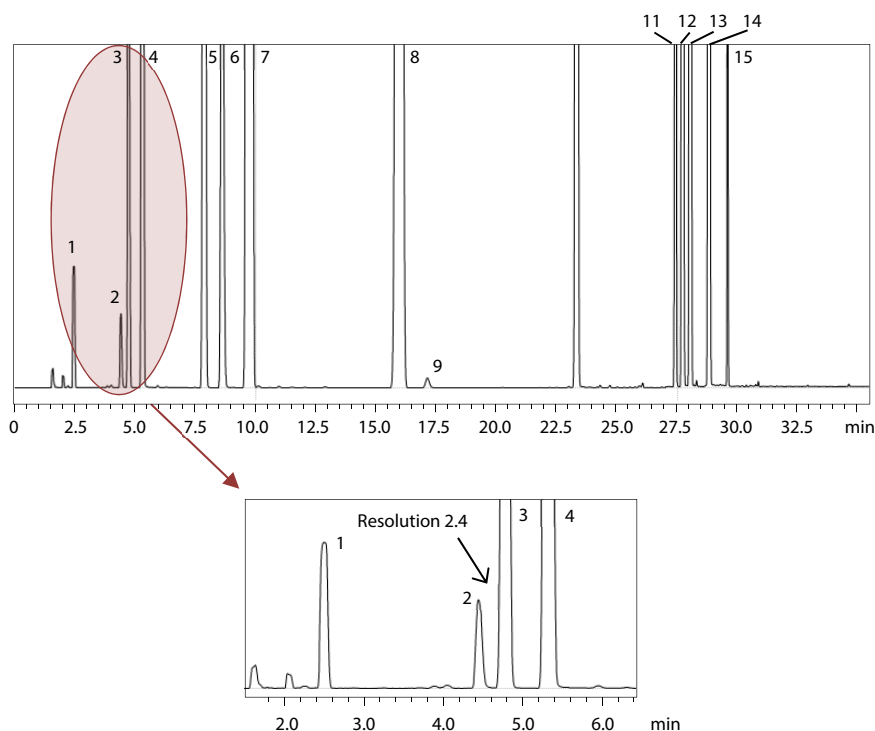


Fig. 3 Chromatogram of WATER-SOLUBLE ARTICLES Class 2A Standard Solution by Procedure A

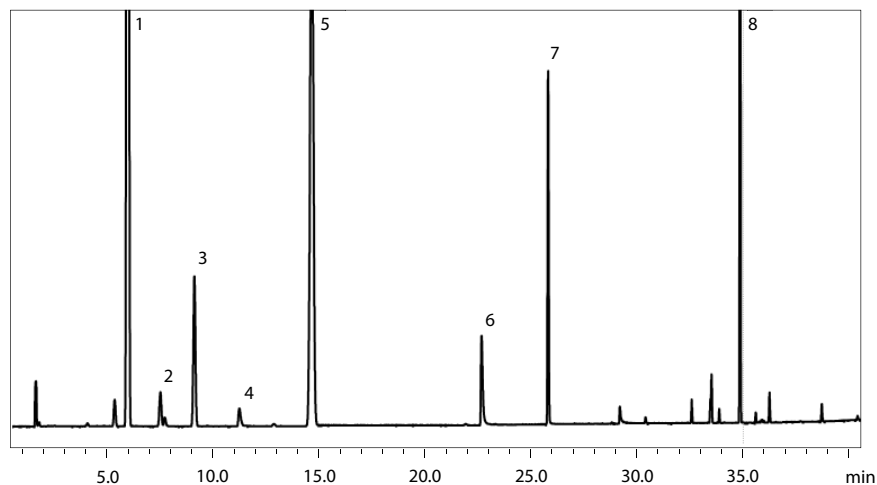


Fig. 4 Chromatogram of WATER-SOLUBLE ARTICLES Class 2B Standard Solution by Procedure A



Shimadzu Corporation  
www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedure.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. Company names, product/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation or its affiliates, whether or not they are used with trademark symbol "TM" or "®". Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services. Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

First Edition: Jun. 2017  
Second Edition: Mar. 2021