

Application Data Sheet

No. 61

System Gas Chromatograph

NGA analysis with ISO6974-4 Nexis GC-2030ISO6974-4 GC-2014ISO6974-4

This method is for determining the chemical composition of natural gases and similar gaseous mixtures within the composition range shown in the specification sheet. This test method provides data for calculating physical properties of the sample, such as heating value and relative density, or for monitoring the concentrations of one or more of the components in a mixture. A total of 1 valves and 2 columns are applied in this GC system. Sample is introduced into the sample loop for determination. Using a pre-column, C6+ components are back-flushed as a single peak. The valve timing then allows C1-C5, N₂, CO₂ to be eluted to TCD through DC-200 in that order. The final analysis time is approximately 22 minutes. The system includes LabSolution workstation software and BTU and Specific Gravity calculation software.

Analyzer Information

System Configuration:

One valve / two packed columns with one TCD detector

Sample Information:

Permanent gas, C1-C6

Concentration Range:

No.	Name of Compound	Concentration Range		Detector
		Low Conc.	High Conc.	
1	N ₂	0.001%	15.0%	TCD-1
2	CH ₄	75.0%	100.0%	TCD-1
3	CO ₂	0.001%	10.0%	TCD-1
4	C ₂ H ₆	0.001%	10.0%	TCD-1
5	C ₃ H ₈	0.001%	3.0%	TCD-1
6	i-C ₄ H ₁₀	0.001%	1.0%	TCD-1
7	n-C ₄ H ₁₀	0.001%	1.0%	TCD-1
8	Neo-C ₅ H ₁₂	0.001%	0.5%	TCD-1
9	i-C ₅ H ₁₂	0.001%	0.5%	TCD-1
10	n-C ₅ H ₁₂	0.001%	0.5%	TCD-1
11	C ₆ +	0.001%	0.2%	TCD-1

Detection limits may vary depending on the sample. Please contact us for more consultation.

System Features

- Versatile software easy GC system operation
- One TCD channel
- Good repeatability

Typical Chromatograms

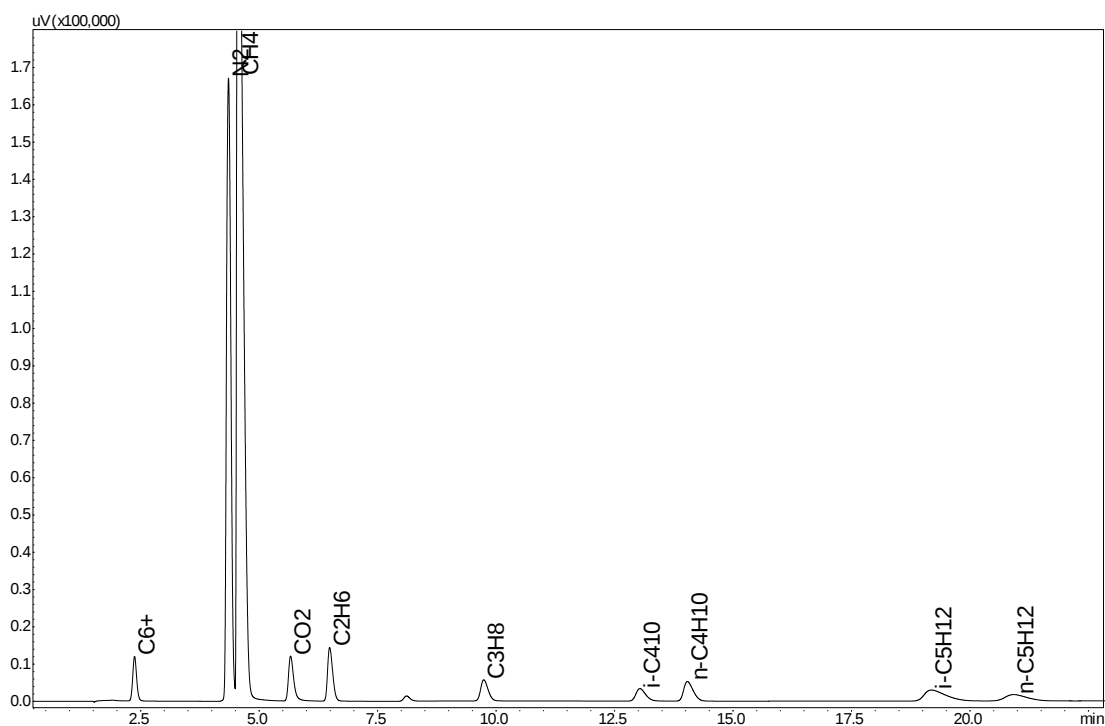


Fig. 2 Chromatogram of TCD

First Edition: November, 2017



Shimadzu Corporation
www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.
The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

© Shimadzu Corporation, 2017