

CoreFocus
Report

No. 122

LC


Reversed-phase

Shim-pack Series

Shim-pack™ GIST C18

High-Speed Analysis of Pramipexole


Keywords: Parkinson's disease, impurity

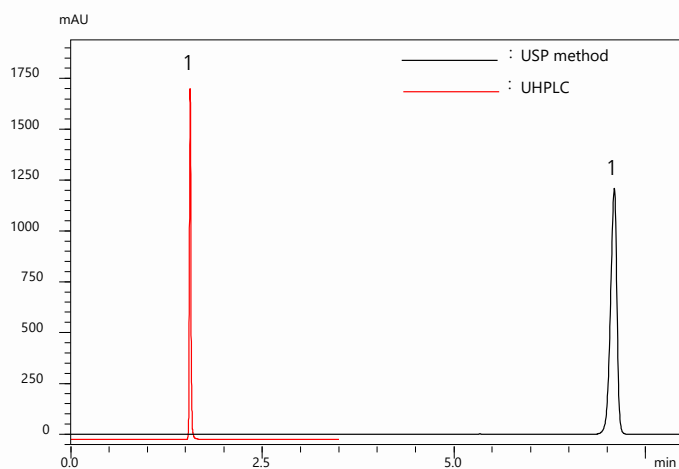

1. Pramipexole


Chromatogram of Pramipexole Dihydrochloride

USP method

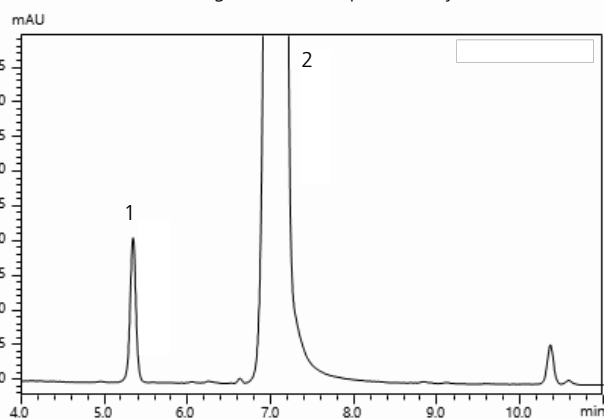

1. Compound A
2. Pramipexole

UHPLC method

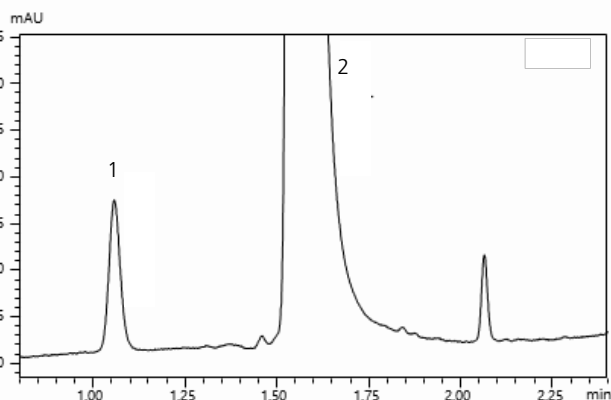

1. Compound A
2. Pramipexole

Enlarged Chromatograms of Vicinity of Pramipexole Dihydrochloride

Analytical Conditions Conforming to USP

System	: Nexera™-i MT
Column	: Shim-pack VP-ODS (150 mm L. × 4.6 mm I.D., 5.0 µm) , P/N : 228-34937-91
Mobile phase	: A) 67 mmol/L (Potassium) Phosphate buffer (pH 3.0) Containing 21 mmol/L 1-Octanesulfonic acid sodium salt B) Solution A / Acetonitrile (1/1)
Flow rate	: 1.5 mL/min
Time program	: B Conc. 40 % (0 min) -> 80 % (15 min) -> 40 % (15.1 - 20 min)
Column temp.	: 40 °C
Injection volume	: 5 µL
Detection	: UV 264 nm
Sample	: Pramipexole Dihydrochloride

Analytical Conditions for UHPLC

System	: Nexera™-i MT
Column	: Shim-pack GIST C18 (50 mm L. × 2.1 mm I.D., 2.0 µm), P/N : 227-30001-02
Mobile phase	: A) 67 mmol/L (Potassium) Phosphate buffer (pH 3.0) Containing 21 mmol/L 1-Octanesulfonic acid sodium salt B) Solution A / Acetonitrile (1/1)
Flow rate	: 0.78 mL/min
Time program	: B Conc. 40 % (0 min) -> 80 % (2.00 min) -> 40 % (2.01 - 3.5 min)
Column temp.	: 40 °C
Injection volume	: 1 µL
Detection	: UV 264 nm
Sample	: Pramipexole Dihydrochloride

Source : Application News L524 ([JP](#), [ENG](#))

CoreFocus, Shim-pack and Nexera are trademarks of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.

Shimadzu Corporation

www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu.

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®".

Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

Related Products

Some products may be updated to newer models.

HPLC Column Solutions – Reversed phase and Normal phase –

Shim-pack

> Shim-pack GIST Series

HPLC Column Solutions

Related Solutions

> Small Molecule Pharmaceutical

> 低分子医薬品

> Price Inquiry

> Product Inquiry

> Technical Service / Support Inquiry

> Other Inquiry