


CoreFocus
Report
No.135

LC SEC (GPC) Shim-pack Series

Shim-pack™ GPC 805+GPC 801

Analysis of Polymer Additives

Keywords: antioxidant, light stabilizer, tinuvin, irganox

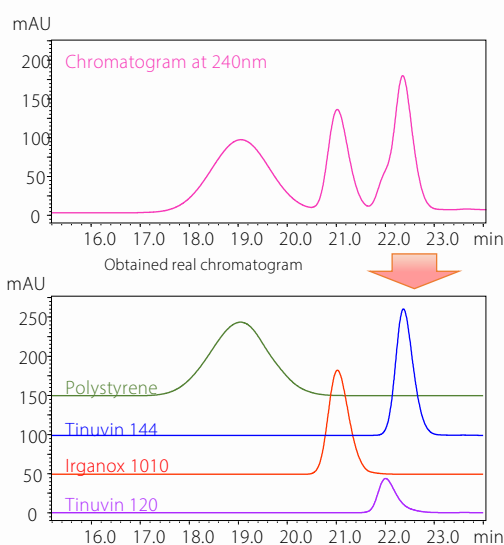

Deconvolution Result of Three Polymer Additives with *i*-PDeA II (PDA)

System	: Nexera™ GPC
Column	: Shim-pack GPC 805+GPC 801 (300 mm L x 8 mm I.D. for each), P/N : 228-20808-91(GPC 805), P/N : 228-20803-91(GPC 801)
Mobile phase	: THF
Flow rate	: 0.8 mL/min
Column temp.	: 40 °C
Injection vol.	: 10 µL
Detection	: RID / PDA (220-400 nm)
Sample	: 0.5 % Polystyrene containing three additives
Cycle time	: 31 min
Overlap time	: 15.5 min

Source : Application News L.537 ([JP](#), [ENG](#))

CoreFocus, Shim-pack and Nexera are trademarks of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.

Shimadzu Corporation
www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu.

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®".

Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

Related Products

Some products may be updated to newer models.

HPLC Column Solutions - Reversed phase and Normal phase -

Shim-pack

> Shim-pack GPC Series

HPLC Column

Related Solutions

> Plastics

> プラスチック-化学

> Price Inquiry

> Product Inquiry

> Technical Service / Support Inquiry

> Other Inquiry