

CoreFocus
Report
No.285


LC

Ion-exclusion

Shim-pack Series

Shim-pack™ Fast-OA and Shim-pack SCR-102H
Analysis of Organic acids by the post-column pH
buffering-electric conductivity detection method

Keywords: biomass fermentation sample, biofuel

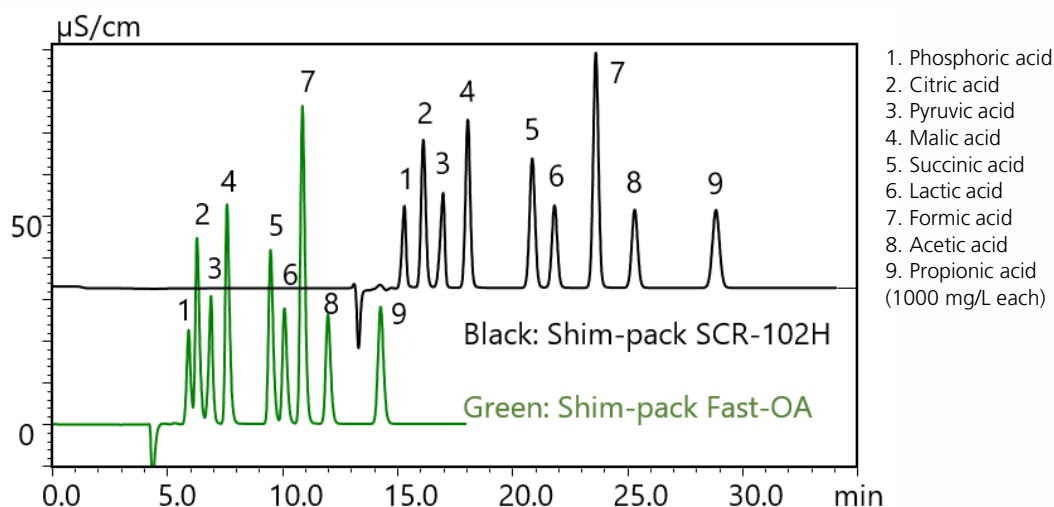

System	: Nexera™ organic acid analysis system
Column	: Shim-pack SCR-102H (300 mm x 8.0 mm I.D., 7 μm) x2, P/N : 228-17893-91
(Conventional)	: Guard column SCR-102H (50 mm x 6.0 mm I.D.), P/N : 228-17924-91
Column	: Shim-pack Fast-OA (100 mm x 7.8 mm I.D., 5 μm) x3, P/N : 228-59942-41
(Fast)	: Shim-pack Fast-OA(G) (10 mm x 4.0 mm I.D.), P/N : 228-59942-42
Mobile phase	: Reagents kit for Organic Acid Analysis System, P/N : 228-61465-91
Flow rate	: 1.0 mL/min
pH Buffering solution	: Reagents kit for Organic Acid Analysis System, P/N : 228-61465-91
Mixer	: MR 20
Column temp.	: 40 °C
Injection vol.	: 10 μL
Vial	: Shimadzu Vials, LC, 1.5 mL, Glass, P/N : S228-15652-92
Detection	: Conductivity detector

Source : Application News L588 ([JP](#), [ENG](#))

CoreFocus, Nexera and Shim-pack are trademarks of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.

Shimadzu Corporation
www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu.

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®". Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

➤ Please fill out the survey

Related Products

Some products may be updated to newer models.

HPLC Column Solutions – Reversed phase and Normal phase –
Shim-pack

➤ Shim-pack Fast-OA
HPLC Column

HPLC Column Solutions – Reversed phase and Normal phase –
Shim-pack

➤ Shim-pack SCR
Series
HPLC Column

Related Solutions

➤ Biomass-SCJ

➤ バイオマス

➤ Price Inquiry

➤ Product Inquiry

➤ Technical Service /
Support Inquiry

➤ Other Inquiry