

**CoreFocus**  
**Report**  
**No.73**

LC

Reversed-phase


Shim-pack Series

**Shim-pack™ GIST C18-AQ**

**Analysis of Caffeic acid in *Dandelion***

**Keywords: Carcinogenic**

Chromatogram of Standard Mix

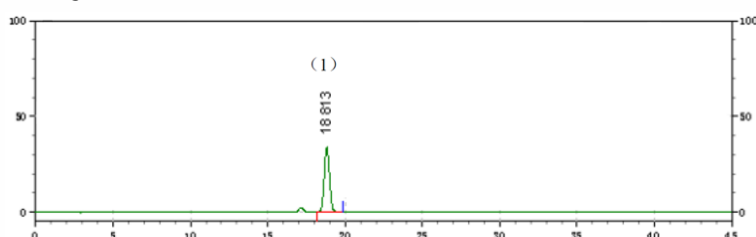


1. Caffeic acid

Standard Mix

Caffeic acid (10 µg/mL in 70% Methanol)

Chromatogram of *Dandelion*

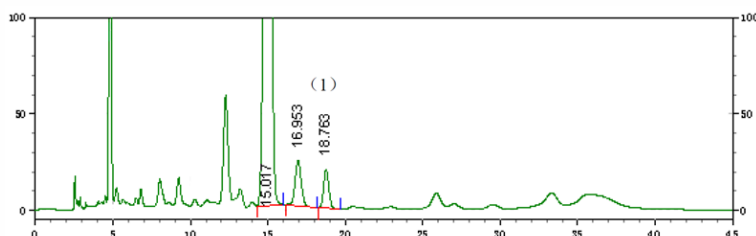

Pretreatment procedure

- Weigh 1 g of sample (commercially available *Dandelion* powder) and transfer into a 50 mL stoppered conical flask.
- Add 10 mL of 5% Formic acid in Methanol solution into flask.
- Close the stopper and reweigh.
- Ultrasonically treat it (power 250 W, frequency 40 kHz) for 30 minutes.
- After cooling, weigh the conical flask.
- Add 5% Formic acid in Methanol to bring the final weight to the pre-reflux weight (Step c).
- Shake well, Centrifuge and collect supernatant.

Column	: Shim-pack GIST C18-AQ (250 mm x 4.6 mm I.D., 5 µm )	P/N: 227-30742-08
Mobile phase	: Phosphate Buffer solution (PH 3.9) : Methanol = 78.5 : 21.5	
Flow rate	: 1.0 mL/min	
Column temp.	: 40°C	
Injection vol.	: 10 µL	
Vial	: LabTotal™ Vial for LC/LCMS, P/N : 227-34001-01	
Detection	: UV at 323 nm	

CoreFocus, Shim-pack and LabTotal are trademarks of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.

**Shimadzu Corporation**  
**www.shimadzu.com/an/**

**For Research Use Only. Not for use in diagnostic procedures.**

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu.

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®". Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".


Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

➤ Please fill out the survey

## Related Products

Some products may be updated to newer models.

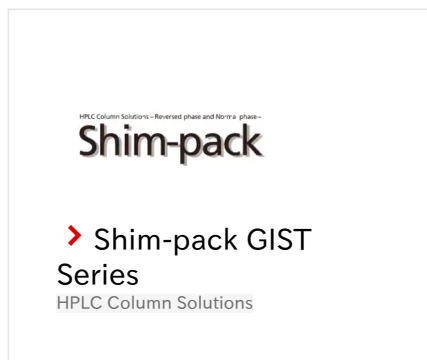

## Related Solutions

➤ Small Molecule  
Pharmaceutical

➤ 低分子医薬品

➤ Price Inquiry

➤ Product Inquiry

➤ Technical Service /  
Support Inquiry

➤ Other Inquiry