

CoreFocus
Report
No.406


GCMS

SH Series

SH-I-5Sil MS

Determine the Phthalate Content in E-liquid

Keywords: the National Food Safety Standards, e-cigarettes

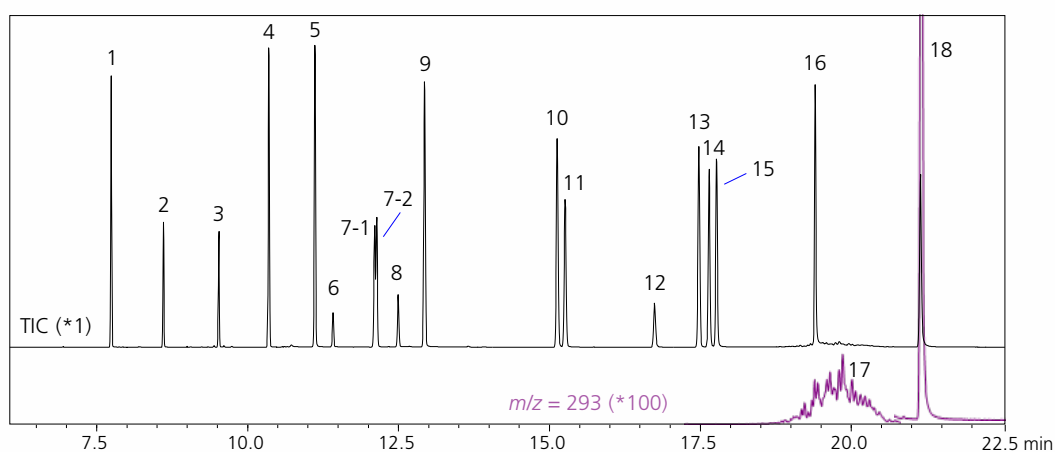


- | | |
|------------------------------------|----------------------------------|
| 1. Dimethyl phthalate | 10. Di-n-hexyl phthalate |
| 2. Diethyl phthalate | 11. Butyl benzyl phthalate |
| 3. Diallyl phthalate | 12. Bis(2-butoxyethyl) phthalate |
| 4. Diisobutyl phthalate | 13. Dicyclohexyl phthalate |
| 5. Dibutyl phthalate | 14. Bis(2-ethylhexyl) phthalate |
| 6. Bis(methylglycol) phthalate | 15. Diphenyl phthalate |
| 7-1. Bis(4-methylpentyl) phthalate | 16. Dioctyl phthalate |
| 7-2. Bis(4-methylpentyl) phthalate | 17. Diisononyl phthalate |
| 8. Bis(2-ethoxyethyl) phthalate | 18. Dinonyl phthalate |
| 9. Di-n-pentyl phthalate | |

TIC Chromatogram for the Phthalate Standards and the m/z 293 Ion Peak for Diisononyl phthalate (No. 17)

Model	: GCMS-QP™2010 NX
Column	: SH-I-5Sil MS (30 m x 0.25 mm I.D., 0.25 μm) P/N : 221-75954-30
Column temp.	: 60 °C (1 min) - 20 °C/min - 220 °C (1 min) - 5 °C/min - 250 °C (1 in) - 20 /min - 290 °C (7.5 min)
Injection temp.	: 260 °C
Carrier gas control mode	: Constant flow (81 mL/min)
Injection mode	: Splitless
Injection volume	: 1.0 μL
Ionization mode	: EI
Ion source temp.	: 230 °C
Interface temp.	: 280 °C
Data acquisition mode	: SIM

Compounds, Retention Times, and Selected Ions

No.	Compound	Abbreviation	Retention Time (min)	CAS No.	Quantitative Ion (m/z)	Qualitative Ion (m/z)
1	Dimethyl phthalate	DMP	7.740	131-11-3	163	77, 194
2	Diethyl phthalate	DEP	8.605	84-66-2	149	177, 105
3	Diallyl phthalate	DAP	9.522	131-17-9	149	189, 132
4	Diisobutyl phthalate	DIBP	10.351	84-69-5	149	223, 163
5	Dibutyl phthalate	DBP	11.116	84-74-2	149	223, 205
6	Bis(methylglycol) phthalate	DMEP	11.415	117-82-8	149	104, 176
7-1	Bis(4-methylpentyl) phthalate	BMPP-1	12.102	146-50-9	149	85, 167
7-2	Bis(4-methylpentyl) phthalate	BMPP-2	12.138	146-50-9	149	85, 167
8	Bis(2-ethoxyethyl) phthalate	DEEP	12.491	605-54-9	149	104, 193
9	Di-n-pentyl phthalate	DPP	12.930	131-18-0	149	237, 219
10	Di-n-hexyl phthalate	DHXP	15.126	84-75-3	149	251, 233
11	Butyl benzyl phthalate	BBP	15.258	85-68-7	149	91, 206
12	Bis(2-butoxyethyl) phthalate	DBEP	16.739	117-83-9	149	101, 193
13	Dicyclohexyl phthalate	DCHP	17.473	84-61-7	149	167, 249
14	Bis(2-ethylhexyl) phthalate	DEHP	17.641	117-81-7	149	167, 279
15	Diphenyl phthalate	DPhP	17.764	84-62-8	225	77, 104
16	Diocetyl phthalate	DNOP	19.402	117-84-0	279	261
17	Diisononyl phthalate	DINP	19.802	28553-12-0	293	127
18	Dinonyl phthalate	DNP	21.141	84-76-4	149	293, 275

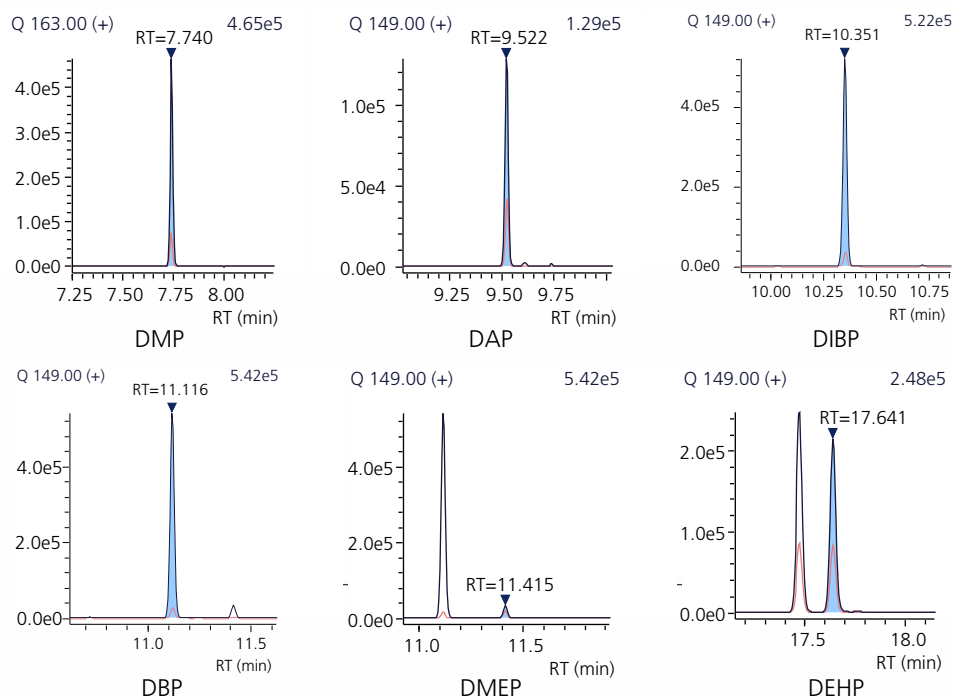

Mass Chromatograms for Some of the Compounds

Source : Application News 03-GCMS-407 ([JP](#), [ENG](#))

CoreFocus and GCMS-QP are trademarks of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.

Shimadzu Corporation

www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu.

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®". Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

➤ Please fill out the survey

Related Products

Some products may be updated to newer models.

HPLC Column Solutions ~ Reversed phase and Normal phase ~
Shim-pack

➤ SH-I-5SiI MS

Related Solutions

➤ Air

➤ 大気

➤ Price Inquiry

➤ Product Inquiry

➤ Technical Service /
Support Inquiry

➤ Other Inquiry