

CoreFocus
Report
No.407


GCMS

AOC

SH Series

SH-PolarWax
Determination of Fatty Acids Methyl
Esters (FAMES) in Aviation Turbine Fuel
as per IP585

Keywords: petroleum product, fuel

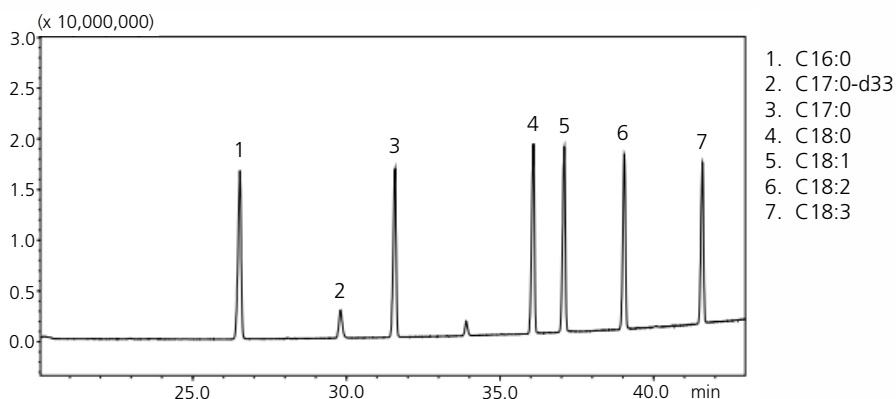

System Configuration

Model : GCMS-QP™2020 NX/AOC™-20i plus + AOC-20s plus
Glass insert : Splitless insert with deactivated glass wool
P/N : 227-35008-01
Column : SH-PolarWax (60m x 0.25 mm I.D., 0.5 µm)
P/N : 227-36248-02

GC

Injection temp. : 260 °C
Flow mode : Constant flow
Column flow : 0.8 mL/min
Purge flow : 3.0 mL/min
Column oven temp. : 150 °C (5 min) – 12 °C/min – 200 °C (17 min) –
3 °C/min – 252 °C (6.5 min)
Injection volume : 1.0 µL

MS

Ion source temp. : 230 °C
Interface temp. : 260 °C
Solvent cut time : 20.00 min
Detector voltage : 1.27 kV
Event time : Scan 0.1 s / SIM 0.3 s
Scan range (m/z) : 32 – 332

Mass table

Analyte	Type	Ions m/z
C16:0	Target	227, 239, 270, 271
C17:0 d33	ISTD	317
C17:0	Target	241, 253, 284
C18:0	Target	255, 267, 298
C18:1	Target	264, 265, 296
C18:2	Target	262, 263, 264, 294, 295
C18:3	Target	236, 263, 292, 293

Source : Application News 07-GC-21007 ([JP](#), [ENG](#))

CoreFocus, GCMS-QP and AOC are trademarks of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.

Shimadzu Corporation
www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu.

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®". Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

➤ Please fill out the survey

Related Products

Some products may be updated to newer models.

HPLC Column Solutions - Reversed phase and Normal phase -
Shim-pack

➤ SH-PolarWax

Related Solutions

Hydrocarbon
➤ Processing Industry
(Petrochemical, Ch

➤ 石油・石油系化学物
質

➤ Price Inquiry

➤ Product Inquiry

➤ Technical Service /
Support Inquiry

➤ Other Inquiry