

CoreFocus Report

No.580

LC-MS


Reversed-phase

Shim-pack™ Series

Shim-pack GISS-HP C18

Analysis of Fumonisinin Livestock Feed Blend

Keywords: Mycotoxins, Grain, Food safety

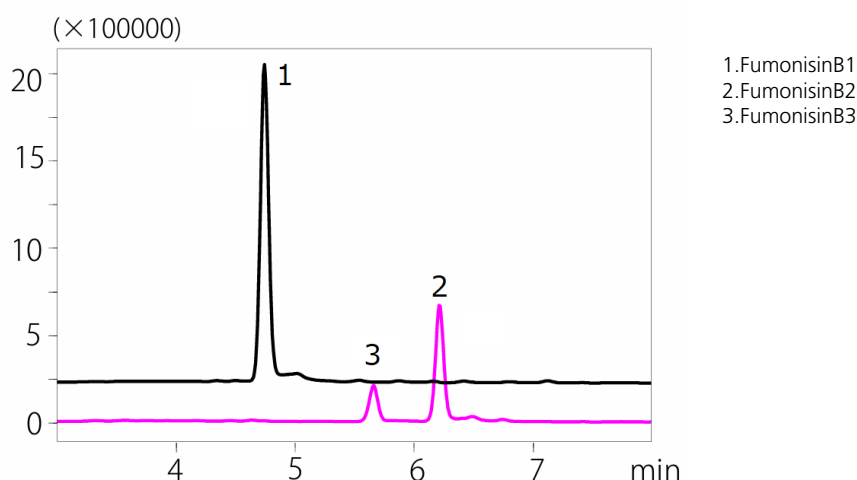

[HPLC conditions] (Nexera™ X3)

Column : Shim-pack GISS-HP C18,
150 mm x 2.1 mm I.D., 3 μm*1
Mobile phase A : 0.1 % formic acid aq.
Mobile phase B : Acetonitrile
Time program : 30 % B(0 min) → 50 % B(5-8 min) →
95 % B(8.01-10 min) → 30 % B(10.01 min)
Flow rate : 0.2 mL/min
Column oven : 40 °C
Injection vol. : 3 μL
Needle rinse
conditions : R0: Acetonitrile/water = 1:1
R1: 10 mmol/L sodium citrate aq.
R2: 1 % formic acid aq./Methanol/
Acetonitrile/isopropyl alcohol = 1:1:1:1
Rinse type: needle inner and outer surface rinse
Needle rinse program:
Needle inner surface rinse: R1 – R2 – R0
Needle outer surface rinse: R2

1*: P/N : 227-30084-03

[MS conditions] (LCMS-2050)

Ionization : ESI/APCI(DUIS™), Positive mode
Mode : SIM
(FB1: *m/z* 722.40, FB2/FB3: *m/z* 706.40)
Nebulizing gas flow : 1.5 L/min
Drying gas flow : 3.0 L/min
Heating gas flow : 5.0 L/min
DL temp. : 200 °C
Desolvation temp. : 500 °C
Interface voltage : +1.0 kV

Source : Application News 01-00649 ([JP,ENG](#))

CoreFocus, Shim-pack, Nexera and DUIS are trademarks of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.

Shimadzu Corporation

www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu.

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®".

Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".


Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

➤ Please fill out the survey

Related Products

Some products may be updated to newer models.

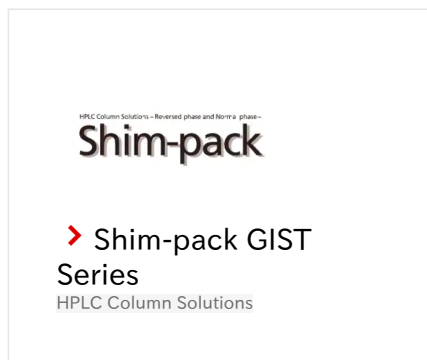

Related Solutions

➤ Natural Toxins

➤ 自然毒

➤ Price Inquiry

➤ Product Inquiry

➤ Technical Service /
Support Inquiry

➤ Other Inquiry