


CoreFocus Report

No.590

LC-MS Reversed-phase Shim-pack™ Series

Shim-pack GIST-HP C18 Analysis of PFAS

Keywords: ASTM

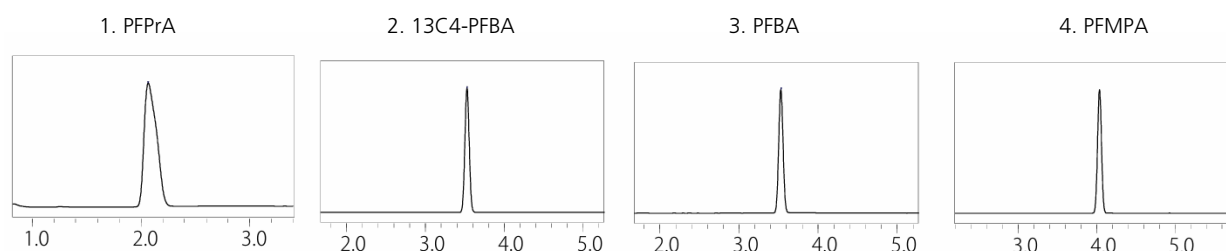

[HPLC conditions] (Nexera™ X3)

Analytical column	: Shim-pack GIST-HP C18, 100 mm x 3.0 mm I.D., 3 μm ^{*1}
Delay column	: Shim-pack Scepter C18-120, 100 mm x 2.1 mm I.D., 3 μm ^{*2}
Mobile phase A	: 2 mmol/L Ammonium acetate in H ₂ O/Acetonitrile = 95/5
Mobile phase B	: Acetonitrile
Gradient program	: B conc. 10 %(0 min) → 22 %(2.3-3.0 min) → 45 %(6.0 min) → 80 %(13.0 min) → 95 %(14.0-16 min) → 10 %(16.01-20.0 min)
Flow rate	: 0.6 mL/min
Interface	: IonFocus™ ESI
Flow rate	: 0.6 mL/min
Column oven temp.	: 40 °C
Injection vol.	: 25 μL
Multiple draw injection program	: Co-injection 25 μL sample → 25 μL 0.1 % acetic acid in H ₂ O

[MS conditions] (LCMS-8060NX)

Ionization	: ESI (Negative mode)
Interface temp.	: 170 °C
Interface voltage	: -0.5 kV
Focus voltage	: -2 kV
ESI Probe position	: +3 mm
Nebulizing gas flow	: 3 L/min
Drying gas flow	: 5 L/min
Heating gas flow	: 15 L/min
DL temp.	: 200 °C
Heat block temp.	: 300 °C

1*: P/N : 227-30040-04, 2*: P/N : 227-31014-05

Source : Application News 01-00150 ([JP](#), [ENG](#))

CoreFocus, Shim-pack, Shim-pack Scepter, Nexera and IonFocus are trademarks of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.

Shimadzu Corporation

www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu.

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®".

Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".


Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

➤ Please fill out the survey

Related Products

Some products may be updated to newer models.

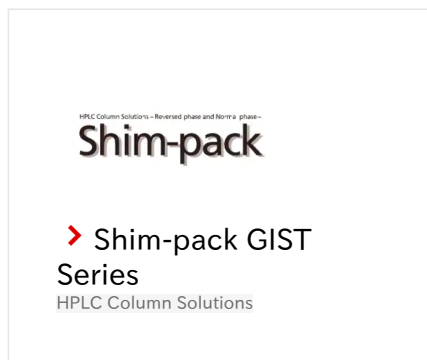

Related Solutions

➤ PFAS

➤ 有機フッ素化合物
(PFAS)

➤ Price Inquiry

➤ Product Inquiry

➤ Technical Service /
Support Inquiry

➤ Other Inquiry