

Application News

Gas Chromatograph-Mass Spectrometry

Ultra-Sensitive Analysis of \(\beta\)-cyclocitral in Water Using Solid Phase Micro Extraction and Gas Chromatography-Mass Spectrometry

Hoonju Kim, Jin Jang Shimadzu Scientific Korea

User Benefits

- ◆ SPME-GC-MS supports effective monitoring of water environments
- lacktriangle Excellent sensitivity, achieving quantification of eta-cyclocitral in water within the ultra-trace level
- SPME provides a solvent-free extraction, efficient, and highly sensitive method for detecting β-cyclocitral

■ Introduction

Cyanobacterial blooms, commonly known as algal blooms, cause rivers and lakes to turn vividly green due to the excessive growth of cyanobacteria. This phenomenon is caused by factors such as global population growth, industrial pollution, and rising water temperatures caused by climate change.¹⁾

Cyanobacterial blooms produce volatile organic compounds (VOCs) like β -cyclocitral and β -ionone, which cause unpleasant odors and reduce the quality of drinking water and aquatic products. This leads to consumer complaints and economic losses in aquaculture and water treatment.

Detecting and quantifying β -cyclocitral (Figure. 1) at trace levels is therefor essential for effective water management. However, conventional solvent extraction methods often fail to detect β -cyclocitral, as it requires acidification or heating to be effectively released from the matrix.²⁾

In contrast, solid phase microextraction (SPME) enables efficient detection by facilitating the thermal desorption of β -cyclocitral from sample matrices onto the fiber. Moreover, SPME reduces solvent use and shortens sample preparation time compared to traditional extraction techniques, making it a suitable approach for the analysis of thermally induced volatile compounds like β -cyclocitral.

This application news describes the suitability of β -cyclocitral analysis using SPME automatic injection (AOCTM-6000 Plus) and GCMS-QP2020 NX of Shimadzu. The developed method enables highly sensitive quantitation of β -cyclocitral, achieving detection at the pg/mL level.

$$O$$
 H H_3C CH_3

Figure 1. Structural Formula of β-cyclocitral

■ Analytical Conditions

β-cyclocitral in water was analyzed using GCMS-QP2020 NX of Shimadzu equipped with the AOC-6000 Plus autosampler as shown in Figure 2. The detailed analytical conditions of instrument was provided in Table 1. For the SPME analysis, 50/30 μm DVB/CAR/PDMS fiber was used, and SH-PolarWax (30 m × 0.25 mm l.D., 0.25 μm) was used as the analytical column.

The AOC-6000 Plus supports various injection techniques, including liquid injection, headspace (HS) injection, and solid-phase microextraction (SPME) injection. Furthermore, sample preparation and preprocessing in the AOC-6000 Plus are automatically carried out during injection into the GC system, improving efficiency and reproducibility.



Figure 2. GCMS-QP2020 NX with AOCTM-6000 Plus

Table 1. Analytical Conditions

AOC-6000 Plus		
SPME Fiber	: 50/30 µm DVB/CAR/PDMS	
Incubation Temp.	: 60 °C	
Incubation Time	: 5 min	
Sample Extract Time	: 20 min	
Sample Desorb Time	: 10 min	
Agitation	: On	
GCMS-QP2020 NX		
Analytical Column	: SH-PolarWax (30 m × 0.25 mm l.D., 0.25 µm)	
Column Temp.	: 50 °C (1 min) \rightarrow 10 °C/min \rightarrow 160 °C \rightarrow	
	$20 ^{\circ}\text{C/min} \rightarrow 240 ^{\circ}\text{C} (2 \text{min})$	
Carrier Gas Flow	: He, 1.0 mL/min	
Injection Mode	: Splitless	
Injector Temp.	: 260 °C	
Ion Source Temp.	: 230 °C	
Interface Temp.	: 250 °C	
Acquisition Mode	: SIM mode (<i>m/z</i> 152, 137)	

■ Materials and Preparation

Materials

 β -cyclocitral standard (97 % purity) was purchased from Sigma Aldrich. tert-Butyl methyl ether (MTBE, 99 % purity) used as a solvent was obtained from Samchun Chemical. Distilled water for the calibration curve was prepared using a Millipore Milli-Q water purification system.

Standard and sample preparation

A stock solution of β -cyclocitral was prepared at a concentration of 1,000 μ g/mL in MTBE. Six standard solutions were obtained by diluting the stock solution with MTBE to final concentrations of 1.0, 5.0, 10.0, 20.0, 50.0, and 100.0 ng/mL.

Each calibration standard was prepared by adding $10~\mu L$ of the corresponding standard solution to 10~mL of distilled water in 20~mL headspace vial, followed by 4~g of NaCl. The resulting final concentrations of calibration standard solutions were 1.0, 5.0, 10.0, 20.0, 50.0, and 100.0~pg/mL in distilled water.

Recovery test

To evaluate accuracy and precision, a sample solution with a concentration of 25 pg/mL was prepared and analyzed. The sample was prepared as follows: 10 mL of distilled water was placed in a 20 mL headspace vial, spiked with 10 μ L of a 25 ng/mL standard solution, and 4 g of NaCl was added. Four replicate measurements were performed, and the recovery percentage were calculated accordingly. Accuracy was assessed based on the mean recovery value, while precision was determined by calculating the relative standard deviation (%RSD) of the recovery values.

■ Results

Calibration curve

The calibration curve for β -cyclocitral was established using the external standard method over the concentration range of 1.0 – 100.0 pg/mL. Excellent linearity was observed, with the coefficient of determination (R²) better than 0.999. The calibration curves and analytical chromatograms at each concentration level are shown in Figure 3.

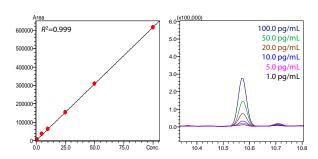


Figure 3. Calibration Curve and Chromatograms of $\beta\mbox{-cyclocitral}$

Accuracy and precision

The analytical method exhibits satisfactory accuracy and excellent precision in recovery test. As summarized in Table 2, the mean recovery was 105.3%, with %RSD of 0.7.

Table 2. Accuracy and Precision of β -cyclocitral (25 pg/mL, n=4)

_					
		Concentration (pg/mL)	Average (pg/mL)	Accuracy (%)	Precision (%RSD)
	1	26.2	26.3	105.3	0.7
	2	26.5			
	3	26.5			
	4	26.1			

■ Conclusions

In this application news, SPME-GC-MS method was developed and validated for the ultra-sensitive quantitation of β -cyclocitral in water using the Shimadzu GCMS-QP2020 NX coupled with the AOC-6000 Plus. The method demonstrated quantification capability down to the picogram-per-milliliter (pg/mL) level, excellent linearity (R² > 0.999). Accuracy and precision were confirmed with a recovery of 105.3% and %RSD of 0.7. These results confirm the reliability and suitability of the SPME-GC-MS method for ultra-trace quantification of β -cyclocitral in environmental water samples.

■ Reference

- Effects of different cultivation conditions on the production of β-cyclocitral and β-ionone in Microcystis aeruginosa, Jessica Aparecide Silva Moretto et al., BMC Microbiology volume 22, Article number: 78 (2022)
- Analytical Technique Optimization on the Detection of βcyclocitral in Microcystis Species: Molecules 2020, 25, 832; doi:10.3390/molecules 25040832

 $AOC\ is\ a\ trademark\ of\ Shimadzu\ Corporation\ or\ its\ affiliated\ companies\ in\ Japan\ and/or\ other\ countries.$



Shimadzu Corporation www.shimadzu.com/an/

Shimadzu Scientific Korea www.shimadzu.co.kr For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. See http://www.shimadzu.com/about/trademarks/index.html for details.

Third party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not

they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

09-SSK-022-EN

First Edition: Jul. 2025

> Please fill out the survey

Related Products Some products may be updated to newer models.



Related Solutions

