



ROI settings: adding attributes

Go to ROI settings

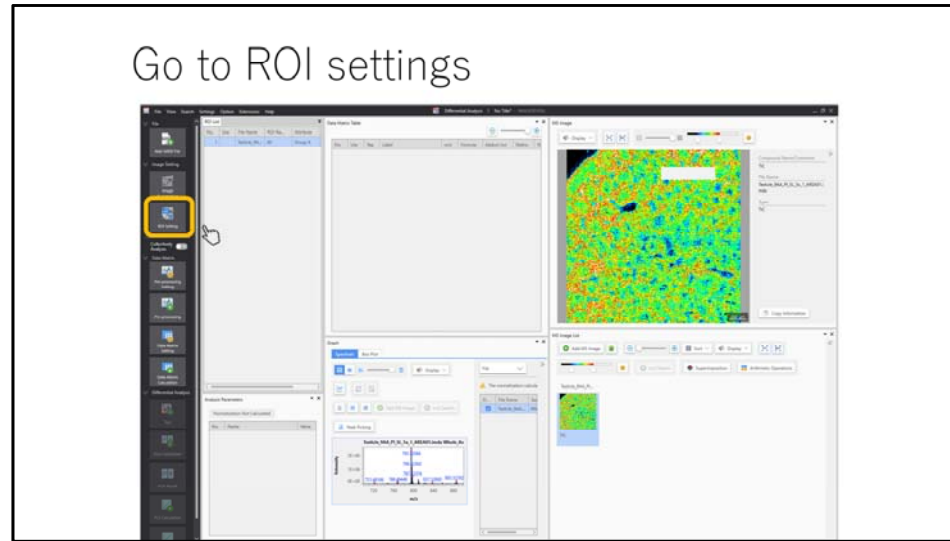


We will apply settings to the relevant ROI (Region Of Interest).

Apply ROI settings

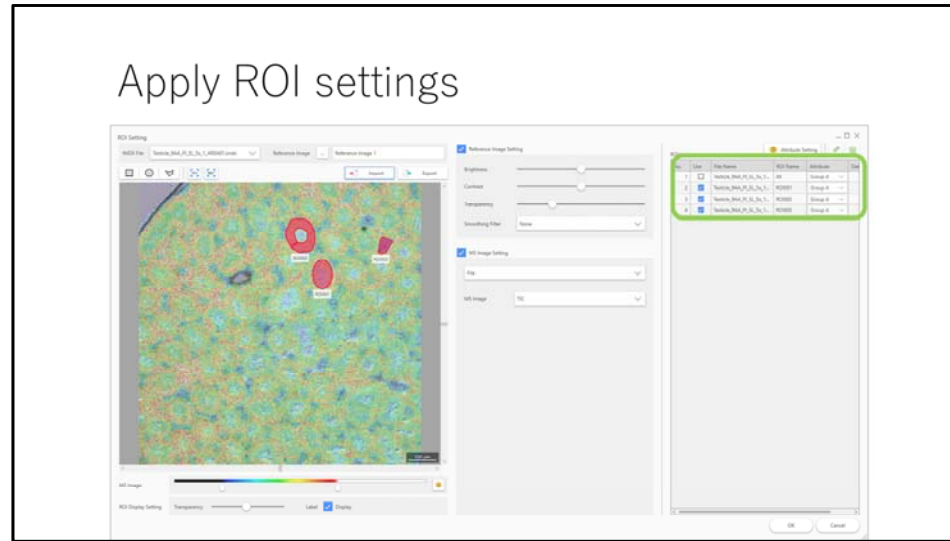

Attributes can be added after drawing the ROI.
*At least 2 attributes must be set for testing.


Add an attribute

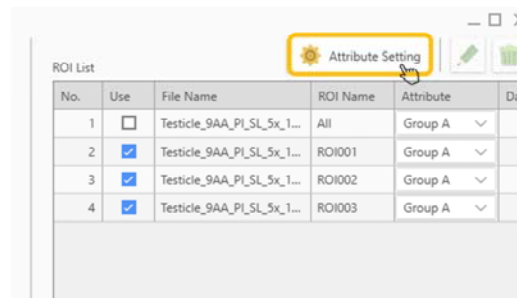

The screenshot shows a window titled "ROI List" with a table of ROI entries. A yellow box highlights the "Attribute Setting" button in the top right corner of the window. The table has columns for "No.", "Use", "File Name", "ROI Name", "Attribute", and "Date".

No.	Use	File Name	ROI Name	Attribute	Date
1	<input type="checkbox"/>	Testicle_9AA_Pi_SL_5x_1...	All	Group A	
2	<input checked="" type="checkbox"/>	Testicle_9AA_Pi_SL_5x_1...	ROI001	Group A	
3	<input checked="" type="checkbox"/>	Testicle_9AA_Pi_SL_5x_1...	ROI002	Group A	
4	<input checked="" type="checkbox"/>	Testicle_9AA_Pi_SL_5x_1...	ROI003	Group A	

Add an attribute

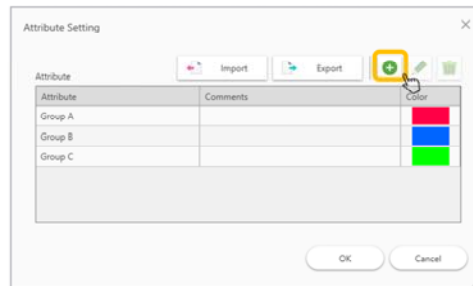

Enter your preferred text in the "Attribute" box

The image shows a dialog box titled "Edit Attribute" with a close button (X) in the top right corner. It contains three main sections: "Attribute" with a text input field highlighted by a green border, "Comments" with a text input field, and "Color" with a red square. At the bottom, there are "OK" and "Cancel" buttons.


Set the colour for the attribute that will be applied to the ROI

The image shows a software dialog box titled "Edit Attribute" with a close button (X) in the top right corner. It contains three input fields: "Attribute" with the text "Center", "Comments" which is empty, and "Color" which displays a red square. The "Color" field is highlighted with a green border. At the bottom of the dialog are "OK" and "Cancel" buttons.

After entering this information, click “OK”

The image shows a dialog box titled "Edit Attribute" with a close button (X) in the top right corner. It contains three input fields: "Attribute" with the text "Center", "Comments" which is empty, and "Color" which shows a purple square. At the bottom, there are two buttons: "OK" and "Cancel". The "OK" button is highlighted with a yellow border, and a mouse cursor is pointing at it.

The new attribute has been added

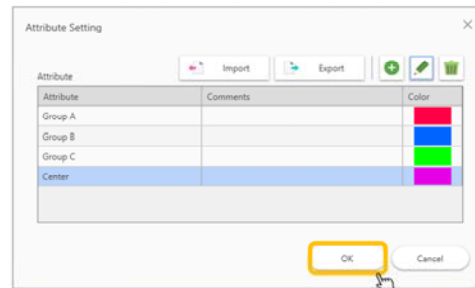

Now the new attribute can be selected from the drop-down list

