CoreFocus

P/N : 227-35032-01

Inlet Liner Removal Tool

INSTRUCTION MANUAL

The Inlet Liner Removal Tool is designed to simplify and efficiently facilitate the removal of gas chromatography (GC) inlet liners.

Applicable liner outer diameter : 5 mm

Recommended operating temperature : Room temperature


CAUTION

Please use this tool only after SPL has cooled down to room temperature.

There is a risk of burns when SPL is at high temperatures. Additionally, using this tool at high temperatures may accelerate its deterioration.

【The procedure for removing the liner】

1. While holding down the injection port, turn the lever counterclockwise.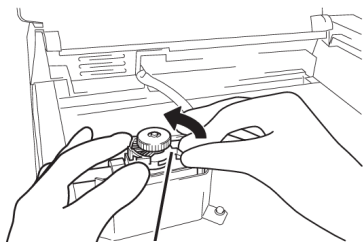


Lever

2. Slowly lift the injection port.

Note

Please lift the injection port slowly and directly upwards with your hand to prevent damage to the glass liner.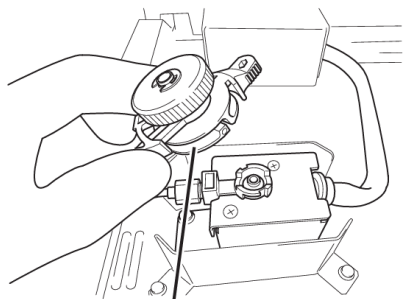

Injection Port


Tips

If the O-ring is stuck on the injection port, please use a pair of tweezers to carefully remove the O-ring before lifting the injection port to prevent any damage to the glass liner.

3. Insert the tool into the liner that is protruding from the SPL.


Note

Please be careful not to damage the sealing surface of the O-ring on the sample injection port.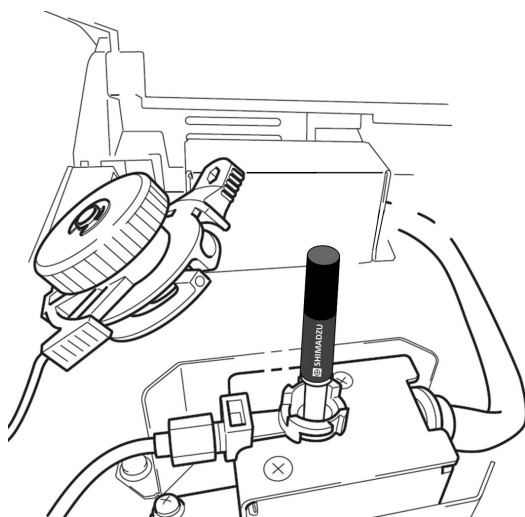

Tips


If the O-ring is stuck due to heat, please use a pair of tweezers to carefully tear the O-ring away from the main body via the two grooves on both sides.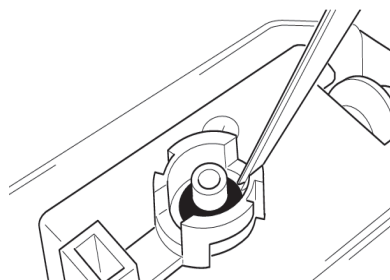

4. Pull it straight up to remove it.

Note

There is a possibility of the liner breaking if you pull it out diagonally.

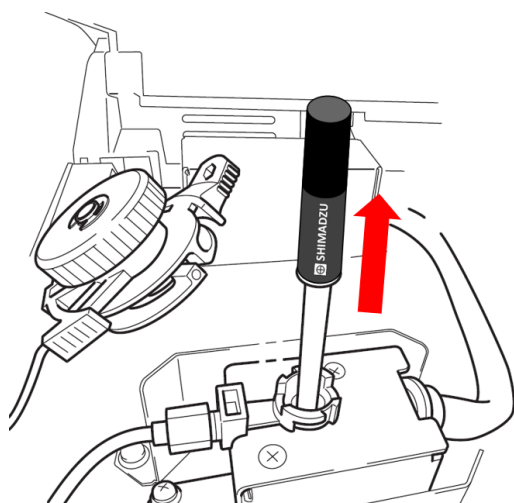5. Remove the tool from the extracted liner.

Note

Please be careful as the liner may be hot.

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