

Instruction Manual For Medium-Sized Electronic Balance

BX-K series

BX32KS, BX52KS, BX12KH, BX22KH, BX32KH

BW-K series

BW32KS, BW52KS, BW12KH, BW22KH, BW32KH

For Basic Operation



Read the instruction manual thoroughly before you use the product.
Keep this instruction manual for future reference.

 **SHIMADZU CORPORATION**
KYOTO JAPAN

ANALYTICAL & MEASURING INSTRUMENTS DIVISION

Requests

- Provide this manual to the next user in the event that the instrument is transferred.
- To ensure safe operation, contact your Shimadzu Balance representative for installation, adjustment, or reinstallation after moving the instrument to a different site.

Notices

- The content of this manual is subject, without notice, to modifications for the sake of improvement.
- Every effort has been made to ensure that the content of this manual was correct at the time of creation. However, in the event that any mistakes or omissions are discovered, it may not be possible to correct them immediately.
- The copyright of this manual is owned by Shimadzu Corporation. Reproduction and duplication of whole or part of the content without permission of the company are strictly prohibited.
- "Microsoft", "Windows", and "Excel" are registered trademarks of Microsoft Corporation of the U.S.A. in the United States and other countries. All other company names and product names that appear in this manual are trademarks or registered trademarks of the companies concerned. Note that ™ and ® indications are not used.
- The company names, organization names and product names in this manual are trademarks or registered trademarks of the companies and organizations concerned.
- Shimadzu does not guarantee that the WindowsDirect communication function will operate without problems on all PCs. Shimadzu will accept no responsibility for any trouble that arises as a result of using this function. You are recommended to back up all important data and programs in advance.



© 1997-2018 Shimadzu Corporation. All rights reserved.

Safety Precautions

To be strictly observed






To ensure that you use the balance safely and correctly, read the following precautions carefully and observe them.

The levels of danger and damage that will arise if the balance is used incorrectly are classified and indicated as shown below.

 WARNING	Indicates a potentially hazardous situation which, if not avoided, could result in serious injury or possibly death.
 CAUTION	Indicates a potentially hazardous situation which, if not avoided, may result in minor to moderate injury or equipment damage.

Precautions are classified and explained by using one of the symbols below, depending on the nature of the precaution.

 Instructions	Indicates an action that must be performed.
 Prohibitions	Indicates an action that must NOT be performed.

 WARNING	
 Prohibitions	Never disassemble, modify or attempt to repair this product or any accessory. You could sustain an electric shock or the product could operate abnormally. If you believe that the balance has failed, contact your Shimadzu representative.
 Instructions	Use the balance with the correct power supply and voltage. Use the balance with the attached AC adapter. Using the balance with an incorrect power supply or voltage will lead to fire or trouble with the balance. Note also that if the power supply or voltage is unstable or if the power supply capacity is insufficient, it will not be possible to obtain satisfactory performance from the balance.
 Instructions	Use the correct weighing units. Using incorrect weighing units can lead to accidents as a result of weighing errors. Check that the weighing units are correct before starting weighing.
 Prohibitions	Do not use the balance outdoors or anywhere where it will be exposed to water. You could sustain an electric shock or the product could operate abnormally.

⚠ CAUTION




Prohibitions

Avoid locations where the balance will be exposed to any of the following.

This could cause accidents or poor performance.

- Air flow from an air conditioner, ventilator, door or window
- Extreme temperature changes
- Vibration
- Direct sunlight
- Corrosive or flammable gases
- Dust, electromagnetic waves or a magnetic field

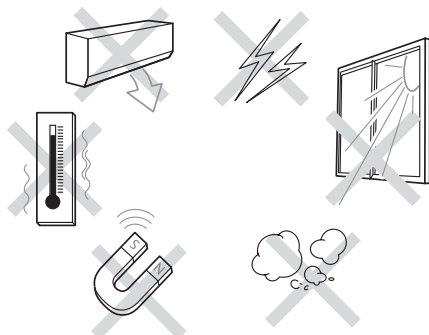

Instructions

Install the balance on a strong and stable flat table or floor in the room.

Placing the balance in an unstable site could lead to injury or trouble with the balance. When selecting the installation site, take into account the combined weight of the balance and the item to be weighed.



Instructions

After a power outage, turn the power back ON.

When a power outage occurs, the power is shut off automatically. Therefore, begin operation from 1.3 "Turning the Power ON" (^ page 7) again.



Instructions

Treat the balance with care and respect.

The balance is a precision instrument. Subjecting it to impacts could cause it to fail. When moving the balance, remove pan and pan supporter. Grasp it firmly with both hands to carry it.

If the balance has to be stored for a long time, store it in the packaging box in which it was delivered.



Prohibitions

Do not connect anything other than peripheral devices specified by Shimadzu to the balance's connector.

If you do, the balance may stop working normally.

In order to avoid trouble, always connect peripheral devices in accordance with the directions in this manual.



Instructions

If you detect anything abnormal (e.g. a burning smell) disconnect the AC adapter immediately.

Continuing to use the balance with an abnormality could lead to fire or an electric shock.



Prohibitions

Do not use balance for other than weighing.

Do not use the balance for other than weighing.

Do not use the attached AC adapter for other equipment.

It may cause fire or malfunction.

In addition, the balance and the accessories are manufactured for weighing and industrial use, not suitable for home use.

Shimadzu Balances and 21 CFR Part 11

21 CFR Part 11

21 CFR Part 11, Electronic Records, Electronic Signatures, Final Rule (often referred to as Part 11) is the United States Food and Drug Administration (FDA) regulation affecting computer resources and electronic records that are used for any document that is required to be kept and maintained by FDA regulations.

Requirements concerning computer resources security are key elements in Part 11.

The controls implemented as a result of security related requirements are intended to result in trusted records.

Shimadzu CLASS-Balance Agent

Shimadzu provides a means for compliance with 21 CFR Part 11 with Shimadzu CLASS-Balance Agent software, part of a comprehensive laboratory data management system, Shimadzu CLASS Agent.

Ask your Shimadzu representative about it.

Shimadzu WindowsDirect

When Shimadzu balances are integrated with laboratory software by means of our WindowsDirect function, no communication software is required or used.

The Shimadzu balance functions as a primary device in the system, just as a keyboard, mouse or other data entry hardware does.

For this reason, system validation and compliance may be greatly simplified with the use of Shimadzu balances.

Two-way Communication

Shimadzu balances have always been computer friendly and they can be set up for bi-directional communication as part of a fully automated production system or LIMS.

This manual includes the command codes and information needed by programmers to integrate Shimadzu balances with their software.

Action for Environment (WEEE)

To all user of Shimadzu equipment in the European Union:

Equipment marked with this symbol indicates that it was sold on or after 13th August 2005, which means it should not be disposed of with general household waste. Note that our equipment is for industrial/professional use only.

Contact Shimadzu service representative when the equipment has reached the end of its life.

They will advise you regarding the equipment take-back.

With your co-operation we are aiming to reduce contamination from waste electronic and electrical equipment and preserve natural resource through re-use and recycling.

Do not hesitate to ask Shimadzu service representative, if you require further information.








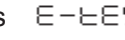










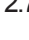

WEEE Mark

INTRODUCTION

Thank you for purchasing this medium-sized electronic balance, the Shimadzu BX-K/ BW-K series.

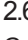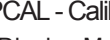In addition to its ability to make rapid mass measurements, the BX-K/ BW-K series is also well-equipped with a four-mode analog display, clock function, various application measurement functions and a unit conversion function. Furthermore, it is a high performance, multi-functional balance which can be used in a great variety of applications when connected to peripherals. The BW-K series features a built-in calibration weight which allows easy sensitivity calibration wherever you are, making it even more convenient to use.

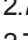Before using the balance read this instruction manual carefully to ensure long and fruitful use of the BX-K/ BW-K series.

1.	BASIC OPERATION.....	1
1.1	Balance Components.....	1
1.1.1	Operations of the Keys on the Main Unit.....	3
1.1.2	Symbol Display.....	4
1.2	Installation.....	5
1.2.1	Choosing the Installation Site.....	5
1.2.2	Balance Installation.....	5
1.3	Turning the Power ON.....	7
1.4	Adjusting the Built-in Clock.....	8
1.5	Span Calibration.....	8
1.6	Weighing.....	8
1.6.1	Changing the Unit Display.....	9
1.6.2	Changing the Minimum Display Digit ( / ).....	9
1.7	Maintaining and Transporting the Balance.....	9
1.7.1	Cleaning the Balance.....	9
1.7.2	Transporting the Balance.....	9
1.8	Specifications.....	10
2	MEMU ITEM SELECTION.....	11
2.1	Introduction.....	11
2.2	Procedure of Menu Item Selection.....	11
2.3	Menu Map.....	13
2.4	General Menu Operations.....	21
2.4.1	Setting Numeric Values.....	21
2.4.2	Setting a Decimal Point.....	21
2.4.3	Menu Lock.....	21
2.4.4	Last Menu Recall.....	22
2.4.5	Returning to the Default Settings (menu reset).....	22
2.5	Calibration Execution Menu.....	23
2.5.1	Calibration.....	23
2.5.2	Span Calibration Using the Built-in Weight. 	24
2.5.3	Calibration Check Using the Built-in Weight 	24
2.5.4	Span Calibration Using External Weights 	25
2.5.5	Calibration Check Using External Weights 	26
2.6	Calibration Setting Menu.....	26
2.6.1	Selecting the Standard Calibration Type ( - ).....	26
2.6.2	PCAL - Calibration of the Built-in Weight  - 	27
2.7	Graphic Display Menu  Std : 	28
2.7.1	Overview.....	28
2.7.2	Full Scale Mode  - 	28
2.7.3	Target Mode  - 	28
2.7.4	Group Mode  - 	29

2.7.5	Level Mode	■ - F t G L n o	29
2.7.6	No Graphic Display	■ - F t G L n o	29
2.8	Environment Menu	■ S t d : E A U S	30
2.8.1	Overview		30
2.8.2	Averaging	E - A b t P 8	30
2.8.3	Stability Detection Band	E - A b t P 8	30
2.8.4	Tracking	E - A b t P 8	31
2.8.5	Pretaring Value	⚡ E - A b t P 8	31
2.8.6	Whole Lighting Mode	E - A b t P 8	32
2.8.7	Taring/Printing	⚡ E - A b t P 8	32
2.9	Application Menu	S t d : E A U S	33
2.9.1	Overview		33
2.9.2	Zero Range		33
2.9.3	Auto Zero		34
2.9.4	Auto Print		34
2.9.5	Peak Hold		35
2.9.6	Interval Timer		36
2.9.7	Auto-Memory and Zeroing		37
2.9.8	Animal Weighing		38
2.10	Unit Registration Menu and Unit Change		39
2.10.1	Unit Registration Menu		39
2.10.2	% Conversion		39
2.10.3	Piece Counting		40
2.10.4	Solid Specific Gravity Measurement		41
2.10.5	Liquid Specific Gravity Measurement		42
2.11	System Setting Menu	■ S t d : E A U S	43
2.11.1	Date for the Built-in Clock		43
2.11.2	Time for the Built-in Clock	S - d t S C r	43
2.11.3	Display During Standby	S - d t S C r	43
2.11.4	Measurement Control System	S - d t S C r	44
2.12	Communication Menu	S t d E A U S (⚡ blinking)	44
2.12.1	Overview		44
2.12.2	Handshaking	⚡ - H F b P S d	45
2.12.3	Format	⚡ - H F b P S d	45
2.12.4	Communication Speed	⚡ - H F b P S d	46
2.12.5	Parity / Bit Length	⚡ - H F b P S d	46
2.12.6	Stop Bit	⚡ - H F b P S d	46
2.12.7	Delimiter	⚡ - H F b P S d	46
3.	Connecting Peripheral Instruments		47
3.1	Printer		47
3.2	Personal Computer - RS-232C -		47
3.2.1	Connecting the Cable		47

3.2.2	Data Format.....	48
3.2.3	Using Command Codes.....	48
3.2.4	Windows® Direct Function.....	54
3.3	Optional AKB-301 Application Keyboard.....	59
3.3.1	Basic Information.....	60
3.3.2	Example of Use.....	61
3.4	Foot Switches	63
3.5	Hook for Below-Balance Weighing.....	64
4.	APPENDIX	65
4.1	Part List.....	65
4.1.1	Standard Accessories and Maintenance Parts.....	65
4.1.2	Options.....	65
4.2	Table of Unit Conversion Constants.....	66
4.3	Specifications for the RS-232C/AUX Connector	67
4.4	Error Display List.....	68
4.4.1	General Display.....	68
4.4.2	Error Display.....	69

1. BASIC OPERATION

1.1 Balance Components

Open the package and sure following items are included.

Balance body, pan, AC adaptor, vinyl cover (*), adhesive tape,





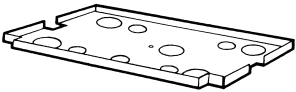

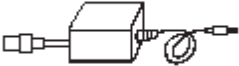
Instruction manual, inspection certificate,



* The adhesive tape may be attached on the vinyl cover.

<BW-K only>


Internal weight, Fix screw of the internal weight, cover of the internal weight

Fix of the screw of the internal cover

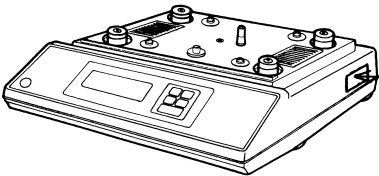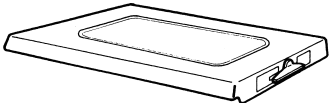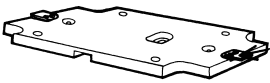[Name]	[Appearance]	[Notes]
Balance unit		The serial number of the balance is stated on this nameplate
Pan		
Built-in calibration mass (BWK series only)		Check that the numerals on this part are the same as the last 5 figures of the serial number on the balance unit
Calibration weight stopper screw (BWK series only)		
Calibration weight cover (BWK series only)		
Calibration weight cover setscrew (BWK series only)		
AC adaptor		※ The shape of the AC adaptor and the plug is subject to change without prior notice.

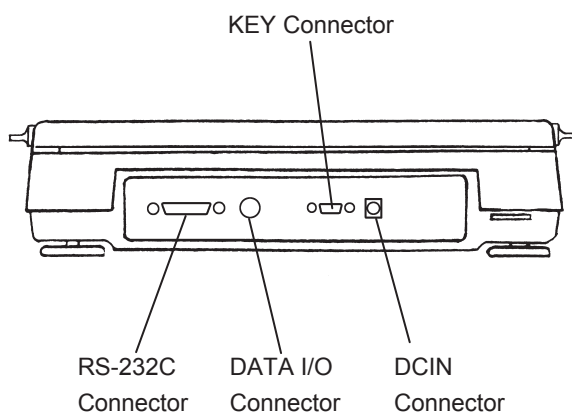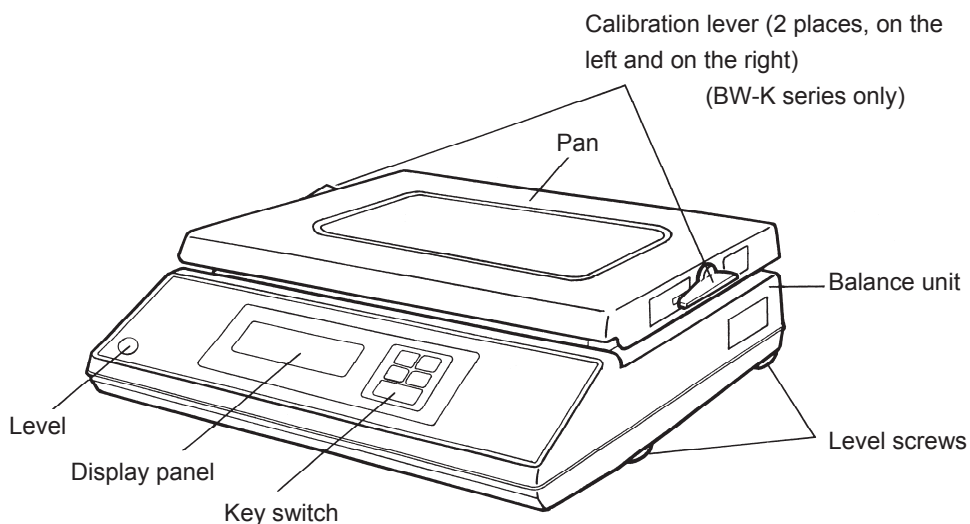

Position of the calibration lever (BW-K series only)

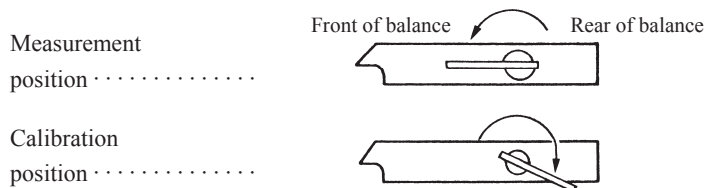

Regardless of whether the calibration lever is in the measurement or calibration position, ensure that both the left and right levers have been turned until they come to a stop.

1.1.1 Operations of the Keys on the Main Unit

The functions of the keys found on the front of the balance are described in the table below.

Key	During Weighing		During Menu Selection (* 1)	
	Pressing Once and Releasing	Pressing and Holding for About 3 Seconds	Pressing Once and Releasing	Pressing and Holding for About 3 Seconds
<u>POWER</u> BRK	Switches between the operation and standby modes.	Exits the Application Measurement function.	Returns to the previous menu.	Returns to the mass display.
<u>CAL</u> MENU	Enters span calibration or menu selection. (*2)	Displays the last Menu Map item that was set. (Last menu recall)	Moves to the next Menu Map item.	Displays the last Menu Map item that was set. (Last menu recall)
$\rightarrow 0/T \leftarrow$ ↵	Tares the balance. (Displays zero.) (*3)	Displays the Pretare value.	Selects and sets the currently displayed Menu Map item.	No operation.
<u>UNIT</u> ▲	Changes the selected unit. (*4)	Switches between the 1d and 10d display. (*5)	Increases the numeric value of the blinking digit by 1.	No operation.
<u>PRINT</u> ▶	Sends the displayed value to a peripheral device.	Sends the date and time to a peripheral device.	Moves to the next digit during numeric value entry.	No operation.

*1 Refer to Section 2. "MENU SELECTION."


*2 This key is used to set values when percent (%), number (PCS), solid specific gravity (▼d), or liquid specific gravity (d) units are displayed.

*3 When a Pretare value is set, zero is not displayed and the [- Pretare value] is displayed.

*4 Units other than "kg" must be registered and set before they can be used for measurement. Kirogram (kg) is the only unit is registered before shipment. To register other units, refer to Section 2.10. "REGISTRATION, RELEASE, AND SELECTION OF MEASUREMENT UNITS".

*5 When the unit is set to 10d, the resolution of the minimum display is decreased by one decimal place.

1.1.2 Symbol Display

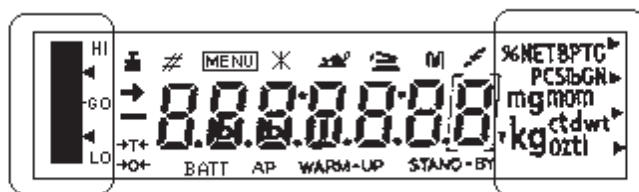

Graphic display

Unit display

An example of “whole lighting” display.

Display	Name	Description
	Stability symbol	Indicates that the balance is stable. (*1)
	Tare symbol	Indicates that a Pretare value has been set.
	Zero symbol	Indicates that Auto Zero is ON for the Application Measurement function.
	Weight symbol	Blinks when span calibration is necessary. This symbol continues to blink until either a manual or automatic span calibration has been made.
	Number symbol	Indicates numeric value entry.
	Menu symbol	Indicates that the menu lock is on.
	Asterisk	Indicates that the displayed numeric value is not a mass value.
	Animal symbol	Illuminates when Animal Weighing is ON for the Application Measurement function.
	Auto-Memory & Zeroing symbol	Illuminates when Auto-Memory and Zeroing are ON for the Application Measurement function.
	Memory symbol	Indicates that Net Total Weighing is ON for the Application Measurement function.
	Communication symbol	Illuminates during communication to external equipment through the RS-232C or DATA I/O connector.
BATT	Battery symbol	This symbol illuminates to indicate that the power voltage has dropped.
AP	Auto Print symbol	Indicates that Auto Print is ON for the Application Measurement function.
STAND-BY	Stand-by symbol	Illuminates when the balance power is in the standby mode. This symbol also illuminates when the Application Measurement function has entered the standby mode.
	Inverse triangle symbol	Illuminates when the solid specific gravity unit is used. This symbol is also used as a substitute for the decimal point.

*1 Stability symbol

The displayed value may change while the stability symbol remains illuminated if the load is changing slowly or if the stability detection band has been set to a large value.

1.2 Installation

1.2.1 Choosing the Installation Site

(1) Power supply



CAUTION

- Select an installation site that is near a power source to ensure that the attached AC adapter is used properly.
- Verify that the supply power voltage conforms to that indicated on the AC adapter.

(2) Installation site

Install the balance on a strong and stable flat table or floor in the room.



CAUTION



Choose an installation site where the balance will be protected from the following:

- corrosive or flammable gasses.
- dust, wind, vibration, electromagnetic waves, or magnetic fields.
- direct sunlight.
- extreme temperature or humidity.
- rain or the possibility of being splashed with water.

Large capacity balances should be installed on a sturdy floor and table that can withstand the total load of the balance AND object to be weighed.

1.2.2 Balance Installation


- 1) Put the balance onto its side and unscrew the 2 transportation screws in accordance with the instruction label on the bottom of the balance. Screw these into the “Release” hole.
- 2) Place the balance unit as shown in the diagram on the right. If intending to fit the vinyl cover, fix it at this stage of the procedure. Refer to “Fitting the Vinyl Cover”, mentioned later.

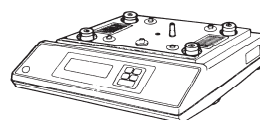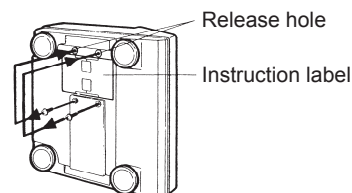


- 3) Fix the calibration weight cover (1) with the calibration mass cover setscrew (2).

Note


Steps (3) to (5) are only required with the BW-K series.

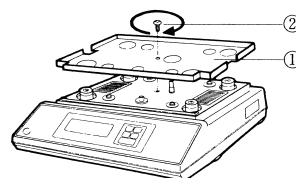


- 4) Put the built-in calibration weight (3) in place and tighten the calibration weight stopper screw (4).

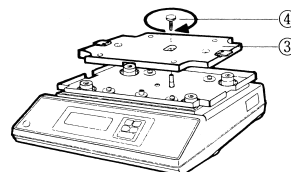

- 5) Turn both the left and right calibration levers of the pan to the Measurement position.

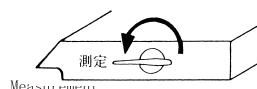

- 6) Place the pan on the balance. For the BW-K series, the pan will be easier to place if it is placed slightly from the rear.

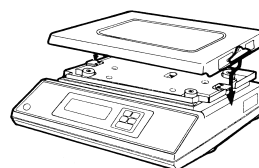

- 7) The level of the balance is adjusted ^{Note}.

Turn the level screws to adjust it so that the air bubble in the level rests inside the red circle. Check that the balance is steady when adjustment is complete.

Note


Adjusting the level can be done with ease if all 4 level screws are at first grounded equally. Next, adjust the front and the back using the 2 front feet, and then adjust left and right using either of the 2 feet on the left and right.

1.2.2.1 Fitting the Vinyl Cover

When using the balance in a location where it is likely to get dirty easily, refer to the diagram below and affix a vinyl cover by following the procedure below. Remove the pan first if fitting the vinyl cover after the balance has been assembled.

- 1) Cut the supplied tape (2) and stick it onto the 4 points on the case (3) (refer to the diagram below). (*)
* The adhesive tape may be attached on the vinyl cover.
- 2) Fix the vinyl cover (1) onto the double-sided tape (2), ensuring that there are no creases.
- 3) Fit the pan.

- 4) Place a weight onto the pan which is close to the weighing capacity to check that the pan and vinyl cover are not touching. Accurate measurements cannot be made if there is contact between the two items. If this is the case, re-fix the vinyl cover.

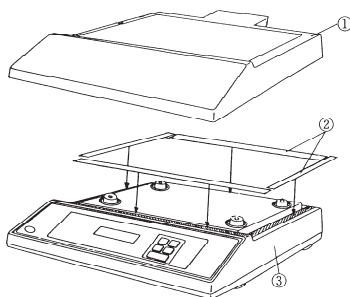

(Note) The vinyl cover (1) is folded up when it is in the packaging.

1.3 Turning the Power ON

- 1) Insert the plug of the AC adapter into the DC IN connector on the rear of the balance, then insert the AC adapter into the power source.
- 2) The balance self-check is activated and the following messages are displayed in the order indicated.

```
HELLO  ➡ CHE 5 ➡ CHE 4 ➡ CHE 3 ➡ CHE 2
➡ CHE 1 ➡ CHE 0 ➡ Whole lighting ➡ OFF
```

(CHE 5 and CHE 4 are not displayed for the BX series.)

- 3) When the $\frac{\text{POWER}}{\text{BRK}}$ key is pressed, the whole display illuminates and then the display changes to indicate the kilogram-display.

Note


If ENVIRONMENT menu 37 E8-STOP has been selected, the

display stops in the whole lighting state. In this case, press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key during whole lighting to change the display to the kilogram-display.

1.4 Adjusting the Built-in Clock

The BX-K and BX series of balances have a built-in battery driven clock. After unpacking and installing the balance, adjust this clock to the appropriate date and time.

 Refer to Section 2.11.1 “Date for the Built in Clock.”


 Refer to Section 2.11.2 “Time for the Built-in Clock.”

1.5 Span Calibration

It is necessary to calibrate the balance after it is moved, or when the ambient temperature has changed.

Verify that the balance is stable before performing the span calibration. To achieve a very stable state, ensure that the balance has been turned on for at least one hour, that the temperature is constant, that there are no breezes or vibrations and that the balance is in an area isolated from the normal traffic flow.

Refer to the following sections for more information on the Span Calibration procedure.

 For **BX-K series** balances, refer to Section 2.5.2 “Span Calibration Using the Built-in Weight.”

 For **BX series** balances refer to Section 2.5.4 “Span Calibration Using External Weights.”

1.6 Weighing




- 1) Place the weigh vessel (tare) on the pan.
- 2) Press the $\rightarrow 0/T \leftarrow$ key to zero the display.

Note

If a Pretare value is set, the  symbol is illuminated and zeroing of the display does not occur.

- 3) Place the object to be weighed on the pan, and read the displayed value after the stability symbol is displayed.

Errors Displayed During Weighing

	<i>Overload:</i> Weighing capacity has been exceeded.
	<i>Negative Overload:</i> The load on the balance is too light. The pan is not adjusted properly.
	<i>Display Overload:</i> This display appears if the mass display exceeds 7 digits due to the choice of unit. May appear after the $\frac{CAL}{MENU}$ key is pressed while the solid specific gravity (∇d) or liquid specific gravity (d) units are used.

1.6.1 Changing the Unit Display

Every time the $\frac{\text{UNIT}}{\blacktriangle}$ key is pressed, the unit display changes sequentially among those registered in the UNIT REGISTRATION menu.

Note: Before a unit can be displayed it must be registered in the UNIT REGISTRATION menu. Refer to “2.10 Unit Registration Menu and Unit Change” for information on how to register units.

The unit is displayed sequentially according to the order of the UNIT REGISTRATION menu.

1.6.2 Changing the Minimum Display Digit ($\text{□} \text{d} / \text{d}$)

It is possible to decrease the resolution of the minimum balance display by one decimal place if necessary.

- 1) Press and hold the $\frac{\text{UNIT}}{\blacktriangle}$ key for approximately three seconds. – $\text{□} \text{d}$ – is displayed and the display is decreased by one decimal place.
- 2) Press and hold the $\frac{\text{UNIT}}{\blacktriangle}$ key for approximately three seconds. – d – is displayed and the display returns to the original number of decimal places.

1.7 Maintaining and Transporting the Balance

1.7.1 Cleaning the Balance

Use a soft damp cloth containing a neutral detergent to clean the balance.

Avoid using organic solvents, chemicals, or dusting sprays as they may damage the coatings of the balance or the display panel.

Attach the in-use balance cover (standard accessory) when the balance is used in an environment where it is susceptible to being soiled.

The pan can be removed and washed with water. Verify that the pan is completely dry before replacing it on the balance.

1.7.2 Transporting the Balance

To carry the balance, hold the balance firmly with both hands.

To transport the balance, use the shipping carton used to deliver the balance.

Before placing the balance into the shipping carton, verify that the transportation screws at the bottom of the balance have been tightened until they stop.

1.8 Specifications

Type	Weighing Capacity (kg)	Minimum Display (g)	Standard Deviation ($\sigma \leq$)g	Linearity \pm g	Range of the External Calibration Weight (kg)
BX32KS BW32KS	32	1	0.6	1	5~32
BX52KS BW52KS	52	1	0.6	1	5~52
BX12KH BW12KH	12	0.1	0.1	0.2	5~12
BX22KH BW22KH	22	0.1	0.1	0.2	5~22
BX32KH BW32KH	32	0.1	0.12	0.2	5~32

The BX-K series is not equipped with a built-in calibration weight.

The BW-K series is equipped with a built-in calibration weight.

Temperature coefficient of sensitivity:

H-type = $\pm 4\text{ppm}/^{\circ}\text{C}$

S-type = $\pm 5\text{ppm}/^{\circ}\text{C}$ (Range of temperature compensation: 10 to 30°C)

Common Specifications

	BW-K	BX-K
Pan size (mm)	Approx.345 W x 250 D	
Dimensions	Approx.360 W x 355 D x 125 H	
Weight (kg)	Approx.16.5kg	Approx 10.5kg
Operating Temperature Range	5 - 40°C	
Rated electric power supply	DC12V, 300mA	
I/O Connector	RS-232C connector DATA I/O connector KEY connector	for RS-232C for Printer for application keyboard AKB-301
Pollution Degree	2	
Overvoltage Category	Category II	
Altitude	Up to 2000m	
Installation Site	device may only used indoors	
AC adaptor (Primary)	AC 100-240V, 480mA, 50/60 Hz	

2 MENU ITEM SELECTION

2.1 Introduction







The BX-W and BW-K series balance has many functions that can be selected to meet the requirements of the user. Menu Item selection is used to program these functions.


The  symbol is displayed during Menu Item selection.


Once the menu items have been set based on the installation environment and weighing purpose, it is not necessary to select the menu items each time the balance is used. Once the contents of the menu are set, they are stored even if the balance is turned OFF or if the power is disconnected.

2.2 Procedure of Menu Item Selection

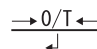The menu of the BX-W / BW-K balance consists of four levels.

- 1) Press the  key to cycle through the items within a menu level.
( in the table below)
- 2) Press the  key to choose the current item or move to the next menu level.
( in the table below)
- 3) Press the  key to move back one menu level.
- 4) Press and hold the  key to return to the mass display.

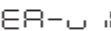

This instruction manual identifies each menu item by a number. For example, , “Anti-Vibration mode” is identified as [25].

 Refer to Section 2.3 "Menu Map" or the attached operation explanatory sheet.



Symbols in the Set : EAUS Display

Press the  key when the following symbols are blinking to enter the associated menu.

- | | |
|---|---|
|  (Calibration) | = Enters the menu related to span calibration. |
| (Graphic display) | = Enters the menu related to the graphic display, target and checkweighing. |
| E (Environment) | = Enters the menu related to the installation environment and general operation of the balance. |
| A (Application) | = Enters the menu related to applications and printing. |
| U (Unit) | = Enters the menu related to unit conversion. |
| S (System) | = Enters the menu related to the system (clock, reports, and passwords). |
|  (Communication) | =Enters the menu related to communication with a computer and software. |

The stability symbol in the Set : EAUS display indicates that EA-AUTO , Eb- I, and Et-on are set.

2.3 Menu Map

To enter the Menu Map:

Turn the balance on.

Press the $\frac{\text{CAL}}{\text{MENU}}$ key during mass display.

(This menu CAN NOT be entered from %, PCS, \blacktriangledown , and \square displays.)

The selected span calibration type is displayed. The possible displays are:

CAL\# (BW only)

EES (BW only)

E-CAL\# (BX or BW)

E-EES (BX or BW)

Press the $\frac{\text{CAL}}{\text{MENU}}$ key again to display the first menu level: Std : EAU

Press the $\frac{\text{CAL}}{\text{MENU}}$ key repeatedly to cycle through the items or press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key to select the blinking item.

Mass Weighing Display


\downarrow
 $\text{CAL\#} \rightarrow$ Performs selected standard span calibration.
 \downarrow

$\text{Std : EAU} \rightarrow$ Selects the blinking item.
 \downarrow

Moves to the next item

Default settings


Graphic Display Menu

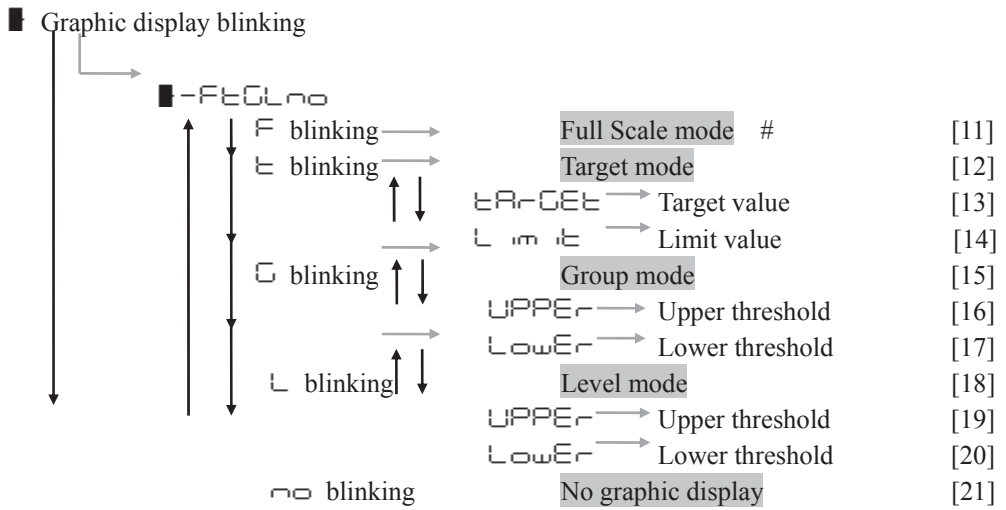

go to next page

Default settings


Environment Menu

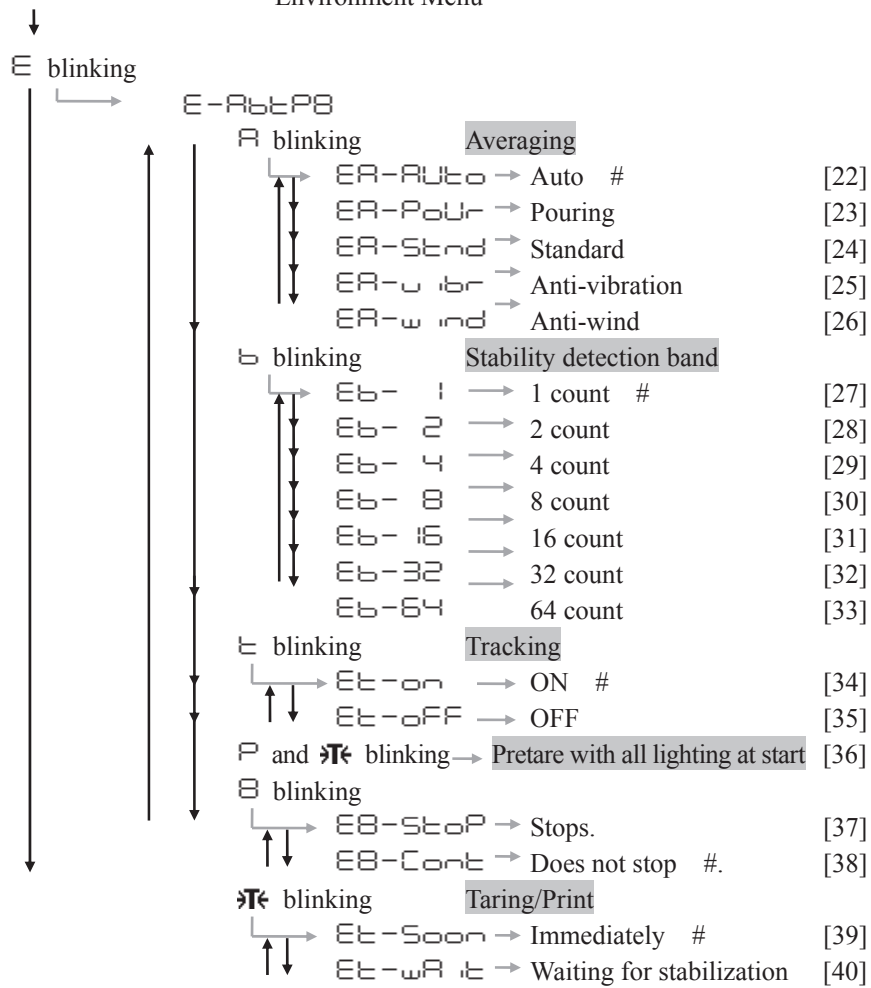

go to next page

Default settings

Applications Menu

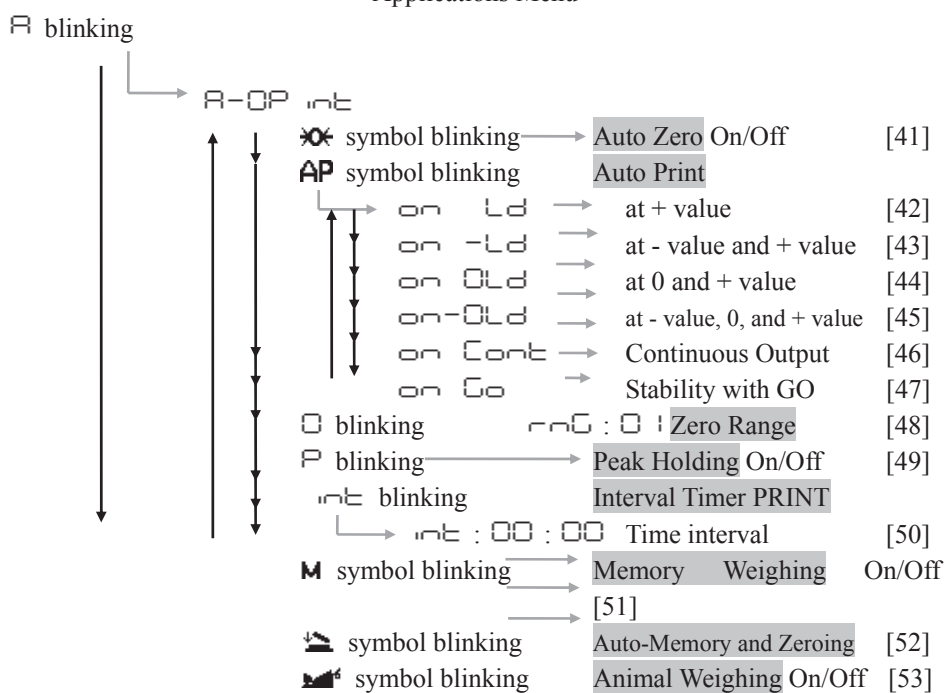


go to next page

Default Setting

Unit Registration Menu

U blinking

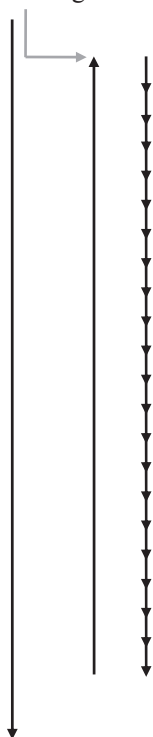


U-g	→	Gram (g)	[54]
U-mg	→	Miligram (mg)	[55]
U-%	→	Percent # (%)	[56]
U-pcs	→	Number # (PC)	[57]
U-ct	→	Carat (CT)	[58]
U-mom	→	Momme (MO)	[59]
U-□d	→	Immersion liq. density Solid density(DS)	[60]
U-d		Reference wt. volume Liquid density(DL)	[61]
U-Lb	→	Pound	[61a]
U-Oz	→	Ounce	[61b]
U-Ozt	→	Troy Ounce	[61c]
U-HK	→	Hong Kong's tael	[61d]
U-SPORE		Singapore tael	[61e]
U-TW		Taiwanese tael	[61f]
U-MAL	→	Malaysia tael	[61g]
U-CH		Chinese tael	[61h]
U-dwt	→	Pennyweight	[61i]
U-GN	→	Grain	[61j]
U-m	→	Mesghal	[61k]
U-b	→	Bahts	[61l]
U-t	→	Tola	[61m]
U-o	→	Parts Pounds	[61n]
U-USER	→	Set the Multiplier for User unit (US)	[62]

During
measurement,
these items are
distinguished
by the position

go to next page


Default settings

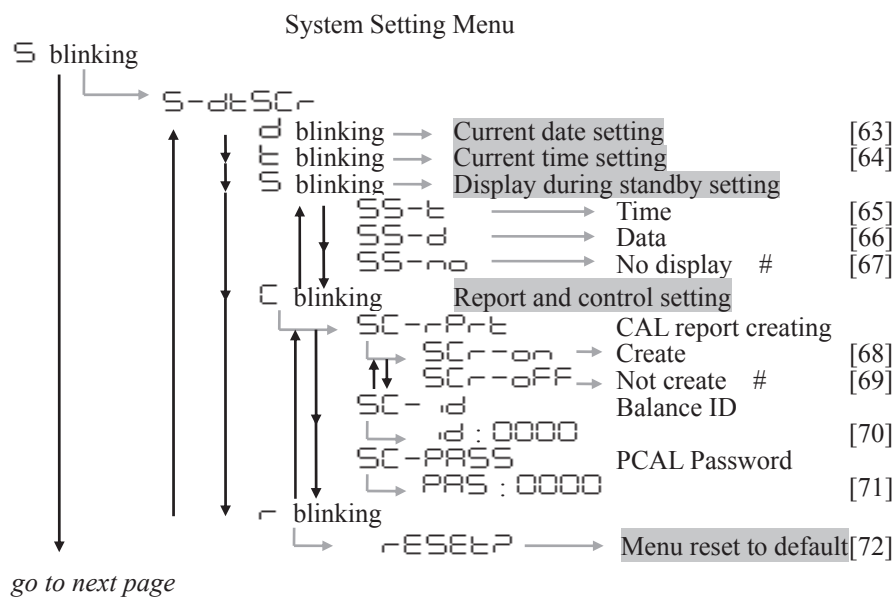

Note: Menu item [71] is not available with the BX-K balances.

Default settings

Communication Settings Menu

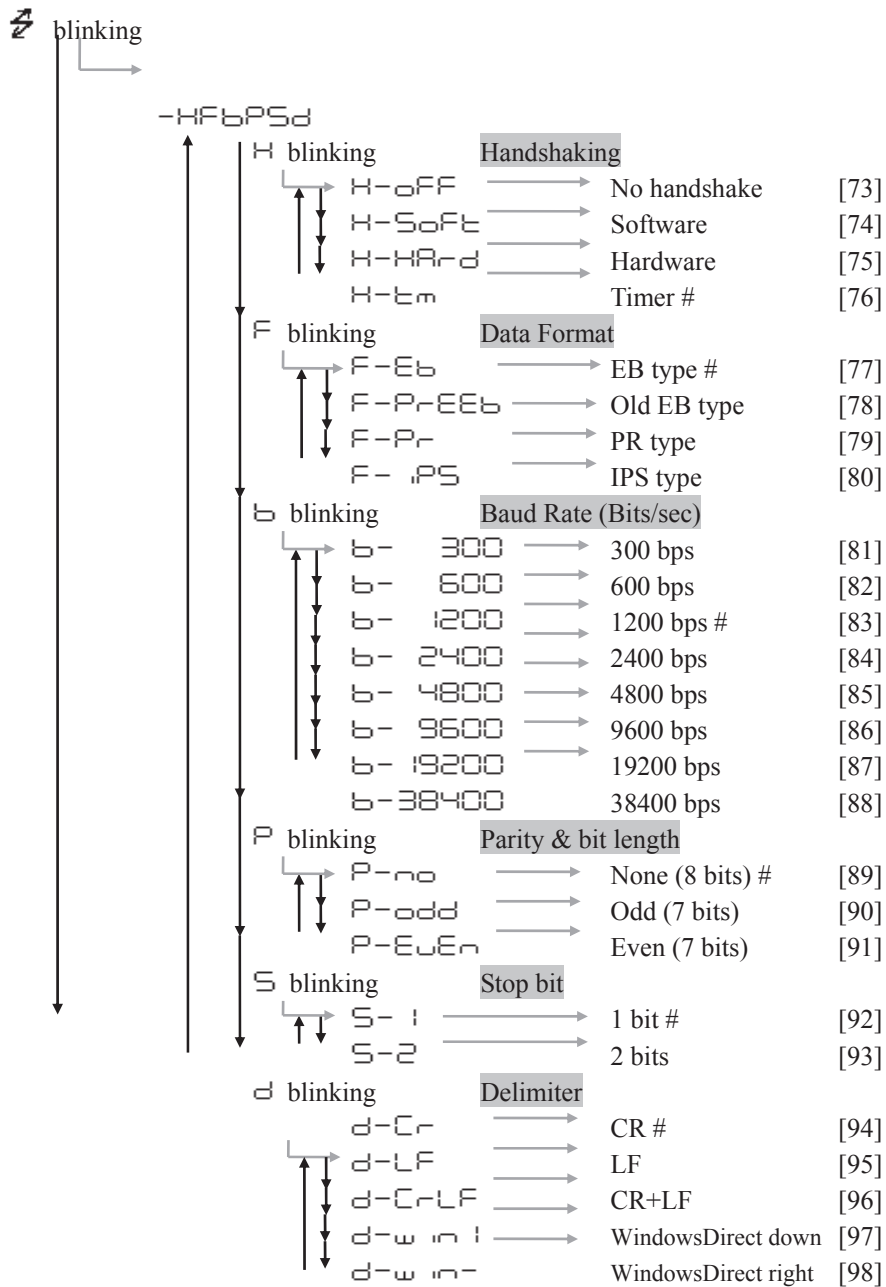

Returns to blinking  symbol.

Default settings

2.4 General Menu Operations

2.4.1 Setting Numeric Values

Numeric values may be used to set the threshold of the comparator with the BX-K and BW-K series balances.

Use the optional AKB-301 Application Keyboard to easily set these values. The values may also be set using only the balance keys.

In a menu used to set numeric values, **MENU** and **#** are both illuminated and the digit to be input blinks.

- Press the **UNIT** key to increase the value of the blinking digit by one. (0...9→
□)
- Press the **PRINT** key to move the blinking digit one place to the right.
- Press the **→0/T←** key to store the displayed value in the balance memory.
SEt is displayed when the value has been successfully saved.
Err is displayed when the balance failed to save the value.
- Press the **POWER** key to stop numeric entry. Abort is displayed briefly and
BRK the display returns to the menu, one level up.

2.4.2 Setting a Decimal Point

A decimal point is only used when setting units for solid density weighing, liquid density weighing or when setting the multiplier for the user defined unit. Set the decimal point while setting numerical values as follows.

- Press the **PRINT** key repeatedly until the last digit is blinking. Press the **PRINT** key
once more to initiate decimal point setting mode.
The ▼ symbol or current decimal point blinks.
- Press the **UNIT** key to move the blinking decimal point one digit at a time to the
desired position.
- Press the **→0/T←** key to set the decimal point position.
SEt is displayed briefly to indicate that the setting is completed.

2.4.3 Menu Lock

The BX-K and BW-K series of balances have a "Menu Lock" function that locks the menu selections to avoid changes being made by mistake._

The menu lock is toggled ON and OFF by pressing the $\frac{\text{CAL}}{\text{MENU}}$ key during OFF display that appears after power is supplied to the balance.

LOCKED is displayed when the menu is locked.

RELEASE is displayed when the menu lock is turned off.

Menu lock is on.

Menu lock is off.



Note: Access is denied and ERR22 is displayed when the user attempts to select a locked menu.

Use the following procedure to turn off the menu lock function and restore access to the menus.

- 1) Disconnect power from the balance and wait 10 seconds. Reconnect power to the balance.
- 2) When OFF is displayed, press the $\frac{\text{CAL}}{\text{MENU}}$ key.
- 3) RELEASE is displayed to indicate that the menu lock has been turned off.

2.4.4 Last Menu Recall

This function is convenient when an application requires frequent changes to a specific Menu Map item.

During mass display or menu selection, press and hold the $\frac{\text{CAL}}{\text{MENU}}$ key for approximately three seconds. The last Menu Map item that was changed or set is displayed.

2.4.5 Returning to the Default Settings (menu reset)

The procedure below describes how to reset the menu and return to the default settings.

Default settings are indicated with the # symbol in the Menu Map. (See Section 2.3 "Menu Map").

Select Menu Map item [72] to reset the menu.

- 1) In the mass display, press the $\frac{\text{CAL}}{\text{MENU}}$ key repeatedly until the S of Std:ERUS blinks.
- 2) Press the $\rightarrow 0/T \leftarrow$ key.
The system setting menu is selected S-dtScr.
- 3) Press the $\frac{\text{CAL}}{\text{MENU}}$ key repeatedly until the r in S-dtScr is blinking.

- 4) Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key to display $\neg E S E \neg$ ("?" means "?").
- 5) Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key again.
 $\neg E S E \neg$ is displayed to indicate menu reset completion.
- 6) Press the $\frac{POWER}{BRK}$ key several times (or hold it for approximately 3 seconds) to return to the mass display.

2.5 Calibration Execution Menu

2.5.1 Calibration

Calibration is required to accurately weigh items with an electronic balance. Calibration should be performed:

- When the location of the balance is changed, even within the same room.
- When the room temperature changes considerably.
- Periodically, according to the quality control plan of the user.


Terms used in this manual:

Span Calibration:	Adjustment of the balance to specifications using two weight values; zero and an appropriate value for the balance capacity.
Calibration Check:	Comparing the current calibration mass reading to the calibration mass reading after the last span calibration.
PCAL:	Procedure to calibrate the built-in calibration mass to an external calibration mass standard. The PCAL procedure is password protected.

Factory Default Settings:

BX-K series	Calibration using the external weight. (E-CAL)
BW-K series	Calibration using built-in weights. (I-CAL)

The settings for the type of calibration can be changed with menu selection.

 Section 2.6.1 "Selecting Standard Calibration Type."

Note: The following displays may appear in the situations listed.

CAL E1	When it takes too much time to stabilize the balance.
CAL E2	When the zero point of the balance appears to have changed greatly.
CAL E3	When the balance calibration appears to have changed greatly or the wrong weight is loaded on the pan during calibration.

In all of these situations, span calibration or calibration check is not performed.

 Refer to Section 4.4.2 "Error Display."

2.5.2 Span Calibration Using the Built-in Weight, \rightarrow CAL

- 1) Verify that the balance is in mass display and that the pan is empty.
- 2) Press the $\frac{\text{CAL}}{\text{MENU}}$ key once, \rightarrow CAL is displayed
(If \rightarrow CAL is not displayed, return to mass display and select Menu Map item [1].)
- 3) Press the $\rightarrow 0/T \leftarrow$ key.
When \rightarrow CAL 2 is displayed, and the stability mark (\rightarrow) illuminates on the left in the display unit, turn both of the calibration levers on the left and right of the pan to the back until they stop. At this point the stability mark will temporarily disappear.
- 4) Press the $\rightarrow 0/T \leftarrow$ key when the stability mark lights up again.
- 5) The display will show \rightarrow CAL 1. Return the calibration levers to the measurement position. The stability mark will disappear temporarily when this is done.
- 6) Press the $\rightarrow 0/T \leftarrow$ key after the stability mark lights up again.
After awhile the SET display will appear, and CALend will be displayed.
Span Calibration will be complete once the balance goes to kg display.

2.5.3 Calibration Check Using the Built-in Weight \rightarrow TEST

- 1) Verify that the balance is in mass display and that the span is empty.
- 2) Press the $\frac{\text{CAL}}{\text{MENU}}$ key once, \rightarrow TEST is displayed.
(If \rightarrow TEST is not displayed, return to mass display and select Menu Map item [2].)
- 3) Press the $\rightarrow 0/T \leftarrow$ key.
When \rightarrow TEST 2 is displayed, and the stability mark (\rightarrow) illuminates on the left in the display unit, turn both of the calibration levers on the left and right of the pan to the back until they stop. When this is done the stability mark will disappear temporarily.
- 4) Press the $\rightarrow 0/T \leftarrow$ key after the stability mark lights up again.
The display will show \rightarrow TEST 1. Return the calibration levers to the measurement position. The stability mark will disappear temporarily when this is done.
- 5) Press the $\rightarrow 0/T \leftarrow$ key after the stability mark lights up again.
After awhile dxxx will be displayed (xxx will be numerals).
This value indicates the extent of the current deviation in the balance sensitivity.

- 6) When correcting the Δ value to zero, press the $\frac{\text{CAL}}{\text{MENU}}$ key.
Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key if not correcting it to zero (although pressing the $\frac{\text{POWER}}{\text{BRK}}$ key will also leave it uncorrected, this will in effect mean “Calibration Aborted”).

Note: Changing the Δ value to zero is equivalent to performing span calibration.

- 7) $\text{CAL} \Delta$ is displayed, indicating the completion of the calibration check.

Note: Examples for interpreting the results of a Calibration Check:

Δ Value	Actual Mass	Displayed Mass
-0.0003	30kg	29.9997 (32kg/0.1g balance)
+0.21	400g	400.21 (420g/0.01g balance)

Error codes that may be displayed:

- $\Delta \text{ OVER}$ indicates that the Δ value is 1000 counts or more.
- $\Delta \text{ UNDER}$ indicates that the Δ value is -1000 counts or less.

2.5.4 Span Calibration Using External Weights E-CAL

- 1) Verify that the balance is in mass display and unload the sample from the pan.
- 2) Press the $\frac{\text{CAL}}{\text{MENU}}$ key once. E-CAL is displayed.
(If E-CAL is not displayed, return to mass display and select Menu Map item [3].)
- 3) Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.

The value of the correct calibration mass to be loaded is displayed and blinks.



Changing the Calibration Mass to be Used

Pressing the $\frac{\text{CAL}}{\text{MENU}}$ key allows changes to the weight value. Modify the value using the $\frac{\text{UNIT}}{\blacktriangle}$ key and $\frac{\text{PRINT}}{\blacktriangleright}$ key, then press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key. To interrupt modification, press the $\frac{\text{POWER}}{\text{BRK}}$ key.

- 4) Load the indicated calibration weight and press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.
- 5) Shortly, zero display blinks. Unload the weight from the pan and press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.
 SEt is displayed briefly to indicate completion of span calibration.

2.5.5 Calibration Check Using External Weights $E-EESE$

- 1) Verify that the balance is in mass display and unload the sample from the pan.
- 2) Press the $\frac{CAL}{MENU}$ key once to display $E-EESE$.
(If $E-EESE$ is not displayed, select Menu Map item [4].)
- 3) Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.

The value of the correct calibration mass to be loaded is displayed and blinks.



Changing the Calibration Mass to be Used

Pressing the $\frac{CAL}{MENU}$ key allows changes to the weight value. Modify the value using the $\frac{UNIT}{\blacktriangle}$ key and $\frac{PRINT}{\blacktriangleright}$ key, then press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key. To interrupt modification, press the $\frac{POWER}{BRK}$ key.

- 4) Load the indicated calibration mass and press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.

The zero display blinks.

- 5) Unload the weight from the pan and press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.

The display changes to the $\square xxx$ display. (xxx indicates a numeric value)

- 6) To change the \square value to zero, press the $\frac{CAL}{MENU}$ key.

Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key to avoid changing the \square value to zero. Pressing the $\frac{POWER}{BRK}$ key interrupts calibration and does not change this value to zero.

Note: Changing the \square value to zero is equivalent to performing span calibration.

2.6 Calibration Setting Menu

2.6.1 Selecting the Standard Calibration Type (\blacksquare - $EEEP$)

Menu Map item [1]	\cdot blinking	Selects "Span calibration using the built-in weight."
Menu Map item [2]	\cdot E blinking	Selects "Calibration check using the built-in weight."
Menu Map item [3]	E blinking	Selects "Span calibration using external weights."

Menu Map item [4]

⏏ blinking

Selects "Calibration check using external weights."

2.6.2 PCAL - Calibration of the Built-in Weight










PCAL is used to calibrate the built-in weight to a standard calibration weight that is correctly adjusted, traceable and/or certified.



CAUTION

Use a correctly controlled, precise calibration mass for this procedure. If it is performed without a correct calibration mass, span calibration and calibration checks using the built-in mass may not be correct in subsequent operations.

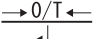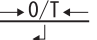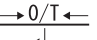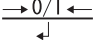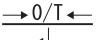- 1) Unload the sample from the pan and verify a zero mass display.
- 2) Select the Menu Map item [10].
PAS : 0000 is displayed.
- 3) Enter the PCAL password using the  and  keys, then press the  key.
The default password is 9999, set at shipment or upon menu reset.
- 4) After PCAL 3 is displayed, the value of the standard weight to be loaded blinks.
Load the standard mass displayed and press the  key.
- 5) Shortly, zero is displayed and blinks. Unload the weight and press the  key.
- 6) From this point the display will proceed automatically. When PCAL 2 is displayed, operate the levers on both sides of the pan to put on the built-in calibration weight.
- 7) Press the  key when the stability symbol re-illuminates after having disappeared for a moment.
- 8) When PCAL 1 is displayed, operate the levers on both sides of the pan to take off the built-in calibration weight.
- 9) Press the  key when the stability symbol re-illuminates after having disappeared for a moment.
- 10) From this point the display will proceed automatically. The procedure is complete when the balance goes to mass display mode.

Note: SEt is displayed during the process. Leave the balance in a stable state until the mass display appears as in step 6.

- In PCAL, the value of the "weight to be loaded" cannot be changed.
- Set the PCAL password using Menu Map item [71].

2.7 Graphic Display Menu ■SEd : ERUS

2.7.1 Overview

The BX-K and BW-K series balance has a bar graph located on the left side of the display. This graph can be conveniently used for liquid weighing and pass or failure judgment (comparator function) of the sample.

The graphic display functions include the Full Scale mode, Target mode, Group mode, Level mode, and no graphic display mode.

Note: Target, limit, upper, and lower values are set as numeric values only. Set the correct numeric value for the unit that will be used for weighing.

- In weighing mode, changing the displayed unit does NOT change the target, limit, upper, or lower numeric values. For example, when the upper limit has been set at 10g, switching the unit from "g" to "kg" does not change the limit to 10kg. It will be 0.01kg.
- The decimal point is invisible in the value setting display. Determine its position based on the resolution of the balance. For example, if the balance's minimum display is 0.01g, 150g must be set by inputting "15000" in the display.
- Numeric values are memorized independently for each graphic display mode.
- The graphic display mode can be selected without setting numeric values.

2.7.2 Full Scale Mode ■-FECLno

The relative amount of the load on the pan is displayed in the bar graph. This feature helps to prevent errors due to OL (overload) status.

Use Menu Map item [11] to select the Full Scale mode.

A bar displayed in the lower areas of the scale indicates that the load on the pan is small.

A bar displayed in the upper areas of the scale indicates that the load on the pan is close to the weighing capacity.

2.7.3 Target Mode ■-FECLno

Target mode is useful for filling applications or for quality check by weight.

The ERCEt value is the numeric value that is the desired amount in the unit that is used for weighing. The L m it value is the numeric amount above or below the target value that is acceptable.

This mode is useful for constant amount weighing of liquid or judgment of excess and shortage.

Menu Map item [12] selects the Target mode.

- Menu Map item [13] sets the $UPPER$ value, which corresponds to the center line of the graphic display.
- Menu Map item [14] sets the $LOWER$ value, which corresponds to the distance between the center line and upper or lower triangles.

2.7.4 Group Mode ■-FtCLno

This is the best mode to determine pass or failure judgment based on the sample weight.

Menu Map item [15] selects the Group mode.

Menu Map item [16] sets the upper threshold value, which corresponds to the upper triangle.

Menu Map item [17] sets the lower threshold value, which corresponds to the lower triangle.

Note:	Determination as follows:	<u>Display</u>
	$UPPER < \text{Sample weight}$	HI
	$LOWER \leq \text{Sample weight} \leq UPPER$	GO
	$\text{Sample weight} < LOWER$	LO

2.7.5 Level Mode ■-FtCLno

Use this mode for classification based on the sample weight. The display looks like a bar graph, but also includes a comparator function.

Menu Map item [18] selects the Level mode.

Menu Map item [19] sets the upper threshold value, which corresponds to the upper triangle.

Menu Map item [20] sets the lower threshold value, which corresponds to the lower triangle.

Note:	Determination as follows:	<u>Display</u>
	$UPPER < \text{Sample weight}$	HI
	$LOWER \leq \text{Sample weight} \leq UPPER$	GO
	$\text{Sample weight} < LOWER$	LO

2.7.6 No Graphic Display ■-FtCLno

Menu Map item [21] turns off the graphic display.

2.8 Environment Menu ■Std : EARS

2.8.1 Overview

Settings on the balance can be changed to compensate for the installation environment such as the degree of vibration etc. or for the purpose of weighing a solid, liquid or powder.

2.8.2 Averaging E-AVEPB

It is possible to match the stability of the display and the degree of response with the requirements of specific applications.

Menu Map item [22] EA-AUTO (Environment/Averaging-AUTOMATIC)
The balance automatically performs optimum averaging dynamically while observing the load data. This is the recommended setting and should be used unless special circumstances exist.

Menu Map item [23] EA-POUR (Environment/Averaging-POURING)
This mode is only suitable for constant volume weighing of liquids etc. This mode is very sensitive to wind and vibration.

Menu Map item [24] EA-STAND (Environment/Averaging-STANDARD)
This mode is suitable for weighing in a normal environment. Averaging is fixed and does not change dynamically as in EA-AUTO.

Menu Map item [25] EA-VIB (Environment/Averaging-VIBRATION)
Use this mode when the balance is used in a location where there are large vibrations and the display fluctuates in the EA-AUTO mode.
Response time is deteriorated by small mass amount changes.

Menu Map item [26] EA-WIND (Environment/Averaging-WIND)
Use this mode when the balance is used in a location where it is exposed to airflow that causes the display to fluctuate in the EA-AUTO mode.
Response time deteriorates further than EA-VIB, but weighing is comparatively stabilized.

Note: If weighing cannot be performed efficiently even with EA-WIND, change the installation site of the balance or use the deluxe windbreak or large pan windbreak.

2.8.3 Stability Detection Band E-AVEPB

Menu items [27] to [33] are used to determine the conditions for indicating balance stability.

EB-1 (Environment/Band-1):

When the display has remained constant (within one display count), the balance is regarded as stable and the stability symbol illuminates. Stability detection band settings can be selected from EB-2 to EB-64.

Note: Use $\text{E} \text{--} \text{B}$ to $\text{E} \text{--} \text{B} \text{--} \text{C}$ only when the Auto Print function is used and it is therefore required that the sensitivity to vibration be reduced to allow the balance to be regarded as stable and print the measurement value.

The balance may not operate properly if these settings are selected under normal use conditions and environment.

2.8.4 Tracking $\text{E} \text{--} \text{A} \text{--} \text{B} \text{--} \text{P} \text{--} \text{B}$

Tracking is the function that will maintain the current displayed value as long as possible.

Menu Map item [34] $\text{E} \text{--} \text{O} \text{--} \text{O}$ Turns ON this function.

Menu Map item [35] $\text{E} \text{--} \text{O} \text{--} \text{F} \text{--} \text{F}$ Turns OFF this function.



Zero Tracking Function

When the display is zero, $\text{E} \text{--} \text{O} \text{--} \text{O}$ menu functions as "zero tracking" to keep the zero display as long as possible. "Zero tracking" automatically cancels small zero drift.

It is recommended to set to $\text{E} \text{--} \text{O} \text{--} \text{F} \text{--} \text{F}$ when measuring slight mass change such as in the process of drop addition or liquid evaporation.

2.8.5 Pretaring Value $\text{P} \text{--} \text{T} \text{--} \text{E}$ $\text{E} \text{--} \text{A} \text{--} \text{B} \text{--} \text{P} \text{--} \text{B}$

This function is used to weigh the mass of a sample packed in a container such as a bottle or bag without opening the container. Pretare function should be used only if the mass of the containers does not vary from sample to sample. Instead of zero the pretare value is displayed (as a negative value) when the $\text{--} \text{O} / \text{T} \text{--}$ key is pressed. The pretare value is then subtracted from the load on the balance pan to determine the weight of the sample.

- 1) Menu Map item [36] opens the Pretare setting screen.
- 2) Set the Pretare value using the UNIT key and PRINT key, and press the

$\text{--} \text{O} / \text{T} \text{--}$ key.



Pretare Value

Cancel the Pretare value by setting the value to zero.

When a Pretare value other than zero has been set, $\text{P} \text{--} \text{T} \text{--} \text{E}$ illuminates.

The Pretare value is set using the "kg" unit. The maximum value is dependant on the weighing capacity of the balance.

Check the Pretare value by pressing and holding the $\text{--} \text{O} / \text{T} \text{--}$ key for approximately 3 seconds during weighing.

To set a pretare value using the optional AKB-301 Application Keyboard, press the [Pretare value] and the $\rightarrow 0/T \leftarrow$ key at the same time. Use the AKB-301 if you have to change the Pretare value frequently.

2.8.6 Whole Lighting Mode $E-ABtP8$

Press the $\frac{POWER}{BRK}$ key to switch from the standby mode to the mass display. The entire display is illuminated. This function sets whether the display automatically proceeds to mass display or waits for a $\rightarrow 0/T \leftarrow$ key command. When the mode has been set to $E8-STOP$, it is easier to check the display because the display remains illuminated.

Menu Map item [37] $E8-STOP$

Display stops at whole lighting. Press the $\rightarrow 0/T \leftarrow$ key to proceed to the mass display.

Menu Map item [38] $E8-CORt$

After the whole lighting display has appeared for approximately 0.5 sec, the display automatically proceeds to the mass display.

2.8.7 Taring/Printing $\overline{I}t E-ABtP8$

Determine if the balance should wait for stability before printing when the $\frac{PRINT}{\blacktriangleright}$ key is pressed or displaying the zero point when the $\rightarrow 0/T \leftarrow$ key is pressed.

When the Application Measurement function is selected, the $\frac{PRINT}{\blacktriangleright}$ key does not wait for stability.

Menu Map item [39] $Et-Soon$

Print and tare operate without waiting for stabilization.


Menu Map item [40] $Et-wA It$

Print and tare operate after the stability symbol illuminates.

Note: When the balance is in the $Et-wA It$ mode and waiting for stability to be determined, ---- is displayed when the $\rightarrow 0/T \leftarrow$ key is pressed.

Press the $\frac{POWER}{BRK}$ key to disable this function. Taring is not executed after the

$\frac{POWER}{BRK}$ key is pressed.

Once the PRINT key is pressed, data is not output until the  and symbol has illuminated and the balance is stable. Data is output only after the stability symbol is lit.

2.9 Application Menu Std : EAUS

2.9.1 Overview

The Application Measurement function is the generic name of the functions used to perform measurement work using the balance in more efficient ways. Only one Application Measurement function can be used at a time. When the power is turned ON, the balance will begin in the function mode that was last set.




Cancellation of Application Measurement

To cancel the Application Measurement, press the POWER
BRK key while in the mass display and hold it until APL End is displayed.

To use the Application Measurement function that was just canceled, use the last menu recall function by pressing and holding the CAL
MENU key for 3 seconds in the mass display.



Using the Application Measurement Function with a Unit Other Than "kg"

Select the Application Measurement function while in the mass display, and select the unit of weighing with the UNIT
 key.

Because the operation of the CAL
MENU differs from the one in normal mass display, the Application Measurement function cannot be selected while %, PCS, or specific gravity units are displayed.

2.9.2 Zero Range

The "Zero Range" value is used in the Application Measurement function as a reference to identify whether the sample is loaded.

For example, if the display is within the Zero Range, the balance determines that a sample is not loaded. If the display is more than five times the Zero Range it determines that a sample is loaded. Application Measurement functions which involves judgment of whether a sample is loaded or not work in accordance with the Zero Range setting.


In Menu Map item [48], the "Zero Range" value is set using the number of counts displayed in kilogram-display.

The setting range is 01 to 99 with 01 being the default value. Even when weighing will be done in another unit, Zero Range setting is made by only gram value.

Note: When a Pretare value is set, the value to determine that there is "no load" becomes "- Pretare \pm Zero Range" during mass display. For example, the Zero Range function works as expected when attempting to weigh and Auto Print the mass of a bottled sample during mass display.


2.9.3 Auto Zero

Select Menu Map item [41].



When the displayed value is within the Zero Range and the stability symbol has illuminated, zeroing occurs automatically. The  symbol appears in the display when the Auto Zero function is active.

Other keys function as expected with the Auto Zero function activated.


2.9.4 Auto Print

Auto Print function allows output of the data automatically without pressing the  key for each sample. The **AP** symbol is illuminated when the Auto Print function is activated.

Six types of Auto Print are possible.


Menu Map item [42]   (on positive load)

Load the sample when the value displayed is within the Zero Range. When the stability symbol has illuminated and the positive displayed value is more than 5 times the Zero Range, data is automatically output.

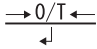The next data output is not performed unless the display has returned to a value within the Zero Range by unloading the sample or pressing the  key.


Menu Map item [43]   (on positive or negative load)

Load or unload the sample when the displayed value is within the Zero Range. When the stability symbol has illuminated and the displayed positive or negative value is more than 5 times the Zero Range, data is automatically output.

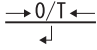The next data output is not performed unless the display has returned to a value within the Zero Range by unloading the sample or pressing the  key.


Menu Map item [44]   (on positive load and on zero)

Load the sample when the value displayed is within the Zero Range. When the stability symbol has illuminated and the positive displayed value is more than 5 times the Zero Range, data is automatically output.

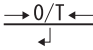

Unload the sample or press the  key. When the displayed value is within the Zero Range and the stability symbol has illuminated, data is output again.

Menu Map item [45]   (on positive or negative load and on zero)

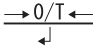
Load the sample when the value displayed is within the Zero Range. When the stability symbol has illuminated and the displayed positive or negative value is more than 5 times the Zero Range, data is automatically output.


Unload the sample or press the  key. When the displayed value is within the Zero Range and the stability symbol has illuminated, data is output again.

Menu Map item [46]   (on /continue)

By pressing the  key while the **AP** and **STAND-BY** symbols are lit, the **STAND-BY** symbol goes out, the  symbol lights and the displayed data is continuously output.

Continuous output stops temporarily when the  key is pressed.

Note: During continuous output, the  symbol may appear to remain lit. If the transfer speed of the data output is slow, the display may flash. Increase the transfer speed as much as possible and set the handshake to **HOFF**.

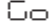Menu Map item [47]   (on / go)

When the graphic display is in Comparator mode (Group mode or Level mode) and the stability symbol lights in the GO range, the data is output once. The next data output is executed after the displayed value is within the Zero Range.

2.9.5 Peak Hold

Menu Map item [49] measures the displayed peak value. The **P** symbol is illuminated when the Peak Hold function is activated.

"Peak value" is the highest or lowest stable value displayed after the display has changed beyond five times the Zero Range.

1) In the peak detection standby state with the **P** and **STAND-BY** symbols illuminated, press the $\xrightarrow{0/T} \leftarrow$ key to tare the display.

2) Press the $\xrightarrow{\text{PRINT}} \leftarrow$ key.

The **STAND-BY** symbol disappears and peak value detection starts.

3) **P** and * are simultaneously displayed after the peak value is detected, and the data is output. This display will not change regardless of the load on the pan.

4) Press the $\xrightarrow{\text{POWER}} \leftarrow$ **BRK** key.

The balance returns to the peak detection standby state in step 1).

Note: Press the $\xrightarrow{\text{POWER}} \leftarrow$ **BRK** key in the peak detection standby state to initiate the power standby state.

Press the $\xrightarrow{\text{POWER}} \leftarrow$ **BRK** key during detection of the peak to return to the peak detection standby state.

Polarity of the peak value displayed is "polarity of the displayed value of the first change by five times or more of Zero Range from the display within Zero Range."

Usually the peak value is easily measured by setting to **EA-POUR** [23]. Depending on the weighting conditions and the sample type, this setting may not always be successful.

2.9.6 Interval Timer

Automatically outputs the displayed value at preset intervals. The **T** symbol is illuminated when the Interval Timer is activated.

The optional AKB-301 Application Keyboard is convenient when the interval is changed frequently.

1) Using Menu Map item [50], set the output interval (00:01 = 1 sec to 99:59 = 99 minutes 59 seconds).

2) In the interval timer standby state when the **T** and **STAND-BY** symbols are both illuminated, press the $\xrightarrow{\text{PRINT}} \leftarrow$ key.

The first data is output. Data is automatically output at the set time intervals.

3) To stop output, press the $\xrightarrow{\text{POWER}} \leftarrow$ **BRK** key.

The balance returns to the interval timer standby state in step 1).


Notes:


- Use the $\xrightarrow{0/T} \leftarrow$ key to erase the tare or zero the balance at any time.

- Pressing the $\frac{\text{POWER}}{\text{BRK}}$ key while in the interval timer standby state causes the power supply standby state to be initiated.
- To release the interval timer function, keep pressing the $\frac{\text{POWER}}{\text{BRK}}$ key. This does not reset the interval timer function to zero.
- Using the interval timer function to record data over a long period may cause data error due to balance drift.
- Some instruments receiving the data may not operate normally if the set time interval is short. To correct this, set the time interval to a longer period. When the set time interval is short and the instrument connected to DATA I/O is unknown, it is recommended to set the handshake to a setting other than H-T-m [76].

2.9.7 Auto-Memory and Zeroing

Use Menu Map item [51].

This menu is used to weigh a large number of individual samples. The  (Auto-Memory and Zeroing) symbol is illuminated when this function is active.

- 1) Load the weighing vessel and press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key in the Auto-Memory and Zeroing standby state ( and **STAND-BY** symbols are lit). Zeroing occurs.
- 2) Press the $\frac{\text{PRINT}}{\blacktriangleright}$ key.
The **STAND-BY** symbol disappears, and Auto-Memory and Zeroing measurement starts.
- 3) Load the first sample. Each time the stability symbol is lit and the display is a value five times or more the Zero Range or the $\frac{\text{PRINT}}{\blacktriangleright}$ key is pressed, the displayed value is output and zeroing occurs.
- 4) For the next sample, additional weighing is performed without pressing the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.
- 5) Press the $\frac{\text{POWER}}{\text{BRK}}$ key.
The balance returns to the Auto-Memory and Zeroing standby state and total mass on the pan without the tare is displayed.
Press the $\frac{\text{PRINT}}{\blacktriangleright}$ key to print this value.


Notes:

- When the stability symbol is illuminated and the displayed value is within the Zero Range, zeroing occurs automatically.

- When the $\frac{\text{PRINT}}{\blacktriangleright}$ key is pressed and the displayed value is five times the Zero Range, zeroing occurs after data output. (Manual loading)
- When the $\frac{\text{POWER}}{\text{BRK}}$ key is pressed in the Auto-Memory and Zeroing standby state, the power supply standby state is initiated.

2.9.8 Animal Weighing

Set with Menu Map item [53].

This function is suitable for weighing animals.  (animal) symbol is illuminated when the Animal Weighing mode is active.

- Note:**
- 1) Load the weighing vessel and press the $\frac{\rightarrow 0/T \leftarrow}{\blacktriangledown}$ key.
 - 2) Load the sample (animal etc.) with a mass more than 50 times the Zero Range.
 - 3) When the weighed value is relatively stable, the value is automatically output.
 - 4) Press the $\frac{\rightarrow 0/T \leftarrow}{\blacktriangledown}$ key or unload the sample.
 - 5) When the displayed value is stable and less than 10 times the Zero Range, automatic zeroing occurs.
- Any residue remaining on the pan (excrement or fur) is automatically canceled and zeroing occurs. If zeroing does not occur, increase the value of the Zero Range.

Notes:

- Standby state is not available in the Animal Weighing mode.
- Press the $\frac{\text{POWER}}{\text{BRK}}$ key to initiate the power supply standby state.
- On the premise of weighing animated objects, the stability detection band is automatically extended in the Animal Weighing mode. Reproducibility of the measurement data is slightly less than with other modes.
- When the animal being weighed cannot be controlled and Auto Print will not activate, press the $\frac{\text{PRINT}}{\blacktriangleright}$ key to output the displayed value. Then unload the animal. Even if the stability symbol lights before the animal is removed, data is not printed again.
- By setting a larger stability detection band in the menu, the stability symbol will light more readily.
- If the balance is slow to return to the zero point set a larger Zero Range value.

2.10 Unit Registration Menu and Unit Change

Std: EAUS

2.10.1 Unit Registration Menu

Press the $\frac{\text{UNIT}}{\blacktriangle}$ key in mass display to sequentially change the registered units. It is possible to display units other than "kg" with the BX-K and BW-K series of balances. Before weighing, set the display units to be used with the unit setting menu [54] to [62]. (Unit kg, %, number (PCS) were set before shipment.)

Note:

- In the unit setting menu, the stability symbol is illuminated to indicate the currently set units.
- Set or release the unit by pressing the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key when the unit is displayed.

Functional Units:

Solid specific gravity unit (\blacktriangledown d): Density of the liquid in which the sample is immersed.

Liquid specific gravity unit (d): Volume of the reference weight to be immersed in the sample.

USE- unit (\blacktriangleright): Numeric value (multiplier) multiplied by the gram (g) weight of the unit.

To register Functional Units, set the constant to something other than 0. Setting the constant to 0 turns the function off.

For the unit names, refer to  Section 2.3 "Menu Map."

2.10.2 % Conversion

- 1) Set the % unit with Menu Map item [56].
The % unit is set before shipment.
- 2) Press the $\frac{\text{UNIT}}{\blacktriangle}$ key several times in the mass display until the % unit is displayed.

Setting the 100% reference

Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key to tare the balance.

Load the reference sample that corresponds to the 100% value. This value must be equivalent to 100 counts or more in the "kg" unit.

When the stability symbol illuminates, press the $\frac{\text{CAL}}{\text{MENU}}$ key.

SEt is displayed briefly and the reference sample weight is displayed as 100%.

Note: The numbers of digits displayed in the % unit and rounding off of the minimum digit vary depending on the mass value of the reference sample and the balance model. It is not possible to obtain resolution greater than that in the "g" unit.

The weight of subsequent samples are displayed as a percentage of the reference sample weight.

2.10.3 Piece Counting

- 1) Register the PCS unit with Menu Map item [57].
(The PCS unit is registered before shipment.)
- 2) Press the $\frac{\text{UNIT}}{\blacktriangle}$ key several times in the mass display until the PCS unit is displayed.
- 3) Load the container and press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key to tare the balance.
- 4) Count exactly five pieces (or 10, 20, 50, 100, or 200 pieces) of sample to be measured and load them on the pan.
- 5) Press the $\frac{\text{CAL}}{\text{MENU}}$ key.
- 6) Every time the $\frac{\text{CAL}}{\text{MENU}}$ key is pressed, the display sequentially changes as
 $\text{Ld } 5\text{pcs} \rightarrow \text{Ld } 200\text{pcs} \rightarrow \text{Ld } 5\text{pcs} \rightarrow \dots$ (Piece menu).
 The default setting is $\text{Ld } 1\text{ pcs.}$
 Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key when the display is equivalent to the number of loaded pieces.

Example: If 50pcs. are loaded, press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key when $\text{Ld } 50\text{pcs}$ is displayed. This determines the unit weight or average weight per piece.
 As sample is added or removed, the piece count (number of pieces) is displayed.

Note: Repeat steps 3 through 6 above when the sample or manufacturing lot is changed.



Piece Count Menu Display a Next Setting

In this example, when the $\frac{\text{CAL}}{\text{MENU}}$ key is pressed in the next PCS menu, display starts from $\text{Ld } 50\text{pcs.}$



Counting Pieces Using the Optional AKB-301 Application Keyboard

By using the optional AKB-301 Application Keyboard, piece count based on an arbitrary number of pieces is possible.

The unit weight itself, if it is known, can be input using the numeric keys.

Section 3.3 "AKB-301 Application Keyboard".

2.10.4 Solid Specific Gravity Measurement

Solid specific gravity measurement refers to the measurement of the sample (solid) weight in the air and in a liquid of known density and the calculation of the sample density.

The ▼d symbol is used to represent the solid density unit in this balance. The data output unit is DS.

- 1) Set Menu Map item [60].

Note: Enter the value for the density (g/cm^3) of the liquid (water, alcohol etc.) in which the sample is immersed.

- 2) Attach the optional weighing hook to the bottom of the balance, attach the hanging pan, and then immerse the hanging pan in the tank filled with the liquid of known density.

- 3) Press the $\frac{\text{UNIT}}{\blacktriangle}$ key until ▼d is displayed.

- 4) Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.

- 5) Load the sample on the balance pan or in the hanging pan in air.

- 6) After the stability symbol illuminates, press the $\frac{\text{CAL}}{\text{MENU}}$ key.
dSPoL may be displayed but does not indicate a malfunction.

- 7) Load the sample on the hanging pan immersed in the liquid. The density of the sample is displayed.

- 8) Repeat steps 4 through 7 for each additional sample.

Note:

- Up to four decimal places are displayed for specific gravity. Since it may not be possible to stabilize the balance using all 4 places, 1d/10d switching is possible.
- When loading the sample on the pan in the liquid, ensure that the entire sample is immersed in the liquid.
- The balance does not re-zero when the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key is pressed in this unit display.
- When the optional Specific Gravity kit is used, refer to the Instructional Manual of the Specific Gravity kit for the procedure.
- Return the balance to kg display mode first before performing the next menu operation.

2.10.5 Liquid Specific Gravity Measurement

Liquid specific gravity measurement refers to the measurement of the weight of a reference solid of a known volume in air and in the sample liquid. Specific gravity for the liquid is calculated from these two values of.

The display unit for liquid specific gravity is "d." The data output unit is DL.

- 1) Set Menu Map item [61].

Note: Enter the value for the volume (cm^3) of the reference weight.

- 2) Attach the optional weighing hook to the bottom of the balance, attach the hanging pan, and then immerse the hanging pan in a tank containing the sample liquid.
- 3) Press the $\frac{\text{UNIT}}{\blacktriangle}$ key until "d" is displayed.
- 4) Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.
- 5) Load the reference weight on the pan of the balance.
- 6) After the stability symbol is illuminated, press the $\frac{\text{CAL}}{\text{MENU}}$ key.
dSPoL may be displayed but does not indicate a malfunction.
- 7) Load the reference weight on the hanging pan and immerse it in the sample liquid. The specific gravity of the sample liquid is displayed.
- 8) Repeat steps 4 through 7 for each additional sample.

Note:

- Up to four decimal places are displayed for specific gravity. Since it may not be possible to stabilize the balance using all 4 places, 1d/10d switching is possible.
- When loading the sample on the pan, immerse the entire sample in the liquid.
- To release the \blacktriangledown d, d or USER units, set the value to zero.
- When the optional Specific Gravity kit is used, refer to the Instructional Manual of the Specific Gravity kit for the procedure.

2.11 System Setting Menu ■ Set : EAUS

The SYSTEM SETTING menu is used to set the items that pertain to or are controlled by the balance.

2.11.1 Date for the Built-in Clock

- 1) Select Menu Map item [63].
- 2) Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.
- 3) Set the last two figures of the year, month and day, using the $\frac{\text{UNIT}}{\blacktriangle}$ and $\frac{\text{PRINT}}{\blacktriangleright}$ keys. Then press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.

Example:

February 1st, 1997	Set as 97. 02. 01
February 29th, 2004	Set as 04. 02. 29

Note:

- The built-in clock corrects for the leap year automatically.
- The moment the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key is pressed in step 2 above, seconds are set to zero.
If the date is set after setting the time, the second value will be incorrect. It is important to set the date first and then the time, or to correct the seconds value using the \pm second correcting function described later.

2.11.2 Time for the Built-in Clock S-dtSCr

- 1) Select Menu Map item [64].
- 2) Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.
- 3) Use the $\frac{\text{UNIT}}{\blacktriangle}$ key and $\frac{\text{PRINT}}{\blacktriangleright}$ keys to set the time in the 24 hour system, then press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key.

Example: 1 o'clock 23 minutes in the afternoon Set as 13 : 23

Note: The moment the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key is pressed is 00 seconds.

2.11.3 Display During Standby S-dtSCr

Determine what is to be displayed during a power supply standby.

Menu Map item [65]	<u>SS</u> - <u>t</u>	Displays the time during standby.
Menu Map item [66]	<u>SS</u> - <u>d</u>	Displays the date during standby.
Menu Map item [67]	<u>SS</u> - <u>no</u>	Displays nothing during standby.



Convenient Functions of Time Display

The following functions are available when the time is displayed during standby.
Seconds display function:

Press the $\frac{\text{UNIT}}{\blacktriangle}$ key to enable the display/non-display of the seconds value.

2.11.4 Measurement Control System S-dtSCr

Items related to the calibration of the balance and those set by the administrator are summarized in this menu.

2.11.4.1 Calibration Report SC-rPrt

Turns the calibration report function on/off. Use this to generate a calibration report as for GLP, GMP, or ISO9000. A printer (special accessory) is required to print the report.

Menu Map item [68] SC-on Calibration report is created.

Menu Map item [69] SC-off Calibration report is not created.

2.11.4.2 Balance ID SC-ID

Individual balances can be identified by the serial number on the main body of the balance. The user can add a four-digit ID number to the calibration report.

Set with Menu Map item [70]. Use a 4-digit number from 0000 to 9999.

2.11.4.3 PCAL Password SC-PASS

Enter a 4-digit number from 0000 to 9999 with Menu Map item [71].

It is recommended that the balance administrator set this password to prevent an unauthorized person from incorrectly calibrating the built-in calibration mass.

 Section 2.6.4 "PCAL - Calibration of the Built-in Weight"

Note: When the menu is reset, the PCAL password is reset to 9999.

2.12 Communication Menu Std EAUS (blinking)

2.12.1 Overview

This menu is used to set the specifications for communication between the balance and a personal computer or printer.

Note: This menu affects both the RS-232C and DATA I/O at the same time. For the instrument to be connected to the DATA I/O connector of a printer set the specifications for communication of the balance to the default settings, which are H-tm [76], F-Eb [77], b-1200 [83], P-no [89], S-1 [92], and d-Cr [94].

2.12.2 Handshaking - HFbPSd

Handshaking determines whether the peripheral equipment can receive communication data from the balance. This function does not relay the status of the balance to the peripheral equipment.

The balance is able to receive as long as there is space in the receiving buffer of the balance. This function operates once OFF is displayed, operation in other states is not guaranteed.

When the balance output is retained by handshaking, the display of the balance is locked. Determine the specifications for handshaking.

- | | |
|---|--|
| Menu Map item [73] H-OFF | Software handshaking is not performed. |
| Menu Map item [74] H-SOFF | Software handshaking is performed. |
| After the balance receives X-OFF (13H), the balance output is retained. | |
| After the balance receives X-ON (11H), the balance output is initiated. | |
| Menu Map item [75] H-HARD | Hardware handshaking is performed. |
| When DTR is OFF, the output from the balance is retained. | |
| When DTR is ON, the output from the balance is initiated. | |
| Menu Map item [76] H-ETM | Timed hardware handshaking is performed. |

2.12.3 Format - HFbPSd

Set the balance output format.

- | | |
|-----------------------------|---|
| Menu Map item [77] F-EB | The standard format for the Shimadzu electronic balance. |
| Menu Map item [78] F-PR-EEB | The old output format for the Shimadzu electronic balance |

The old output format is employed in the following models.

EB-500, 5000, 280, 2800, AEL-1600, EB-50K (except -15)

Note: In this format, the number of the lowest place of Menu Map item [70] SC-ID is assigned to identify the balance.

- | | |
|--------------------------|--|
| Menu Map item [79] F-PR | Compatible format for the PR and SR series of Shimadzu electronic balances. Available commands, functions and responses are limited. |
| Menu Map item [80] F-IPS | Compatible format for the IPS series of Shimadzu electronic balances. Available commands, functions and responses are limited. |

2.12.4 Communication Speed -HFbPSd

Select the communication speed (300, 600, 1200, 2400, 4800, 9600, 19200, or 38400 bps).

Number of b-xxx shows bps (bits/second). Baud rate and bps are the same value.

Set with Menu Map items [81] to [88].

2.12.5 Parity / Bit Length -HFbPSd

Select the parity and bit length.

Menu Map item [89]	P-no	No parity, 8-bit length
Menu Map item [90]	P-odd	Odd number parity, 7-bit length
Menu Map item [91]	P-even	Even number parity, 7-bit length

2.12.6 Stop Bit -HFbPSd

Select the number of stop bits.

Menu Map item [92]	S-1	Stop bit 1.
Menu Map item [93]	S-2	Stop bit 2.

2.12.7 Delimiter -HFbPSd

The "delimiter" is used to separate individual pieces of data or commands. Set the delimiter as follows:

Menu Map item [94]	d-CR	Set to CR(0DH).
Menu Map item [95]	d-LF	Set to LF(0AH).
Menu Map item [96]	d-CR-LF	Set to CR+LF(0D0AH).
Menu Map item [97]	d-win	Transfers the data directly to Microsoft® Windows®. This is equivalent to pressing the Enter key of the personal computer.
Menu Map item [98]	d-win-	Transfers the data directly to Microsoft® Windows®. This is equivalent to pressing the right cursor key of the personal computer.

Note: d-win • d-win-

 Section 3.2.4 "Windows® Direct".

When these Menu Map items are set, it is not possible to send commands to the balance from the peripheral instruments. The personal computer and printer cannot be used at the same time.

3. Connecting Peripheral Instruments

A variety of peripheral instruments are available for use with the BW and BX series balance, such as a printer, keyboard or personal computer.

This chapter describes how to connect and use peripheral instruments.

3.1 Printer

Shimadzu Printer EP-100, EP-110 can be connected.

- 1) Use the default communication settings for the balance listed below.
 Menu [76] [77] [83] [89] [92] [94]
 H-tm F-Eb b-1200 P-no S-1 d-Cr
- 2) Disconnect the power to the balance and Printer.
- 3) Connect the DATA I/O connector of the balance to the Printer using the attached cable of the Printer.
- 4) First turn ON the power to the balance, then to the Printer.

3.2 Personal Computer - RS-232C -

3.2.1 Connecting the Cable



Caution

Signals other than RS-232C are also output by the BW and BX series balances through the RS-232C/AUX connector. If these signal lines are incorrectly connected, damaged may occur to the personal computer or balance. Correctly connect an appropriate cable for communication between the balance and personal computer.

All types of personal computers may not operate normally when connected with the optional RS-232C cable according to the diagram.

When using WindowsDirect Function, follow 3.2.4 for setting.

(1) For DOS/V computers (D-sub9 pin) (Null-modem)

Computer side			Balance side		
RXD	2	-----	2		TXD
TXD	3	-----	3		RXD
DTR	4	-----	6		DSR
SG	5	-----	7		SG
DSR	6	-----	20		DTR
RTS	7	┌	5	┐	CTS
CTS	8	└	4	┘	RTS

(2) IEEE Standard (D-sub 25-pin)

Personal computer side		Balance side
1	-----	1
2	-----	3
3	-----	2
4	-----	5
5	-----	4
6	-----	20
7	-----	7
20	-----	6

3.2.2 Data Format

The following explanation is applied when F-E is selected.

For other formats, refer to the data corresponding to the compatible machine.

Note: \square indicates space code and <delimiter> indicates delimiter code.

(1) For the measured value`

First character: Minus: '-' Non-minus: space

2nd to 11th characters: Numeric values or "[", "]" are flushed right. The position of the decimal point varies depending on the type of instrument.

12th to 13th characters: Unit such as g \square or kg

14th to 15th characters: Delimiter

Note: When the delimiter is d-C or d-LF , the 13th character is not present.

When printing stability information, the following characters are put in front of the first character mentioned above.

Stable time: S

Unstable time: D

When the delimiter is d-w or d-w , the special format is used instead of that described above.

(2) For of OL or -OL

OL $\square\square\square\square$ OL $\square\square\square$ <Delimiter>

-OL $\text{-}\square\square\square\square$ OL $\square\square\square$ <Delimiter>

3.2.3 Using Command Codes

Note: If communication conditions are incorrectly set, a communication error Error is displayed.

(1) Commands that end with a number, character, or symbol other than [=]: Transmit to the balance with a delimiter for each command code.

Example 1: PRINT<CR> ... The same operation as pressing the

PRINT
▶ key

(2) Commands that end with a [=]: Transmit the number to the balance with a delimiter.

Example 2: TIME=1234 <CR> .. 12:34 is set as the current time.

Example 3: P.TARE=1.23 <CR> (example of type of the second decimal place).

...1.23g is set as a Pretare value.

Example 4: P.TARE=0.00 <CR> (example of the second decimal place)

...Clears (cancels) the Pretare value.

Note: Number of digits, decimal point, position of decimal point of the numeral transmitted succeeding to '=' are the same as the case of setting the numeric value using the AKB-301 Application Keyboard.

Number of digits below the decimal point is as follows.

S-type and **H-type:** Use the same number of decimal places as in the gram-display.

This restriction does not apply to USER=, SOLID=, and LIQUID=.

Example 5: MENU=4630 <CR> (4-digit number after [=])

This is the same in the following operation.

... from mass display, press the $\frac{\text{CAL}}{\text{MENU}}$ key four times, the

press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key once and the press the $\frac{\text{CAL}}{\text{MENU}}$ key

six times. This returns to the same display as if the $\frac{\text{CAL}}{\text{MENU}}$ key was never pressed.

Press the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key once, the $\frac{\text{CAL}}{\text{MENU}}$ key three times, and

the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key once more.

This example indicates that $\overline{\text{EA}}-\overline{\text{B}}$ is selected.

Note: If there is 0 in the four-digit number, the setting is complete at that point and menu selection is ended.

The result of this command varies depending on the type of the balance.

Example 6: #=2.56 <CR>

Example 7: #=12.345.67 <CR>

A personal computer can instruct the weighing and display a specific number on the balance.

With the commands in Example 6 &7, [#2.56] and [#12.345.67] are displayed on the balance. When the

operator presses the $\frac{\text{PRINT}}{\blacktriangleright}$ key, the character string '2-56<CR>' and '12-345-67<CR>' are output from the balance.

(3) Echo back command

The balance again transmits the character strings of N pieces included between an echo back command '{' or '}' and the delimiter.

An unprocessed echo back command is not left in the receiving buffer of the balance, $N \leq 30$.

Example 8: ABCDEFG12345<CR>

... After receiving this command, the balance outputs ABCDEFG12345<CR>. The printer can print this character string.

Note: Only capital alphabets and a part of symbols (decimal point, symbol etc.) can be used when printing with a printer. A maximum of 15 characters per line.

(4) Command codes for F-EB and F-PEEB

(i) Commands related to output

D01	Continuous output
D03	Continuous output with stability information
D05	Single output
D06	Auto Print setting (type of Auto Print is set separately)
D07	Single output with stability information
D09	Release of continuous output and Auto Print

(ii) Commands related to operation keys

POWER	Equivalent to the $\frac{\text{POWER}}{\text{BRK}}$ key
Q	Equivalent to the $\frac{\text{POWER}}{\text{BRK}}$ key
MENU	Equivalent to the $\frac{\text{CAL}}{\text{MENU}}$ key
TARE	Equivalent to the $\frac{\Rightarrow 0/T \Leftarrow}{\downarrow}$ key
T	Equivalent to the $\frac{\Rightarrow 0/T \Leftarrow}{\downarrow}$ key
UNIT	Equivalent to the $\frac{\text{UNIT}}{\blacktriangle}$ key
PRINT	Equivalent to the $\frac{\text{PRINT}}{\blacktriangleright}$ key

POWER+	Equivalent to holding the $\frac{\text{POWER}}{\text{BRK}}$ key for approximately 3 seconds.
MENU+	Equivalent to holding the $\boxed{\text{MENU}}$ key for approximately 3 seconds.
UNIT+	Equivalent to holding the $\frac{\text{UNIT}}{\blacktriangle}$ key for approximately 3 seconds.
PRINT+	Equivalent to holding the $\frac{\text{PRINT}}{\blacktriangleright}$ key for approximately 3 seconds.
RECALC	Equivalent to the [RECALC] key of the AKB-301 Application Keyboard
C	Equivalent to the [C] key of the AKB-301 Application Keyboard

(iii) Commands related to Application Measurement

PEAK	Sets the Peak Hold mode.
AZERO	Sets the Auto Zero mode ON.
INTERVAL	Sets the Interval Timer mode.
MEMORY	Sets the Memory mode.
M	Immediately operates after setting the Memory mode,.
ADDON	Sets the Auto-Memory and Zeroing mode.
+	Immediately operates after setting the Auto-Memory and Zeroing mode,.
A	Sets the Animal Weighing mode.
ANIMAL	Sets the Animal Weighing mode.
R	Releases the Application weighing mode.

(iv) Commands related to unit conversion

g	Switches to "g" unit
kg	Registration of "kg" unit and switching
PERCENT	Registration of "%" unit and switching
%	Sets 100% when display is in "%" unit.
G	kg <- -> % switching
PCS	Registration of "PCS" unit and switching
CT	Registration of "karat" unit and switching
MOM	Registration of "monme" unit and switching

CU Switches to "user" unit (Set the conversion coefficient beforehand.)

RSTUNIT Returns the default units.

(v) Readout commands of set value

TARGET Readout of target set value

LIMIT Readout of limit set value

G.LO Readout of lower limit set value in Group mode

G.LO Readout of upper limit set value in Group mode

L.LO Readout of lower limit set value in Level mode

L.UP Readout of upper limit set value in Level mode

UW Readout of unit weight set value

G/PCS Equivalent to g/PCS key

CALWT Readout of external weights set value for span calibration

P.TARE Readout of Pretare set value

ZRNG Readout of Zero Range set value

USER Readout of user unit conversion coefficient

VOL Readout of reference weight set value

DENSE Readout of surrounding liquid density set value

I.TIME Readout of Interval Timer set value

Note: The command to read out the numerical value settings can only be given by the personal computer.

(vi) Commands for numeric value setting

CALWT= Sets external weights value for span calibration.

P.TARE= Sets Pretare value.

ZRNG= Sets Zero Range value.

UW= Sets unit weight.

USER= Sets user unit conversion coefficient.

VOL= Sets volume of reference weight.

I.TIME= Sets interval timer value.

DATE= Sets the date.

TIME= Sets the time.

TARGET= Sets the target value.

LIMIT= Sets the limit value.

G.LO= Sets the lower limit value of Group Display mode.

G.UP= Sets the upper limit value of Group Display mode.

L.LO=	Sets the lower limit value of Level Display mode.
L.UP=	Sets the upper limit value of Level Display mode.
PCS=	Sets the arbitrary loading piece.
#=	Corresponds to numeral keys of AKB-301 Application Keyboard.
ID=	Sets ID.
PASSSET=	Sets PCAL password.
PASS=	Inputs PCAL password.

(vii) Commands of special functions

CAL	Enters Span Calibration mode.
C18	Enters Span Calibration mode.
LOCK	Sets menu lock.
RELEASE	Releases menu lock.
TIME	Readout of date and time
RSTMN	Menu reset
MENU=	Sets arbitrary menu.
{	Echo back
}	Echo back

(5) Compatible commands with Mettler® Electronic Balances

S	One time output at a stable state
SI	Immediate one time output
SIR	Continuous output
SR	Continuous output at a stable state
T	Taring after stabilized
TI	Immediate taring
Z	Zero setting (same as immediate taring)

(6) Compatible command with Sartorius® Electronic Balances

<ESC>P	One time output
<ESC>T	Taring

Note: <ESC> indicates escape code (1BH).

3.2.4 WindowsDirect Function

3.2.4.1 Introduction: Experience it!

The BW-K / BX-K series balance can transfer data directly to a computer running MS-Excel, mass input window of analytical instrument software or other applications on Windows® OS, as if the displayed value were typed from the keyboard to the cursor position. This function is called WindowsDirect. As this function directly accesses the Windows® OS, communication software-installation troubles are eliminated. A cable and a few simple settings are all that is needed to enable data transmission from the balance.

For bi-directional communication between the balance and the computer, software is required. WindowsDirect does not send commands from the computer.

- Only numerical values can be transferred through WindowsDirect.

3.2.4.2 Set Up WindowsDirect

Simple settings are made for the balance and the computer. Connection is by RS-232C cable (3.2.4.2.2) specified by Shimadzu.

If bi-directional communication software is used: WindowsDirect function should be turned off. Set up the optimal communication parameters for the software according to "2.12 Communication Menu".


3.2.4.2.1 Setting Up the Balance

1. Supply power to the balance and set the condition for communication as follows.

Menu	[74]	[77]	[81]	[89]	[92]
	H-SOFT	F-Eb	b-300	P-no	S-1
Menu	[97]	[98]			
	d-w in	d-w in			


2. Go to "STAND-BY" by pressing the [POWER] key several times and unplug the AC adapter from the balance. Unplugging the balance once is necessary after the above setting.

3.2.4.2.2 Cable Connection

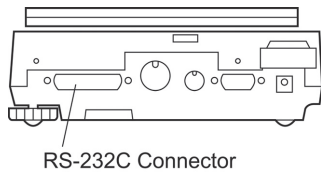

1. Verify the balance display is "STAND-BY".
2. Turn off the computer and remove power from the balance.
3. Connect the RS-232C cable to the balance.
4. Connect the RS-232C cable to the computer.

When using WindowsDirect, use a Null modem cable of one of the below wirings.

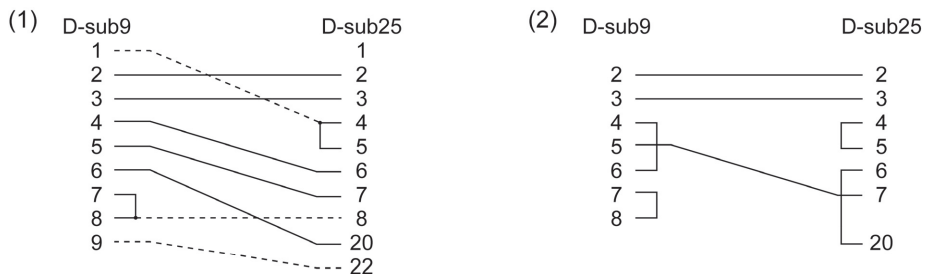


--- is not necessarily required.


A cable of the (1) wiring (including --- wiring) is available as an optional accessory.

RS-232C Cable 25P-9P (1.5m) P/N 321-60754-01

3.2.4.2.3 Setting Up the Computer (leave the balance unplugged)

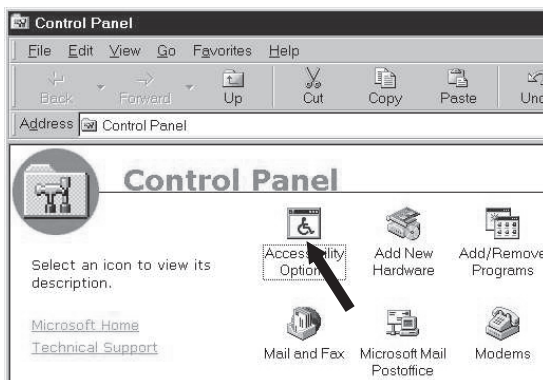


1. Turn ON the power to the computer and start Windows®.
2. Click "Start", choose "Settings", and "Control Panel".
3. Select "Accessibility Options."
4. Verify that there are no check marks for any items on all five tabs including "General."

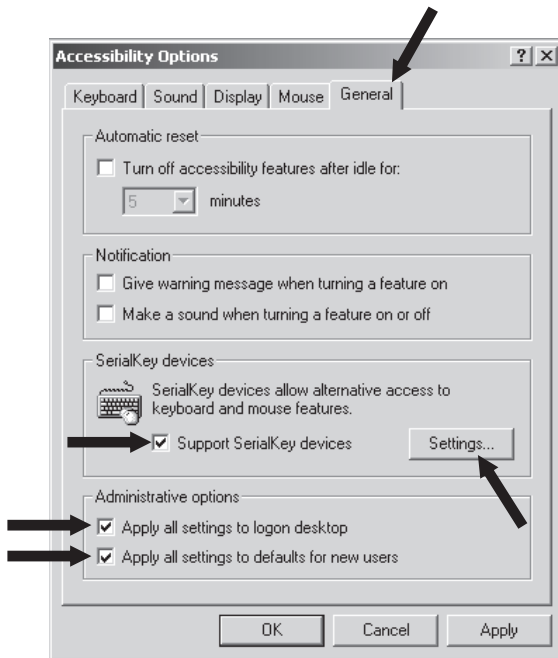

5. Put a check mark at "Support Serialkey device" in the "General" tab. This should be the only check mark on all the tabs of Accessibility Options unless "Administrative options" appears in the "General" tab. Put check marks at both the items of "Administrative options" to maintain the settings even after restarting Windows®.

6. Open "Settings".

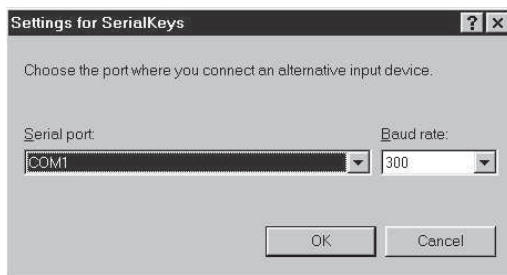


7. Select the serial port corresponding to the RS-232C port of your personal computer. (Serial port: any one of COM1 to 4. Usually, COM1)

8. Select a "Baud rate" of 300.

9. Click "OK".

10. Click "Apply" and wait.

11. Click "OK".

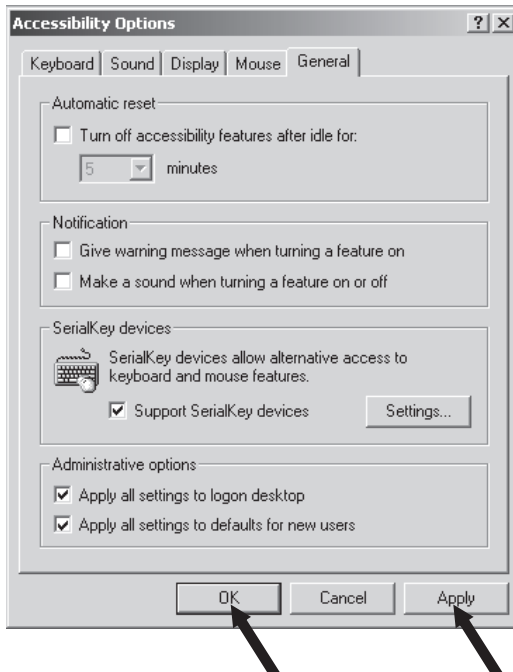

3.2.4.2.4 Start and Checking Operation

1. Confirm Windows® is free from any application.
2. After Windows® has completely started, connect power cord from the AC adapter to the balance, when "oFF" is displayed, press the [POWER] key. The mass display appears.

Note: Turning ON the balance before Windows®* is completely activated may cause incorrect operation.

3. Open the "Note pad" accessory in Windows®* (or start the application you wish to use).
4. Press the [PRINT] key of the balance.
Verify that the numeric value displayed on the balance appears at the cursor position on the screen of computer. The effect is the same as typing the value from the computer keyboard and pressing the ENTER key. Characteristics indicating the unit of measure are not sent to the computer.

Note: To output date and time from the balance, press and hold the [PRINT] key for about 3 seconds.

5. Test combination with Auto Print function, if you wish to use it.
(Refer to 2.9.4)
6. End the operation using the standard close or exit procedure.

3.2.4.3 Troubleshooting the WindowsDirect Communication Function

If the WindowsDirect communication function doesn't run properly, check the following points.

If this doesn't resolve the problem, contact your Shimadzu representative.

Q1 WindowsDirect communication has been set but it is not operating at all.

A1

- Check the type of communications cable used for the connection (Shimadzu authorized part or another part available on the general market) and the soundness of the connection.
- If a USB serial converter is used, depending on the circumstances at the setup there is a possibility that it has been automatically set to a COM port number higher than 4, and in this case you should reassign it to a COM port number that can be used by serial key devices (COM1 to COM4).
- It is possible that the driver used as an accessory with the USB serial converter has not been set up properly. Try uninstalling the driver and installing it again.
- Some notebook PCs feature a setting for disabling RS-232C ports as a power-saving measure. Before trying to use the WindowsDirect communication function, make the setting that enables the use of RS-232C ports.

- Communications with other applications and PCs via a LAN may interfere with the serial key device settings. Try using WindowsDirect communication without using the LAN.

Q2 The WindowsDirect communication function won't work after I restart the PC.

A2

- Some PCs don't recognize that a serial key device has been set when they start up. For details on how to deal with this, contact your Shimadzu representative.

Q3 I want to use the WindowsDirect communication function with Windows Vista or later OS.

A3

- Windows Vista or later OS doesn't have the serial device setting screen that is required to set the WindowsDirect communication function. For details on the setting, contact your Shimadzu representative.

Q4 Data is input to the PC as garbled characters.

A4

- Either the balance or the PC is not set for the WindowsDirect connection function. Make the settings again.

Q5 When data is input into Excel, the cursor doesn't move to another cell.

A5

- If a function for conversion to 2-byte characters is available in Windows, turn the setting for this function off.
- Click the [Edit] tab under [Options] in Excel and check [Move selection after Enter] (if cells move in response to keyboard input there is no problem).
- Check the input data in another application (e.g. Notepad).

Q6 The operation is sometimes abnormal.


A6

- Depending on the processing capability of the PC, malfunctions may occur if the communications speed is high. Set 300 bps for the communication speed. Malfunctions may also occur if the interval for data transmission from the balance is too short. Ensure that one batch of data is displayed on the screen before the next batch of data is sent. And if there is limited processing capability, don't use the continuous output function.
- When data is sent from the balance, don't touch the PC's keyboard or mouse.

3.3 Optional AKB-301 Application Keyboard

The optional AKB-301 Application Keyboard is provides numeric keys which are very convenient for inputting numbers. It is possible to perform various numeric settings such as, piece count and threshold value of graphic display (comparator), as if you were operating a pocket calculator.

Use of the AKB-301 Application Keyboard is recommended for performing piece count and pass or failure judgment in daily operation.

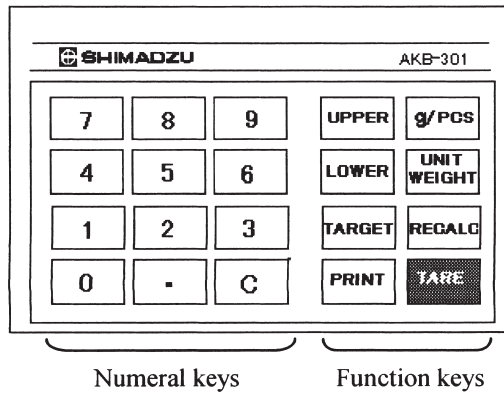

3.3.1 Basic Information

3.3.1.1 Connection

- 1) Disconnect the power to the balance.
- 2) Connect the AKB-301 cable to the KEY connector on the back of the balance.
- 3) Turn on the power to the balance.

3.3.1.2 Overview of Use

- 1) To set numeric values ➡ Press a function keys after the number keys.
 - 2) To check the set numeric values ➡ Press only a function key, not a number key.
 - 3) To return to the mass display ➡ Press the C key.
- If numeric keys other than C are pressed, the # symbol illuminates to indicate that a number is being entered.

3.3.1.3 Summary of Function Keys

Function key	Function	
	When this key is pressed after the numeric keys	If this key is pressed while the numeric keys are not pressed *1
UPPER key *2	Sets UPPER value of graphic display.	Displays UPPER value of graphic display.
LOWER key *3	Sets LOWER value of graphic display or sets L m t value.	Displays LOWER value of graphic display or displays L m t value.
TARGET key *4	For target display, sets TARGET value.	For target display, displays TARGET value.
g/PCS key	Recalculates the weight with regard to the present load as the input value (pieces) and saves.	Switches gram-display unit and displays in piece unit.
UNIT WEIGHT key	Stores the numeric values entered as the unit weight.	Displays the unit weight presently stored.
RECALC key	Nothing is done.	When piece is displayed, divides the present load by the present displayed piece count to recalculate the unit weight.
PRINT key *5	Outputs the input numerals.	Outputs the displayed value.
TARE key *5	Sets Pretare value.	Performs taring.

*1 When these values are displayed, the * and # symbols are illuminated on the screen.

Pressing the C key returns these displays to the mass display.

These displays can be output (printed) with the PRINT key.

*2 This key functions only when a comparator is used in Group or Level mode. ERR is displayed for other modes.

*3 When a comparator is used in Group or Level mode, LOWER is the target value. L m t is the target value in the Target mode. ERR is displayed for other modes.

*4 This operation is performed only in Target mode. $\Xi \text{---} \text{---}$ is displayed for other modes.

*5 Operation of the TARE and PRINT keys are the same as keys on the main balance body.

3.3.2 Example of Use

3.3.2.1 Using with a Comparator

Note: The decimal key of the AKB is not used for setting comparator values. The decimal place is set automatically according to the display capabilities of the balance.

Preparation

Set the kind of graphic display you desire to use with menu selection.

1) From the mass display, press the $\frac{\text{CAL}}{\text{MENU}}$ key three times and the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key once.

2) Use the $\frac{\text{CAL}}{\text{MENU}}$ and $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ keys to set the desired graphic display mode.

(1) Target display mode

Setting the target value

Example: Press the numeric keys for 10000 and then the [TARGET]key.

➡ The Target value is set to 10000.

Setting the limit value

Example: Press the numeric keys for 5000 and then the [LOWER] key.

➡ The Limit value is set to 5000.

Checking the target value

Press the [TARGET] key to display the set value.

Checking the limit value

Press the [LOWER] key to display the set value.

After checking, press the g/PCS key to return to the mass display.

Note: $\Xi \text{---} \text{---}$ is displayed if the [TARGET] key is pressed without being in the Target display mode.

(2) In Comparator mode (group display and level display)

Setting the upper limit value

Example: Press the numeric keys for 20000 and then the [UPPER] key.

➡ The upper limit value is set to 20000.

Setting the lower limit value

Example: Press the numeric keys for 10000 and then the [LOWER] key.

➡ The lower limit value is set to 10000.

Press the [UPPER] key to display the set value.

Press the [UPPER] key to display the set value.

Press the [LOWER] key to display the set value.

Press the [LOWER] key to display the set value.

After checking, press the g/PCS key to return to the mass display.

Use the [g/PCS], [UNITWT], and [RECALC] keys.

Use the following procedure to count pieces based on a sample with an arbitrary number of pieces. For example, to recalculate the unit weight based on 25 pieces and perform piece counting.

- 1) Load the container and press [TARE] key or zero the balance.
- 2) Count 25 pieces and load them on the pan.

Note: If in gram-display, the gram value of the 25 piece sample is displayed. If in PCS display, the value is converted into piece display using the previously stored unit weight. The display may not be 25 pieces.

- 3) Press the numeric keys for 25 and then the [g/PCS] key.
- 4) Piece count is now displayed in the new unit weight.

To recalculate the unit weight with the currently displayed piece count, press the [RECALC] key when the correct piece count is displayed.

To input a unit weight directly, (for example, 1.234). Press 1.234 and then press the [UNIT WEIGHT] key. Set the unit weight in "g". A value using the maximum number of digits displayed by the balance can be set. Use the decimal point key on the AKB when entering this value.

Press the [UNIT WEIGHT] key.

Press the [g/PCS] key to switch between the "g" unit and "piece count" unit.

Setting the Pretare Value

Example: Enter 1.234 and press the [TARE] key.

1.234 g is set as the Pretare value.

Note: Number of decimal places is set as follows:

H-type and S-type ... with number of decimal places at 1d gram-display unit

If not correct, Ξ_{rr} is displayed.

Example: To cancel the Pretare value for a balance with three decimal places:

X wrong example: 0 [TARE]

O correct example: 0.00 [TARE]

Checking the Pretare value

Press and hold the [TARE] key for approximately 3 seconds. Press the [TARE] key again to return to the mass display.

Cancel a Pretare Value

Cancel a pretare value by setting the value to zero using the correct decimal placement for the balance.

Output Keyboard Numbers to Peripheral Devices

Arbitrary numbers with a maximum of seven digits can be output using the numeric keys and the [PRINT] key.

The sample number can be input at the balance and transferred with the balance output to the personal computer.

Example: To input 123-456:

Enter 123.456 and press the [PRINT] key.

Setting a Numeric Menu Value

Use the number keys of the AKB to enter numeric values during menu selection.

Numbers are set sequentially starting from the blinking digit. The decimal point can be set using the decimal point on the AKB.



Last Menu Recall function with AKB-301

Last Menu Recall function is helpful when changing the numeric values for the same

Menu Map item. To use this function, press and hold the $\frac{\text{CAL}}{\text{MENU}}$ key until the menu is displayed.

3.4 Foot Switches

A foot switch is convenient when the operator must work with both hands. Two types of foot switches are available.

FSB-102T has the same operation as the $\frac{\rightarrow 0/T \leftarrow}{\downarrow}$ key of the balance

FSB-102P has the same operation as the $\frac{\text{PRINT}}{\blacktriangleright}$ key of the balance

Note: Either one of the foot switches can be connected at a time. The foot switches cannot be used together with the AKB-301 Application Keyboard.

3.5 Hook for Below-Balance Weighing

This hook is useful for conducting measurements with the sample suspended beneath the balance, such as in specific gravity measurement.


Attachment method

Attach the below-balance weighing fitting by screwing the supplied screws into the screw holes for the transportation screws on the bottom of the balance (2 places), as shown in the diagram on the right, until the screws cannot be tightened anymore.

Note:

Be careful to ensure that no undue force is applied to the fitting when putting down the balance while the fitting is attached.

Otherwise, the balance may not operate correctly. In particular, do not put down the balance in a location where there is no hole for below-balance weighing.

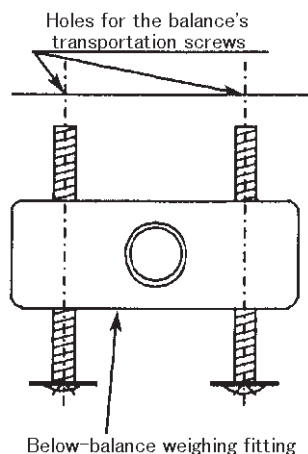

4. APPENDIX

4.1 Part List

4.1.1 Standard Accessories and Maintenance Parts

Name	Part No.	Remarks
Pan ASSY, KW	S321-41074	For the BW-K
Pan ASSY	S321-40919-01	For the BX-K
Pan support cap	S321-40965-04	
Vinyl cover	S321-40924	
Vinyl cover with double-sided tape	S321-73267-01	
Double-sided tape for fixing the vinyl cover	S320-02138	
Calibration weight stopper screw	S321-41046	For the BW-K
Calibration mass cover setscrew (SUS bind M4x8)	S020-37512	For the BW-K
Transportation screw (SUS Sems P3BK M4x16)	S020-46551	
Caster ASSY	S321-30070-03	

4.1.2 Options

Item	Part No.	Description
EP-100	S321-73900-11	
EP-110	S321-73900-12	
AKB-301 Application Keyboard	S321-53382-01	For single/multiple value setting, etc. in piece counting mode. For setting thresholds in pass/fail evaluation. For setting other numerical values.
FSB-102TK Foot Switch	S321-60110-12(D9)	For "TARE"
FSB-102PK Foot Switch	S321-60110-11(D9)	For "PRINT"
Hook for Below-Balance Weighing	S321-40962(D9)	This hook is useful for conducting measurements with the sample suspended beneath the balance, such as in specific gravity measurement.
RS-232C Cable (1.5m)	S321-60754-01(D9)	
RS-232C Cable (3.0m)	S321-60754-02(D9)	
RS-232C Cable (4.5m)	S321-60754-03(D9)	
USB-Serial Conversion Kit	S321-62520-01(D9)	The item includes an RS-232C cable.

Note:

- Part number and specification are subject to change without notice.
- It is not guaranteed that RS-232C cable will conform to all computers.

4.2 Table of Unit Conversion Constants

SCROLL ORDER	MENU	UNIT	Conversion coefficient (g=)	Note (display)
[54]	U-kg	kg	0.1	
[55]	U-mg	mg	1000	
[56]	U-%	%		
[57]	U-pcs	Pcs		
[58]	U-ct	ct	5	
[59]	U-mom	Mom	0.266667	
[60]	U-□d	DS		
[61]	U-d	DL		
[61a]	U-Lb	Lb	0.00220462	
[61b]	U-Oz	Oz	0.0352740	
[61c]	U-Ozt	Ozt	0.0321507	
[61d]	U-HK	TL-HK	0.0267165	A triangle symbol displays
[61e]	U-HK	TL-HK (Jwl)	0.0267173	Three triangle symbols display
[61f]	U-SporE	TL-S'pore	0.0264554	
[61g]	U-tiwAn	TL-Taiwan	0.0266667	A triangle symbol displays
[61h]	U-tiwAn	TL-Taiwan	0.0266667	Three triangle symbols display
[61i]	U-mal	TL-Malaysia	0.0264600	
[61j]	U-ChinA	TL-China	0.0266071	A triangle symbol displays
[61k]	U-ChinA	TL-China	0.0266071	Three triangle symbols display
[61l]	U-dwt	dwt	0.643015	
[61m]	U-GN	GN	15.4324	
[61n]	U-t	t	0.0857339	
62	U-USEr			Set by user

4.3 Specifications for the RS-232C/AUX Connector

Pin number	Name	Function	Remarks
1	FG	Frame ground	
2	TXD	Data output	
3	RXD	Data input	
4	RTS	Internal connection with CTS	
5	CTS	Internal connection with RTS	
6	DSR	Handshake (receiving)	
7	SG	Signal grounding	
8	NC	Blank	
9	NC	Blank	
10	NC	Blank	
11	NC	Blank	
12	NC	Blank	
13	NC	Blank	
14	NC	Blank	
15	NC	Blank	
16	NC	Blank	
17	NC	Blank	
18	NC	Blank	
19	NC	Blank	
20	DTR	Handshake (receiving)	
21	NC	Blank	
22	NC	Blank	
23	NC	Blank	
24	NC	Blank	
25	NC	Blank	


Note: NC= No connection within the balance.

4.4 Error Display List

4.4.1 General Display

Display	Description of message
---	Wait for next display.
- 10d-	Minimum display resolution is decreased by one decimal place.
- 1d-	Minimum display digit is returned to original state.
-t ime-	Date and time are being output.
Abort	Operation was aborted.
APL End	Application Measurement was released.
d over	Calibration check detects too large error.
d Under	Calibration check detects too large error.
LoCKEd	Menu lock is applied.
rELEASE	Menu lock is released.
rESEt	Menu was reset.
SEt	Contents of new setting and coefficient were stored.
oFF	Reset by power failure.
wA it	Built-in weight is moving. Wait.
All numerals blinking	Place the displayed calibration mass.

4.4.2 Error Display

Error display	Description	Countermeasure
CAL E1	The load on the pan is unstable at calibration. Only one of the calibration levers is working.	Avoid wind and vibration.
CAL E2	The drift of zero point is large at calibration.	Empty the pan.
CAL E3	The drift is large at the time of PCAL.	Use correct weight.
CAL E4	The drift is large at span calibration.	Use correct weight.
CAL E5	Weight is wrong.	Use correct weight.
CHE x	Stops when this is displayed.	*
ComErr	Received command code is not correct.	Check delimiter etc.
dSP oL	Integer of the displayed unit exceeded 7 digits.	Decrease the load.
Err 0x	Failure in the balance.	*
Err 10	PCAL password error. “—” is displayed after the error.	Check the password. Press the <u>POWER</u> BRK key and return the balance to mass display mode.
Err 20	Attempted to set improper numeric value.	Review the numeric value and decimal places.
Err 21	Necessary conditions and numeric values are not set.	Check graphic display mode.
Err 22	 key was pressed during menu lock.	Release menu lock.
Err 23	Mode for operation is different.	Check graphic display mode.
Err 24	Cannot be stored, as the supply voltage is abnormal.	Check the supply voltage.

*Contact your Shimadzu representative.

**TOKYO OFFICE**

3, Kanda-Nishikicho 1-chome, Chiyoda-ku, Tokyo 101-8448, Japan
Phone: 81(3)3219-5641 Fax: 81(3)3219-5710
Cable Add.:SHIMADZU TOKYO
Overseas Telex No.: 0232-3291 (SHMDT J)

KYOTO OFFICE

1, Nishinokyo-Kuwabaracho, Nakagyo-ku, Kyoto 604-8511, Japan
Phone: 81(75)823-1200 Fax: 81(75)812-3438
Cable Add.:SHIMADZU KYOTO
Overseas Telex No.: 05422-166 (SHMDS J)

SHIMADZU(CHINA)CO.,LTD. PUXI BRANCH

Block E, No.570 West Huaihai Road, Shanghai, 200052 P.R. China
Phone: 86(21)2201-3888 FAX: 86(21)2201-3666

SHIMADZU SCIENTIFIC INSTRUMENTS, INC.

7102, Riverwood Drive, Columbia, Maryland 21046, U.S.A
Phone: 1(410)381-1227 Fax: 1(410)381-1222

SHIMADZU EUROPA GmbH

Albert-Hahn-Strasse 6-10, D-47269 Duisburg, F.R. Germany
Phone: 49(203)7687-0 Fax: 49(203)7666-25

SHIMADZU MIDDLE EAST & AFRICA FZE UAE HEAD OFFICE

Warehouse No. RA08UC02, Jebel Ali, Dubai, P.O. Box 262081 United Arab Emirates
Phone: 971(4)883-6668 FAX: 971(4)883-6808

SHIMADZU MIDDLE EAST & AFRICA FZE ISTANBUL BRANCH

Kosuyolu Mah., Salih Omurtak Cad., No:57
34718, Kadikoy Istanbul, Turkey
Phone: 90(216)545-8940 FAX: 90(216)545-8941

SHIMADZU (ASIA PACIFIC) PTE LTD.

79 Science Park Drive, #02-01/08 Cintech IV, Singapore Science Park I, Singapore 118264
Phone: 65(6778)6280 Fax: 65(6779)2935

SHIMADZU DO BRASIL COMERCIO LTDA.

Avenida Marques de Sao Vicente, 1771-Barra Funda,
CEP 01139-003, Sao Paulo, SP, Brazil
Phone: 55(11)2134-1688 Fax: 55(11)3611-1635

SHIMADZU SCIENTIFIC INSTRUMENTS (OCEANIA) PTY. LTD.

Unit F, 10-16 South Street Rydalmere N.S.W. 2116
Phone: 61(2)9684-4200 Fax: 61(2)9684-4055