High Performance Packed Column

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
Shim-pack Guard column SCR(H)

Instruction Manual

- “Guard column SCR(N)” and “Guard column SCR(H)” are guard columns (4 mm I.D., 50 mm L.) packed with a cation exchange resin consisting of styrene-divinylbenzene copolymer.
- Guard column SCR(N) is exclusively used in combination with Shim-pack SCR-101N separation column. Refer to the manuals of Shim-pack SCR-101N as well.
- Guard column SCR(H) is exclusively used in combination with Shim-pack SCR-101H separation column. Refer to the manuals of Shim-pack SCR-101H as well.
- The tubing kit available for column connection is listed below.

Item Name	P/N	Comment
Column connector kit	228-17943-92	Ferrule 1.6 F x 2 Male nut 1.6 MN x 2

- When installing the guard column, please refer to the diagram below to connect the joints to both ends of the column. Be careful not to over-tighten, as excessive tightening may lead to damage at the connection part.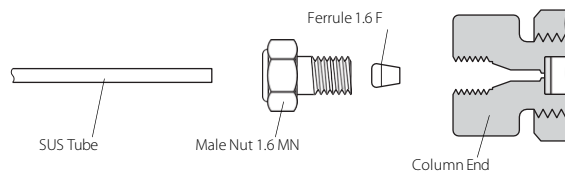

Column Connection diagram

■ Disposal Precautions

Separate the column from general industrial or household waste and dispose of it in accordance with local government regulations.

■ Technical Support

If you encounter any performance defects, please contact your local Shimadzu sales office for assistance. However, please note that we cannot guarantee the lifetime of the columns, nor can we accept any claims when performance has degraded due to non-compliance with the operation procedures outlined above, or as a result of normal aging.

*Please note that the appearance, specifications, and the contents of this instruction sheet are subject to change without notice.