High Performance Packed Column for HPLC

CoreFocus

Shim-pack™ SR-C30 Guard column (for analysis)

INSTRUCTION MANUAL

■ Introduction

Shim-pack SR-C30 Guard column (for analysis) is cartridge-type guard column filled with high pressure packing materials. Please read the following instructions before use.

■ Operating Precautions

Check if anything is missing or damaged. If there are any signs of damage, notify your local Shimadzu representative at once.

■ Specifications


The product specifications are as follows.

Size	Particle size	Product name	Chemical bonding group
I.D. : 4 mm Length : 10 mm, 20 mm	3 μ m	Shim-pack SR-C30 (G)	Triacetyl (C30)
	5 μ m		

■ Installing the Column

Please note the following when you install the cartridge column.

- The optional cartridge holder is required when you install the column.
- The cartridge column can be replaced by hand.
- Do not remove the black end caps at both ends of the cartridge column.
- The cartridge holder can be used repeatedly. In case of degradation, replace the cartridge column only.
- The cartridge holder and the column coupler can be removed with a wrench (1/4 inch).

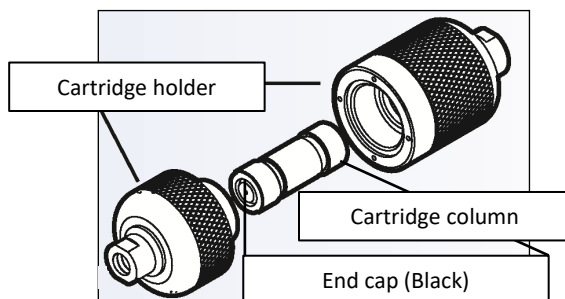■ Use Method

- 1) Loosen the cartridge holder by hand and place the cartridge column into the cartridge holder.
- 2) Tighten the cartridge holder by hand. Then additionally tighten it up to 30° using 1/4-inch wrench. At this time, normally, there is a little gap between the cartridge holders (figure below). Tightening more than requirement may cause damage.
- 3) Connect the assembled guard cartridge column in front of the main-column (the injector side) using tools such as a column coupler. Use a short tube to avoid diffusion of the injection sample solution.

Cartridge holder

■ Safety Guideline

- Select the cartridge columns whose functional groups are basically the same as those of the main column.
- The cartridge holder is sold separately. Select the cartridge holder according to the length of the cartridge column. Use cartridge holder set for Length 10 mm (P/N: 227-31172-03) / 20 mm (P/N: 227-31172-04).
- If the cartridge column is attached and detached repeatedly, the airtightness may be lost, and fluid may leak.
- The timing of replacement of the cartridge column should be determined by the increase of pressure, the change of peak shapes and the separations, etc.
- There is no specified direction for connecting cartridge columns, but do not change the flow direction of a cartridge column once it has been used. Reversing the flow may cause insoluble material adsorbed at the cartridge inlet to be carried into the main column.
- If you inject a large amount of a sample repeatedly, it may cause impurities to be adsorbed in the main column. Therefore, it may deteriorate the main column. Early replacement of the cartridge column is recommended to prevent contamination of main column.

- Please use tubing that has the same or smaller diameter as the autosampler outlet for the connection between the guard column and the analytical column.

■ Disposal Precautions

When disposing of the column, do so in accordance with the processing standards determined by law, separately from general industrial waste and household garbage.

■ Technical Support

Shim-pack SR-C30 series columns are manufactured, inspected, packaged and shipped under strict standards of quality control. Should you find any defect in performance, please contact your local Shimadzu representative, who will ensure your complete satisfaction.

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